

Feb 25, 2023

Measuring neutral lipids in fixed diatom cells using BODIPY 505/515

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l2427jg1y/v1>

Phoebe Argyle^{1,2}, Jana Hinners³, Nathan G. Walworth⁴, Sinéad ollins⁵, Naomi M. Levine⁴, Martina A. Doblin^{1,6}

¹Climate Change Cluster, University of Technology Sydney; ²Ministry of Marine Resources, Cook Islands;

³Institute of Coastal Ocean Dynamics, Helmholtz-Zentrum Hereon, 21502, Geesthacht, Germany;

⁴Department of Biological Sciences, University of Southern California, Los Angeles, CA, 90089-0371, USA;

⁵Institute of Evolutionary Biology, University of Edinburgh, Edinburgh, EH9 3JF, UK;

⁶Sydney Institute of Marine Science, Mosman, NSW, 2088, Australia



Phoebe Argyle

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l2427jg1y/v1>

Protocol Citation: Phoebe Argyle, Jana Hinners, Nathan G. Walworth, Sinéad ollins, Naomi M. Levine, Martina A. Doblin 2023. Measuring neutral lipids in fixed diatom cells using BODIPY 505/515. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.4r3l2427jg1y/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 30, 2020

Last Modified: February 25, 2023

Protocol Integer ID: 45067

Keywords: centric diatom thalassiosira spp, diatom cell, fixed diatom cell, measuring neutral lipid, phenotypic traits of microalgae, diatom plasticity, neutral lipid, frontiers in microbiology, throughput assay, microalgae, quantifying phenotypic trait, microbiology, cell

Funders Acknowledgements:

Gordon and Betty Moore Foundation Marine Microbes Initiative

Grant ID: MMI 7397

Abstract

This protocol is for the measuring neutral lipids in fixed diatom cells using BODIPY 505/515, as developed on the centric diatom *Thalassiosira* spp.

This protocol has been used in the following publications:

Argyle, P. A., Walworth, N. G., Hinnners, J., Collins, S., Levine, N. M., & Doblin, M. A. (2021). Multivariate trait analysis reveals diatom plasticity constrained to a reduced set of biological axes. *ISME Communications*, 1(1), 59.

Argyle, P. A., Hinnners, J., Walworth, N. G., Collins, S., Levine, N. M., & Doblin, M. A. (2021). A high-throughput assay for quantifying phenotypic traits of microalgae. *Frontiers in microbiology*, 12, 706235.

Materials



BODIPY[®] 505/515 (4,4-Difluoro-1,3,5,7-Tetramethyl-4-Bora-3a,4a-Diaza-s-Indacene) Thermo Fisher Catalog #D3921



Paraformaldehyde fixative: 4% paraformaldehyde in phosphate buffered saline (PBS)



Protocol materials



BODIPY[®]; 505/515 (4,4-Difluoro-1,3,5,7-Tetramethyl-4-Bora-3a,4a-Diaza-s-Indacene) **Thermo Fisher Catalog #D3921**



Paraformaldehyde fixative: 4% paraformaldehyde in phosphate buffered saline (PBS)



Liquid nitrogen



DMSO **Merck MilliporeSigma (Sigma-Aldrich) Catalog #D1435**



Preparation of BODIPY stock solution

- 1 Add  2 mg of  BODIPY[®]; 505/515 (4,4-Difluoro-1,3,5,7-Tetramethyl-4-Bora-3a,4a-Diaza-s-Indacene) **Thermo Fisher Catalog #D3921** in powder form to  1 mL of  DMSO **Merck MilliporeSigma (Sigma-Aldrich) Catalog #D1435** to create the stock solution of  2 mg/mL .

Store this in dark glass at  -20 °C to prevent degradation of the fluorescent dye.

Diatom sample preparation

10m

- 2 Sample the diatom culture of interest. For 'tube mode' flow cytometry take minimum  500 μ L into an eppendorf tube or other flow cytometry tube. For plate mode take  200 μ L into a round-bottom flow cytometry plate.

- 3 If using live cells, go to step 6

If using fixed cells:

Add  8 % volume

 Paraformaldehyde fixative: 4% paraformaldehyde in phosphate buffered saline (PBS)

at  10 % (v/v) to reach a final concentration of approx  0.8 % volume . For example, add  100 μ L of

 Paraformaldehyde fixative: 4% paraformaldehyde in phosphate buffered saline (PBS)

to  1000 μ L of microalgae sample.

- 4 Agitate to mix, then leave to fix for  00:10:00 to ensure fixation

10m



- 5 If samples are to be analyzed at a later date, samples should be stored in the fridge (max 48 hours), or flash frozen in  Liquid nitrogen then stored at  -80 °C .

Flow cytometry analysis

- 6 Measure the background fluorescence of unstained cells using a flow cytometer. This protocol was developed using a CytoFlex LX.

Equipment

CytoFLEX LX

NAME

Flow cytometer

TYPE

Beckman Coulter

BRAND

C40312

SKU

<https://www.beckman.com>

LINK

CytoFLEX LX N3-V5-B3-Y5-R3-I2 Flow Cytometer (21 Detectors, 6 Lasers)

SPECIFICATIONS

Fluorescence is measured using 488 nm excitation and 525/40 nm detection. Measure the background fluorescence of at least 200 cells but ideally 2000 or more.

- 7 Add the BODIPY stain in a  1 % (v/v) to final concentration of approx  0.002 mg/mL . For example add  2 µL to  200 µL of microalgae sample, or  2.2 µL to a fixed microalgae sample ( 200 µL sample with  20 µL



Paraformaldehyde fixative: 4% paraformaldehyde in phosphate buffered saline (PBS)

. Remember to take into account volume removed due to flow cytometry analysis.

- 8 Re-read the sample on the flow cytometer using the same laser parameters. Record the median fluorescence on the x channel for at least 200 cells



- 9 Calculate the change in median fluorescence between the stained and unstained sample. This is indicative of the level of neutral lipids present in the cell
- 10 If comparing between taxa of different sizes, size correction may be advisable. This may be done by dividing the change in median fluorescence by the equivalent spherical size, approximated from forward scatter which is measured in tandem during the flow cytometry analysis.