

Jan 15, 2025

Version 2

Meat preparation for meat identification V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov1961plr2/v2>

openbioscience Adrian¹

¹OBS



openbioscience Adrian

OBS

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1961plr2/v2>

Protocol Citation: openbioscience Adrian 2025. Meat preparation for meat identification. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.ewov1961plr2/v2> Version created by **openbioscience Adrian**

Manuscript citation:

Based on <https://www.ncbi.nlm.nih.gov/pmc/articles/PMC3550954/>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: November 28, 2024

Last Modified: January 15, 2025

Protocol Integer ID: 113108

Keywords: alkaline, LAMP, pork, identification, DNA, pork dna, absence of pork dna, meat preparation for meat identification, meat identification, alkaline method for dna extraction, meat sample, dna extraction, meat preparation, identifying unique dna sequence, lamp method for dna amplification, dna amplification, dna, unique dna sequence, identifying species, alkaline method, extraction, detection

Disclaimer

This protocol is work in progress for now.

Abstract

Identifying species by identifying unique DNA sequences is common practice.

In this case we will identify the presence or absence of pork DNA in a meat sample.

We will use the alkaline method for DNA extraction and LAMP method for DNA amplification and detection.

This is not a new protocol but rather the detailing of a previously published protocol listed in the manuscript citation.

Guidelines

Follow the country specific guidelines.

Materials

Consumables: 15 ml falcon tube, 1.5 ml centrifuge tube or equivalent, 1 PCR tube 0.1 mL PCR tubes with flat caps, commonly known as low-profile PCR tubes that matches the device well size.

Labware: graduated cylinder

Devices: Mortar and pestle, LAMP/PCR/Water bath and thermometer, scale, 1- 10 µl pipette, 100- 1000 µl pipette and tips. Optional: vortex, LAMP or PCR device or water bath, scale and boat.

Reagents: NaOH, Tris, HCl, DNase free water, meat sample to be tested

Others: stir rod, spatula, nitrile gloves, pipette rack , tube rack, labeling pen, tube rack.



Safety warnings

- ! Use appropriate PPE and follow Safe Work Practices.
NaOH (sodium hydroxide or caustic soda) needs to be handled with gloves because it's a highly corrosive alkali.
Seek and follow Safety Data Sheets for NaOH and HCl

Before start

Pork meat identification protocol. Meat Preparation protocol based on

<https://www.ncbi.nlm.nih.gov/pmc/articles/PMC3550954/>

"About 500 mg of meat was triturated in pestle & mortar with four volumes (4 ml) of 0.2N NaOH. 5 µl of the extract was again mixed with eight volumes (40 µl) of 0.2N NaOH in a sterile micro centrifuge tube. Mix was heated at 75 °C in water bath for 20 min and then added eight volumes (360 µl) of 0.04 M Tris HCl (pH 7.75) for neutralization of pH. 1 µl of final mix containing about 100 ng DNA was used for."



- 1 Prepare 10 ml of NaOH 0.2 mol/L (sodium hydroxide or caustic soda). This will be used twice, once for 4 mL and once for 40µL.
- 1.1 **Label a new 15 ml falcon tube "NaOH 0.2 mol/L" and add 10 mL DNAfree water and 0.080 g of NaOH. Mix to dissolve,** manually or using a vortex.

*For general molarity calculations see **Molarity calculation**. Molar concentration is the number of moles contained in one Liter of a substance. The concentration of a solution is expressed in mol/L $C=(m/M)/V$ where V is the volume in liters, m is mass in grams and M is molar weight? in grams/mol*

<https://www.alloprof.qc.ca/en/students/vl/sciences/calculating-molar-concentration-s1622>

So:

Molecular Weight of sodium hydroxide (NaOH) =40 g/mol: Na=22.989g/mol

O=15.999g/mol H=1.008g/mol ([http://chembook.org/page-nonav.php?chnum=1\\$=5](http://chembook.org/page-nonav.php?chnum=1$=5))

so: 22.989g/mol+15.999g/mol+1.008g/mol=39.996g/mol so about 40 g/mol.

so $m=C*M*V=0.2\text{mol/L}*40\text{g/mol}*10\text{ mL}=8\text{ g/L}*10\text{ mL}=80\text{ mg of NaOH}$

Meat Preparation

- 2 **Add 0.5 g fresh pork meat without fat in a pestle and mortar and add 4 ml NaOH from the tube labelled "NaOH 0.2 mol/L".**
- 3 **triturate (grind) the mix.** This will break the structure and results in a paste/liquid.
- 4 **Aliquot 5 µL from the mix and add it to a new sterile PCR (100 µl) tube labeled "prepared sample".** This is done with a 1- 10 µl pipette, preferably using filtered tips.

Meat Preparation

- 5 Optionally filter the result from previous step using a syringe and Luer filter.
- 6 **Add 40µL from the tube labelled "NaOH 0.2 mol/L" to the "prepared sample" tube.**



- 7 **Heat the resulting mix (45 µL) at 75 Celsius for 20 minutes** in either water bath, LAMP or PCR device.
- 8 **Move the content of the "prepared sample" to a new 1.5 ml centrifuge tube labeled final prepared sample and add 360 µl of 0.04 mol/L Tris HCl (pH 7.75)** prepared using the <https://www.protocols.io/view/prepare-tris-buffer-for-meat-identification-protocol-dndu5a6w> protocol. !!! Check for the latest version. *This is done for for neutralization of pH.*
Note and warning. The paper specifies pH=7 while the referenced paper specifies pH=7.75
Note: We assume pH is measured at 25 Celsius.
Store at -20.
- 9 Optional, before freezing at -20, aliquot 1 µl measure the result using a NanoDrop or equivalent device. That should normally contain about 100 ng of DNA according to the reference paper.

Protocol references

1 Girish et al: A rapid method for authentication of Buffalo (*Bubalus bubalis*) meat by Alkaline Lysis method of DNA extraction <https://www.ncbi.nlm.nih.gov/pmc/articles/PMC3550954/>

2 Girish et al: Rapid detection of pork using alkaline lysis- Loop Mediated Isothermal Amplification (AL-LAMP) technique. Published, journal pre-proof can be downloaded from Academia https://www.academia.edu/114013841/Rapid_detection_of_pork_using_alkaline_lysiss_Loop_Mediated_Isothermal_Amplification_AL_LAMP_technique