


Oct 24, 2019

Version 1

METHOD FOR THE METABOLIC PROFILE OF PLANT TISSUES BY GAS CHROMATOGRAPHY COUPLED TO MASS SPECTROMETRY (GC/MS) V.1

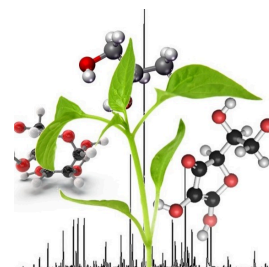

DOI

<https://dx.doi.org/10.17504/protocols.io.wzvff66>

Nívea . Vieira¹, Camilo E. Vital¹, Claudia S.L. Pontes¹, Pedro M. Vidigal¹, Jenny D. Gómez², Edvaldo Barros¹, Humberto Ramos¹

¹Center of Analysis of Biomolecules (NuBioMol), Universidade Federal de Viçosa - UFV, Viçosa-MG, Brazil;

²Department of Biochemistry and Molecular Biology, Universidade Federal de Viçosa - UFV, BIOAGRO/INCT-IPP, Viçosa-MG, Brazil



Humberto J O Ramos

Universidade Federal de Viçosa

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.wzvff66>

Protocol Citation: Nívea . Vieira, Camilo E. Vital, Claudia S.L. Pontes, Pedro M. Vidigal, Jenny D. Gómez, Edvaldo Barros, Humberto Ramos 2019. METHOD FOR THE METABOLIC PROFILE OF PLANT TISSUES BY GAS CHROMATOGRAPHY COUPLED TO MASS SPECTROMETRY (GC/MS). **protocols.io** <https://dx.doi.org/10.17504/protocols.io.wzvff66>

Manuscript citation:

Coutinho, FS; Santos, DS; Lima, LL; Vital, CE; Santos, LA; Pimenta, MR; Silva, JC; Soares-Ramos, JRL; Mehta, A; Fontes, EPB; Ramos, HJO. Mechanism of the Drought Tolerance of a Transgenic Soybean Overexpressing the Molecular Chaperone BiP. Physiology and Molecular Biology of Plants 2019. DOI : 10.1007/s12298-019-00643-x

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 11, 2019

Last Modified: October 24, 2019

Protocol Integer ID: 19221

Keywords: generation of the metabolite profile, metabolite profile, global analysis of the metabolome, metabolome, plant tissues by gas chromatography, metabolite, metabolites for each run, method for the metabolic profile, metabolic profile, mass spectrometry, gas chromatography, plant tissue, plant

Abstract

Plants produce a variety of compounds with diverse structure and abundance that play an important role in the development, growth and response to the environment. Metabolite profiles can be obtained from hundreds of samples by CG/MS (gas chromatography coupled to mass spectrometry). However, the large number of identify metabolites for each run and of treatments and replicates require automatic processing for both spectra (metabolite identification and quantification) and global analysis of the metabolome. Thus, this protocol describes in detail, step-by-step, the generation of the metabolite profiles from GC/MS data, as well as their Exploratory Statistical and Functional Enrichment analyzes.

Materials

REAGENTS

- Methanol (LiChrosolv, Merck, Order n° 1.06001)
- Chloroform (LiChrosolv, Merck , Order n° 1.0244)
- Ribitol (Sigma)
- High pure water (18.2M Ω cm⁻¹) provided by a Milli-Q system (Burlington, Massachusetts, USA)
- Methoxyamine hydrochloride (Sigma Part Number 226904)
- Pyridine (Sigma Part Number 270970)
- N-Methyl-N-(trimethylsilyl)trifluoroacetamide (MSTFA) (Sigma Part number 69479-10 X 1 mL)
- FAME ?????(retention time standard)
- Liquid Nitrogen.

EQUIPMENTS AND SUPPLIES

- Gaseous Chromatography System coupled to mass spectrometry (GC-MS) with the following specifications:
Agilent 7890A GC System coupled to Mass Spectrometric TruTOF® HT TOFMS, Leco
- Capillary column 30 m (MDN-35)
- Thermomixer (Benchtop laboratory incubator)
- Benchtop centrifuge.
- Sppedvac (vacuum centrifugation drier)
- Ultra-freezer.
- Benchtop balance
- Mortar and pestle.
- Vials, caps and septa (PTFE).
- Microtubes e tips (Axygem ou Sigma)

METABOLITE EXTRACTION

- 1
 - 1) Collect samples of plant tissues, immediately freeze in liquid nitrogen and store them in freezer -80°C until use.
 - 2) Macerate the samples in liquid nitrogen using mortar and pestle. Do not allow to thaw. Weigh approximately 100mg of each sample into microtubes (2ml) and annotate the weight (used for normalization).
 - Note: Always use microtubes and tips of good quality.
 - 3) Add 700 µl of methanol and vortex for 10 seconds to stop the enzymatic activity
 - 4) Add 30 µl ribitol from stock solution (0.2 mg/ml in water) as an internal quantitative standard for the polar phase. Vortex for 10 seconds.
 - 5) Use a termomixer to shake 15 minutes at 70°C, 1000 rpm . After 2 minutes of incubation, open the microtubes to relieve internal pressure.
 - 6) Centrifuge the sample for 10 min at 14000 rpm and transfer supernatant to new microtubes.
 - 7) Add 375µl of chloroform and 750µl ultrapure water. Vortex for 15 seconds and centrifuge at 4000rpm for 15 minutes.
 - 8) Collect the upper phase (polar phase) to new microtubes (2ml).
 - Note: It is recommended to choose 2 contrasting samples for initial GC/MS tests. The goal is select the best sample volume for derivatization and GC injection. Test 3 aliquots for each sample (50, 100, 150 µl), evaluate the results in the GC/MS and define the best aliquote to derivatize the remaining samples.
 - 9) Dry the aliquots in vacuum centrifugation drier (speedvac) overnight without heating.

DERIVATIZATION

- 2
 - 1) Dissolve each dried aliquote with 40 µl methoxyamine hydrochloride (20 mg/ml in pyridin) and vortex for 30 seconds. Derivatize also one empty microtube as a 'blank'.
 - 2) Shake in thermomixer for 2 hours at 37°C, 1000 rpm and then centrifuge fastly ("spin") to remove the drops of cover.
 - 3) Add 70µl MSTFA containing retention time standard (Fatty Acid Methyl Esters, FAME) (20µl/ml) and shake in thermomixer for 30 minutes at 37°C.
 - 4) Centrifuge fastly ("spin") to remove the drops of cover and transfer the sample into vials.

METABOLIC PROFILE ANALYSIS BY GC/MS

- 3
 - 1) The method were optimized for a GC/MS system equipped with a capillary column of 30m ((see specifications above).
 - 2) Inject the samples (1µl) in splitless mode at 230 °C using a continuous flow of helium (2 mL/min).

- 3) Maintain the oven temperature initially at 80°C and increase 15°C per minute until it reaches 330°C. Keep this last temperature for 5 minutes.
- 4) Get the mass spectrum by full-scan method ranging from 33 to 600 m/z and use n-alkanes series (retention time standard) to calculate retention index.

DATA PROCESSING AND METABOLITE IDENTIFICATION

- 4 **Note:** For GC/MS TruTOF (Leco), the raw spectra are processed and converted to the CDF format (NetCDF) using the ChromaTof package and analyzed by TargetSearch algorithms (Cuadros-Inostroza et al. 2009). We use a script (Supplementary Material) for identification and quantification of metabolites designed to run on R package (Cuadros-Inostroza et al. 2009). The processing parameters and alignment utilized were optimized for our GC/TOF platform. The parameters used in the TargetSearch were: mass range of 85-500 Da; threshold 50; TopMasses 10; r thresh of 0.05; quality index threshold 600. The compound identification were performed by database searches based on the combination of the mass spectrum and chromatographic retention indices (Cuadros-Inostroza et al. 2009). In this search were used a metabolites fragmentation library by electron impact (EI) and retention index (RI) GMDB_FAMELib_TS_20110228_IS.txt, generated by GMD Mass Spectrum Reference Library (<http://gmd.mpimp-golm.mpg.de/download/>). This library was generated using the same configuration of our GC/MS platform. A filtered identification table containing the identified compounds and their intensities were normalized by fresh mass of the leaves and internal standard ribitol and was used as input data for processing and statistical analysis by MetaboAnalyst platform (<http://www.metaboanalyst.ca/>).

- 1) Install the **R package** for the your operational system (**32 bits ou 64 bits**):

<https://cran.r-project.org/bin/windows/base/>

- 2) Install **OpenOffice**

<https://www.openoffice.org/download/>

- 3) Paste **all the spectra** exported as **.cdf files** from ChromaTof (Leco TruTof users) in your work directory. **C:/myworkdirectory/**

a) Download the files my.work.directory.rar from www.nubiomol.ufv/???

b) Unpack in the "C:/myworkdirectory/". You will see the **R script ("MainScrip.r")**, the library downloaded from GMD Mass Spectrum Reference Library ("**GMDB_FAMELib_TS_20110228_IS.txt**") and the txt files containing information for corrections of the FAMES retention times ("**FAME_limist.txt**" and "**rimLimits.txt**"). The file "**FAME_limist.txt**" and "**rimLimits.txt**" **must to be edited by the user** in accordance with the lower and upper limits of the **TICs for each alkanes**. This information is obtained using ChromaTof package. This file must to be edited for each new analysis cycle.

- Exemple of the Information from **FAME_limist.txt** file:

LowerLimit UpperLimit Standard Mass

267 278 262320 87

328 330 323120 87

382 391 381020 87

486 495 487220 87

575 586 582620 87

662 666 668720 87

739 745 747420 87

810 817 819620 87

876 882 886620 87

937 945 948820 87

994 1000 1006900 87

1048 1054 1061700 87

4) The file "*MainScript.r*" must also be edited:

Open the file in **RStudio** [File > Open file...] (or type Ctrl+O). Edit the following lines of the script:

[Windows]

#Set working directory

setwd("C:\\my.work.directory/")

[Linux]

Set working directory

setwd("~/my.working.directory")

Save the modifications! [File > Save] (or type Ctrl+S)

5) Install the TargetSearch package: Select the lines 9 to 11, then click on "Run" or type "Ctrl+Enter". Wait the installation process to finish

6) Load the TargetSearch and **execute** the script: Select the lines 14 to 102, then click on "Run" or type "Ctrl+Enter"

7) Export data to MetaboAnalyst: Select the lines 107 and 108 of the script to install and load the package **dplyr**.

Edit the lines 110 and 111 of the script, including the correct names of "**profile**" and "**quantmatrix**" files which were created as **output of TargetSearch**:

```
profile_info<-read.delim("TargetSearch-XXXX-XX-XX.profile.info.txt", sep="\t",  
header=TRUE)
```

```
quant_matrix<-read.delim("TargetSearch-XXXX-XX-XX.profile.quantmatrix.txt", sep="\t",  
header=TRUE)
```


8) Select the lines 110 to 121, then click on **"Run"** or type "Ctrl+Enter"

9) Open the file **"Metaboanalyst_data.txt"** in a Text Editor or Spreadsheet Editor such as OpenOffice. Install from <https://www.openoffice.org/download/index.html>

Note: This file contain the **intensity** for each compound identify from spectra GC/MS with score higher than **600** and maintain only the compounds with **high confidence identification**.

Use the parameters below:

Click in **ok!!!**

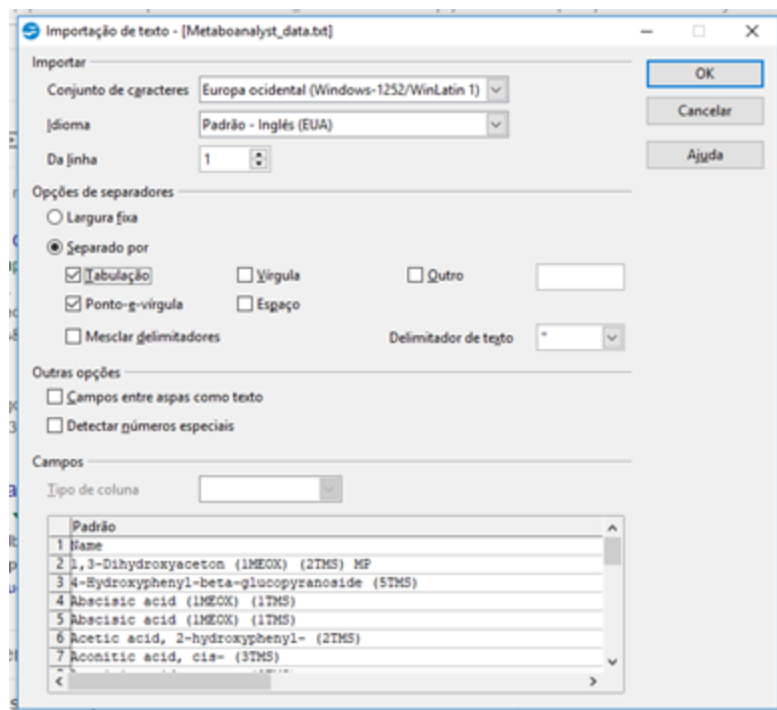

Reorganize the columns to **identify the treatments and replicates**. Insert lines as "name". "sample" and "label" as in the figure below:

Before edition:

Name	Azo_14_R3	Azo_25_R2	Azo_36_R1	CTR_14_R1	CTR_25_R2	CTR_36_R3	Herb_14_R2	Herb_25_R1	Herb_36_R3
1,3-Dihydroxyacetone (1MEOX) (2TMS) MP	NA	70	NA	68	51	58	64	NA	66
4-Hydroxyphenyl-beta-glucopyranoside (3TMS)	1212	782	1561	488	643	662	520	515	1253
Abscisic acid (1MEOX) (1TMS)	NA	NA	61	58	54	NA	54	NA	67
Abscisic acid (1MEOX) (1TMS)	161	77	144	NA	71	87	1067	NA	404
Acetic acid, 2-hydroxyphenyl- (2TMS)	NA	55	73	60	63	78	54	69	NA
Aconitic acid, cis- (3TMS)	303464	309599	322020	255659	311819	319158	301315	243369	296809
Aconitic acid, trans- (3TMS)	19478	19376	20837	NA	11969	NA	NA	NA	18248
Acrylic acid, 4-imidazole- (2TMS)	65	NA	NA	60	65	70	60	NA	NA

After edition:

name	Azo_36_R1	Azo_25_R2	Azo_14_R3	CTR_14_R1	CTR_25_R2	CTR_36_R3	Herb_25_R1	Herb_14_R2	Herb_36_R3
1	1	2	3	1	2	3	1	2	3
4	868146	794995	799669	695700	840948	699469	700343	862915	837617
5	1534	1972	1942	878	1903	841	895	1723	1860
6	322020	309599	303464	255659	311819	319158	243369	301315	296809
7	354466	323384	319754	116400	245922	145650	220858	256255	322623
8	289571	196266	189848	91836	150543	96678	146917	190988	226461
9	55333	46911	62057	29115	37929	53146	36689	46878	62731
10	113490	116243	145055	66731	84669	113444	73906	98891	120072
11	33824	22165	28455	19039	23802	23035	17643	31121	29710
12	30900	28396	27489	22189	25931	25868	11651	27155	24806
13	145	124	142	123	85	128	85	144	138

Save the TXT edited file, such as **"Metaboanalyst_data.txt"**.

Save as **text CSV** "Metaboanalyst_data.csv">>> **select filter configuration**

Score_cor_masses	quantMass	isSelectiveMass	Azo_14_R3	Azo_25_R2	Azo_36_R1
960	87	VERDADEIRO	799669	794995	868146
949	130	VERDADEIRO	1942	1972	1534
949	229	VERDADEIRO	303464	309599	322020
948	217	VERDADEIRO	319754	323384	354466
945	217	VERDADEIRO	189848	196266	289571
945	117	VERDADEIRO	62057	46911	55333
113490	116243	145055	66731	84669	113444
33824	22165	28455	19039	23802	23035
30900	28396	27489	22189	25931	25868
145	124	142	123	85	128
85	128	85	144	138	

Note: This file contain the **intensity** for each compound identify from spectra GC/MS , that could be used for the **statistical** and **pathways** analyzes, compatible by the

MetaboAnalyst platform. It provide a user-friendly, web-based analytical pipeline for high-throughput metabolomics studies. In particular, MetaboAnalyst aims to offer a variety of commonly used procedures for metabolomic data processing, normalization, multivariate statistical analysis, as well as data annotation.

METABOLITE DATA ANALYSIS

5

1) Go to <https://www.metaboanalyst.ca/faces/ModuleView.xhtml>

Click in **Statistical Analysis** >>> **"Select "Peak Intensity Table" and format as "sample in columns (unpaired)".**

Select the file exported from OpenOffice such as **"TargetSearch-2016-06-06.profile.quantmatrixYourData.txt"**

Click **in submit** You will see **error message** such as:

ErrorFailed to read in the CSV file.

Possible causes of error (last one being the most relevant):

Duplicate feature names are not allowed!

Abscisic acid (1MEOX) (1TMS)

The duplicate name **must be renamed**. Go back to the spreadsheet **"TargetSearch-2016-06-06.profile.quantmatrixYourData.txt"** and **edit** using the OpenOffice:

Use CTRL+F

Find Abscisic acid (1MEOX) (1TMS)

Replace Abscisic acid₂ (1MEOX) (1TMS)

Repeat for **all duplicate names** indicated in the error message!!!

Save again as before.

Repeat the process in the **MetaboAnalyst** and **verify** if it was **fixed**.

2) The next module is **"Data Normalization"**.

Note: The sample normalization allows general-purpose adjustment for differences among your sample; data transformation and scaling are two different approaches to make individual features more comparable. You can use one or combine them to achieve better results. For details, see documentation on the MetaboAnalyst site.

Here is also possible normalize the data using the fresh mass of the tissues used for metabolite extraction (Sample-specific normalization). However, for better results proceed it in the spreadsheet "TargetSearch-2016-06-06.profile.quantmatrixYourData.txt". Use the table containing the fresh tissue masses to generate a correction factor for each sample. Localize the sample showing highest fresh weight and call 1.0 (100%), then generated the relative factor for the fresh weight of the others samples that will be relative and higher than 1.0. Thus, differences in the weight between samples are corrected without modified the magnitude of the peak intensity from the GC/MS data. Now, the "corrected" data by fresh weight can be normalized using different statistical methods.

During normalization process the "Data transformation" and "data scaling" have been recommended when some requirement of the parametric statistic, such as normality of the error distribution, homogeneity of variance and etc., are not met. The MetaboAnalyst pipeline enable to **choice the better normalization options**, indicating the result after and before the procedure. Thus, it is necessary that you data be in normal distribution.


The normalization using an internal control is also possible. For this, use "**Normalization by reference feature**".

Find and select as normalizer, in the **tab of the compound list**, the internal control **Ribitol** that was added in the extraction procedure. Data normalization by media or median also show satisfactory results and could be used when normalization by Ribitol is not efficient. See the results before and after normalization to verify if the data is under normal distribution and if the variability were reduced.


After the selecting of normalization method, click in "**Normalize**" and in "**view results**". If the normalization was efficient, click in "**proceed**".

3) Analysis by MetaboAnalyst


Now, you can choice the analysis according with your experimental design and hypothesis. It convenient start with "**Cluster Analysis**" such as PCA enable view the general behaviors of the treatments and replicates, identifying the effect of specific treatments or genotype relative to the metabolite abundances.

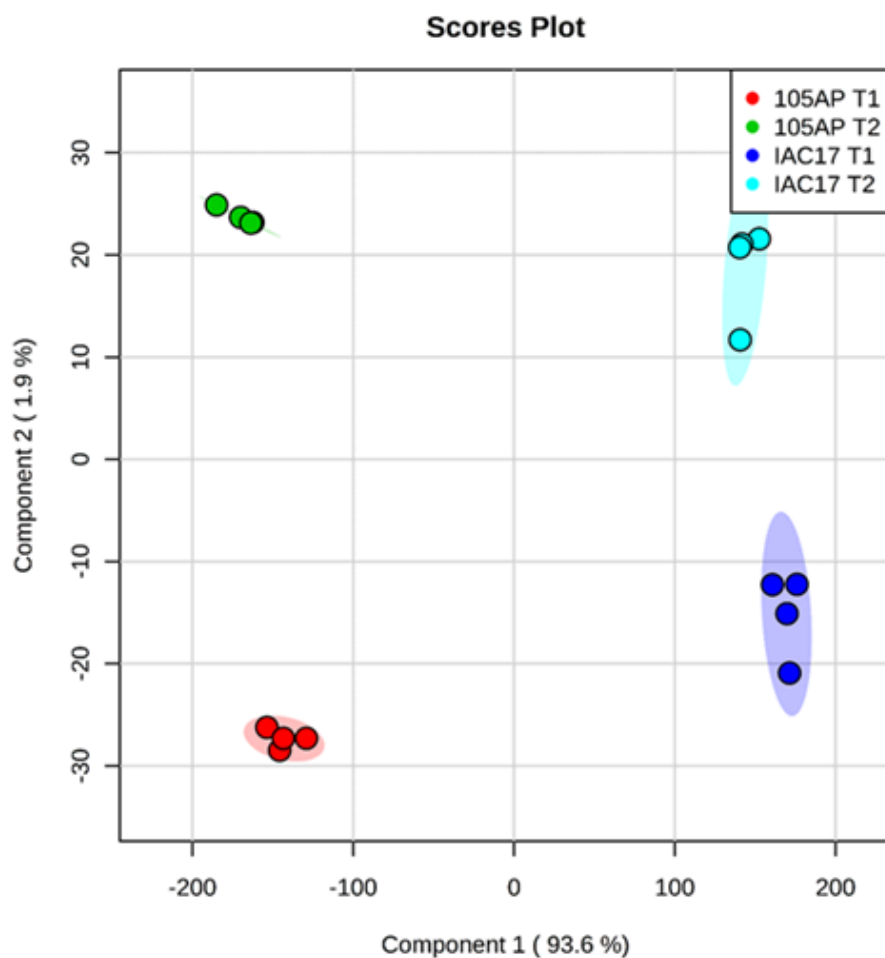

“Cluster Analysis” by PSDLA enable also identify which **metabolites** showed **higher variations** for each group (the output indicate the relative concentrations of the corresponding metabolite in each group under study).

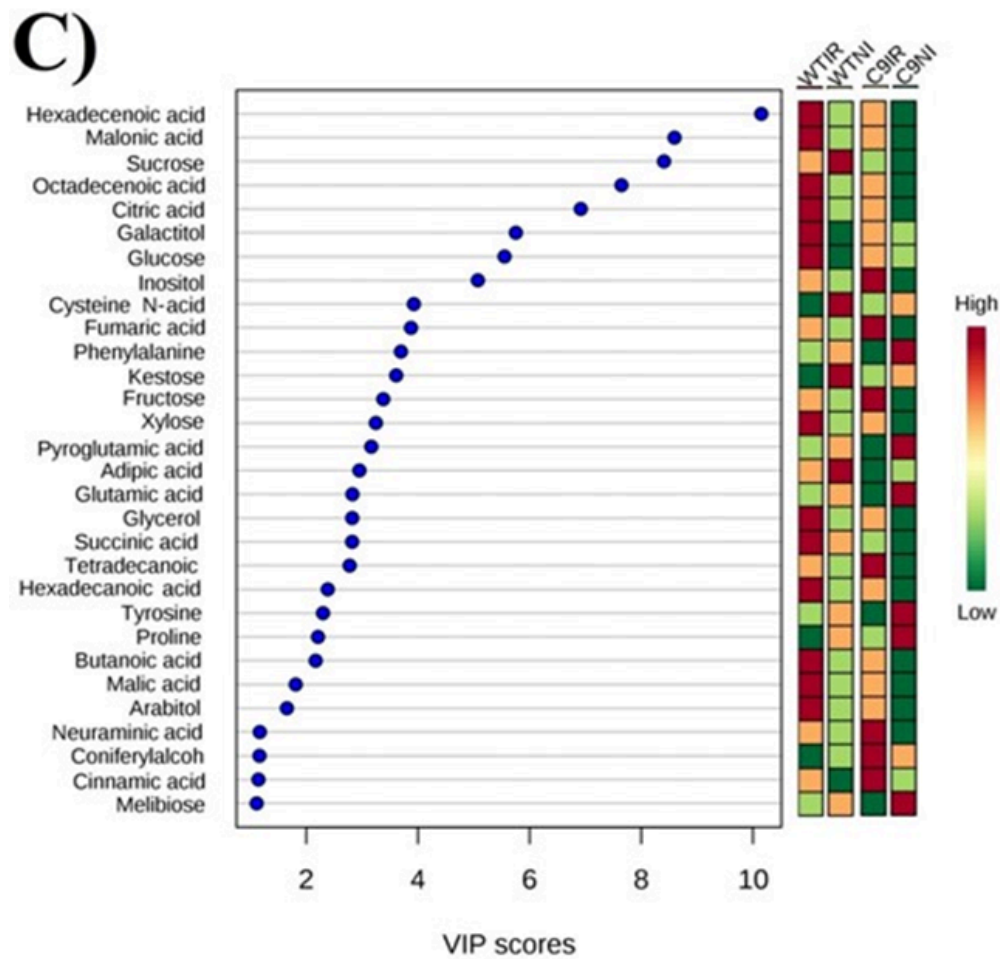


Analysis of variance (**ANOVA**) is very useful to detect which compounds showed significantly variation related to treatments.

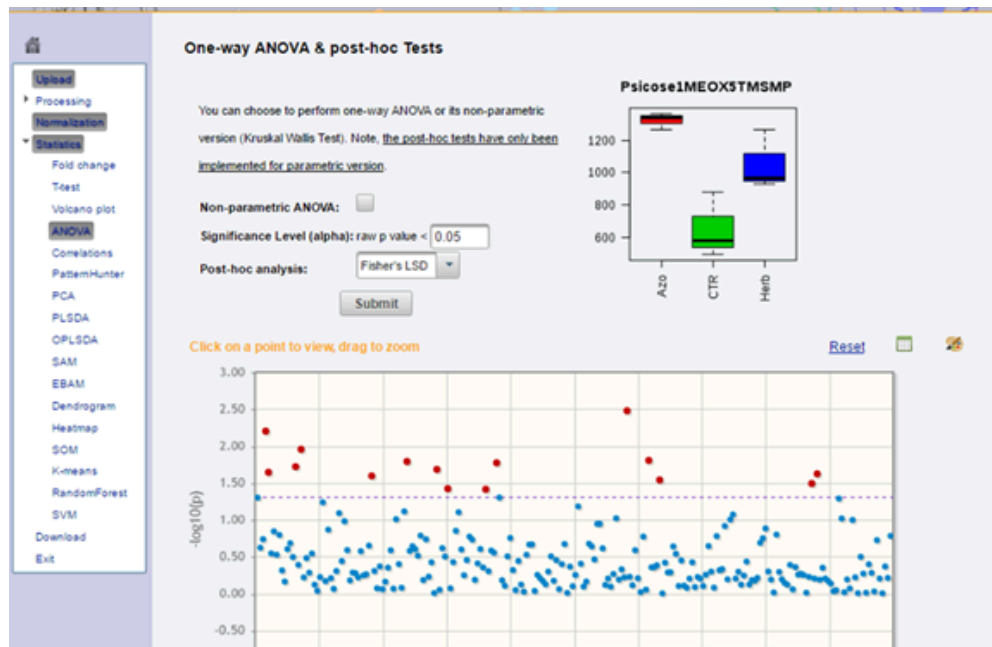

Note: MetaboAnalyst containing a total of 12 modules organized into four general categories: **Exploratory statistical analysis, Functional enrichment analysis, Data integration & systems biology and Data processing & utility functions.**

Pathway Analysis generate the **metabolome views** from a list of compounds for each treatment (with or without intensity). The list is submitted to pipeline, however is necessary altered the compound synonymous in your compound list extracted from the *"TargetSearch-2016-06-06.profile.quantmatrixYourData.txt"*, in order to use the names compatible with the pathway library. The pipeline will suggest synonymous, for example, from KEGG and PubChen. This analyzes shows all matched pathways according to the p values from the pathway enrichment analysis and pathway impact values from the pathway topology analysis.

Overview of Pathway Analysis

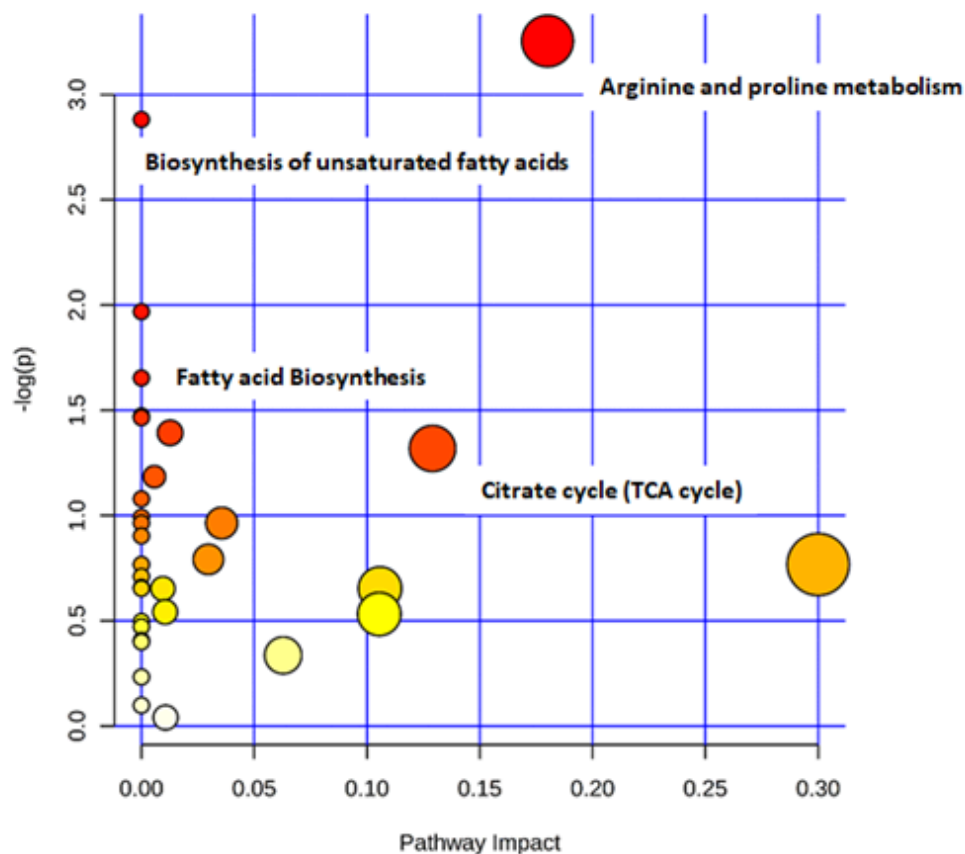

Finally. The result tables and the figures can be **downloaded** with high **resolution**.

ACKNOWLEDGMENTS

- The authors would like to thank the Núcleo de Análises de Biomoléculas (NuBioMol) of the Universidade Federal de Viçosa for providing the facilities for the data analysis. The authors also acknowledge the financial support provided by the following Brazilian agencies: Fundação de Amparo à Pesquisa do Estado de Minas Gerais (Fapemig), Coordenação de Aperfeiçoamento de Pessoal de Nível Superior (CAPES – Finance code 001), Conselho Nacional de Desenvolvimento Científico e Tecnológico (CNPq), Financiadora de Estudos e Projetos (Finep), and Sistema Nacional de Laboratórios em Nanotecnologias (SisNANO)/Ministério da Ciência, Tecnologia e Informação (MCTI).

REFERENCES

- 7 Coutinho, FS; Santos, DS; Lima, LL; Vital, CE; Santos, LA; Pimenta, MR; Silva, JC; Soares-Ramos, JRL; Mehta, A; Fontes, EPB; Ramos, HJO (2019) Mechanism of the Drought Tolerance of a Transgenic Soybean Overexpressing the Molecular Chaperone BiP. *Physiology and Molecular Biology of Plants*. DOI : 10.1007/s12298-019-00643-x.

Lucas Leal Lima, Bruno Pereira Balbi, Rosilene Oliveira Mesquita, José Cleydson Ferreira da Silva, Flaviane Silva Coutinho, Flavia Maria Silva Carmo, Camilo Elber Vital, Angela Mehta, Marcelo Elher Loureiro, Elizabeth Pacheco Batista Fontes, Everaldo Gonçalves Barros, Humberto Josué Oliveira Ramos (2019) Proteomic and Metabolomic Analysis of a Drought Tolerant Soybean Cultivar from Brazilian Savanna. *Crop Breeding, Genetics and Genomics* 2019;1(2):e190022; DOI:10.20900/cbagg20190022

Lisec J, Schauer N, Kopka J, Willmitzer L, Fernie A R (2006). Gas chromatography mass spectrometry-based metabolite profiling in plants. *Nature Protocols*, 1: 387–396.

Cuadros-Inostroza Á, Caldana C, Redestig H et al (2009) TargetSearch-a Bioconductor package for the efficient preprocessing of GC-MS metabolite profiling data. *BMC bioinformatics* 10(1): 428. <https://doi.org/10.1186/1471-2105-10-428>

Vital CE, Giordano A, de Almeida Soares E, Rhys Williams TC, Mesquita RO, Vidigal PMP, de Santana Lopes A, Pacheco TG, Rogalski M, de Oliveira Ramos HJ, Loureiro ME (2017). An integrative overview of the molecular and physiological responses of sugarcane under drought conditions. *Plant Mol Biol*. 6:577-594. DOI: 10.1007/s11103-017-0611-y.