




Sep 27, 2023

🌐 MicroCT Imaging of the Fascicular Structure in the Porcine Right and Left Cervical Vagus Nerve

DOI

<https://dx.doi.org/10.17504/protocols.io.ewov1qd2ygr2/v1>

MicroCT Imaging of the Fascicular Structure in the Porcine Right and Left Cervical Vagus Nerve

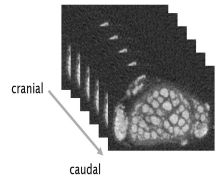

Bettina Kronsteiner¹, Max Haberbush², Francesco Moscato²

¹Center of Anatomy and Cell Biology, Medical University of Vienna;

²Center for Medical Physics and Biomedical Engineering, Medical University of Vienna

SPARC

Tech. support email: info@neuinfo.org



Max Haberbush

Center for Medical Physics and Biomedical Engineering, Medic...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.ewov1qd2ygr2/v1>



Protocol Citation: Bettina Kronsteiner, Max Haberbush, Francesco Moscato 2023. MicroCT Imaging of the Fascicular Structure in the Porcine Right and Left Cervical Vagus Nerve. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.ewov1qd2ygr2/v1>

Manuscript citation:

Kronsteiner, Bettina, et al. "Mapping the functional anatomy and topography of the cardiac autonomic innervation for selective cardiac neuromodulation using MicroCT." *Frontiers in Cell and Developmental Biology* 10 (2022): 968870.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: September 14, 2023

Last Modified: September 27, 2023

Protocol Integer ID: 87769

Keywords: vagus nerve in domestic pig, microct imaging of the fascicular structure, microct imaging, left cervical vagus nerve, fascicular anatomy, vagus nerve, left cervical vagus nerve this protocol, domestic pig, fascicular structure, porcine

Funders Acknowledgements:

European Project H2020

Grant ID: 824071

Austrian Research Promotion Agency R&D Infrastructure Project

Grant ID: 858060

Abstract

This protocol outlines the procedure to dissect, prepare, and perform micro-CT imaging of the fascicular anatomy of the right and left vagus nerve in domestic pigs.

Materials

- Fresh cadavers of pigs (domestic pigs)
- Isotonic sodium chloride solution (NaCl, 0.9%)
- Surgical sutures (5-0 silk suture, Silkam® B-Braun Aesculap, Tutlingen, Germany)
- Polystyrene plate with V-shaped incisions
- PBS (Phosphate-Buffered Saline)
- 4% phosphate-buffered 4% Paraformaldehyde (Roti-Histofix, Roth, Karlsruhe, Germany)
- 1% Lugol's solution
- GIMP open-source software (GIMP 2.10.24, revision 2)
- SCANCO μ CT 50 (SCANCO Medical AG, Brütistellen, Switzerland) specimen μ CT scanner

Surgical Dissection and Macroscopic Examination of Nerve Samples

- 1 Acquire fresh cadavers of male and/or female domestic pigs.
- 2 Place cadavers in the dorsal recumbence position.
- 3 Make a ventral skin incision (15–20 cm) on both sides next to the trachea starting at the level of the mandibles extending to the sternal notch.
- 4 Dissect tissue and open the carotid sheath to expose the common carotid artery, internal jugular vein, the vagus nerve, and the sympathetic trunk.
- 5 Begin nerve dissection at the level of the nodose ganglion (NG) and extend caudally to the superior cardiac branch, which originates approximately 2–5 cm cranial to the subclavian artery on both sides.
- 6 Track branches emerging from and terminating in the vagus nerve (VN) to their origin structures.
- 7 Identify both recurrent laryngeal nerves running close to the trachea looping back around the subclavian arteries.
- 8 Open the chest by median sternotomy followed by gentle exposure of the heart.
- 9 Continue surgical preparation of the VN towards the heart. Dissect the superior cardiac branch from the cardiac branching points at the VN towards the insertion points at the heart, ensuring not to cut any of the cross-connections between the VN and the sympathetic trunk (ST).
- 10 During dissection, moisten the nerves with isotonic sodium chloride solution to prevent drying.
- 11 Measure dimensions of the VN using a ruler and capture a photo for documentation.

Vagal Nerve Preparation and Harvesting

- 12 Fix surgical sutures onto the epineurium of the VN below the nodose ganglion to maintain the anatomical in-situ orientation. Nerves should be harvested 1-2 hours post-mortem.
- 13 After complete dissection, place and fix nerves onto a polystyrene plate, ensuring the anatomical in-situ orientation is maintained.
- 14 Wash nerve specimens in phosphate-buffered saline (PBS)
- 15 Fixate in 4% phosphate-buffered 4% Paraformaldehyde (PFA) at 4°C for 16–18 hours.
- 16 After fixation, wash the nerve samples again and store in freshly made PBS at 4°C until staining for micro-computed tomography imaging.

Micro-computed Tomography

- 17 Stain nerves in 14 mL of 1% Lugol's solution for 24 hours.
- 18 Place nerve specimens in 15 ml conical centrifuge tubes with residual staining solution and the polystyrene plate used for fixation.
- 19 Perform low-resolution scanning using the SCANCO μ CT 50 scanner with 70 kVp (200 μ A, 0.5 mm Al filter, 550 projections, 100 ms integration time) at an isotropic resolution of 20 μ m.
- 20 Divide the VNs into two nerve specimens for scanning, ensuring proper anatomical orientation for alignment.
- 21 Perform high-resolution scans at the cardiac branching point with 70 kVp (85 μ A, 0.5 mm Al filter, 850 projections, 350 ms integration time) at an isotropic resolution of 8.9 μ m over 1 cm of nerve length.
- 22 Re-stain nerves 24 hours prior to high-resolution scans to maintain full saturation with Lugol's solution.
- 23 Analyze high-resolution data at the cardiac branching point.



Protocol references

Kronsteiner, Bettina, et al. "Mapping the functional anatomy and topography of the cardiac autonomic innervation for selective cardiac neuromodulation using MicroCT." *Frontiers in Cell and Developmental Biology* 10 (2022): 968870. doi: <https://doi.org/10.3389/fcell.2022.968870>