



Sep 23, 2023

Microscopy-based bead assay

 [Nature Structural & Molecular Biology](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.14egn38pzl5d/v1>

Elias Adriaenssens¹

¹Sascha Martens lab, University of Vienna, Max Perutz Labs - Vienna



Elias Adriaenssens

Sascha Martens lab, University of Vienna, Max Perutz Labs - ...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn38pzl5d/v1>

External link: <https://doi.org/10.1038/s41594-024-01338-y>

Protocol Citation: Elias Adriaenssens 2023. Microscopy-based bead assay. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.14egn38pzl5d/v1>

**Manuscript citation:**

Elias Adriaenssens, Thanh Ngoc Nguyen, Justyna Sawa-Makarska, Grace Khuu, Martina Schuschnig, Stephen Shoebridge, Marvin Skulsuppaisarn, Emily Maria Watts, Kitti Dora Csalyi, Benjamin Scott Padman, Michael Lazarou, Sascha Martens (2024) Control of mitophagy initiation and progression by the TBK1 adaptors NAP1 and SINTBAD. Nature Structural & Molecular Biology doi: [10.1038/s41594-024-01338-y](https://doi.org/10.1038/s41594-024-01338-y)

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 07, 2023

Last Modified: June 01, 2024

Protocol Integer ID: 84651

Keywords: Microscopy-based bead assay, ASAPCRN, based bead assay, bead assay, based bead, microscopy, assay

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000350

Marie Skłodowska-Curie MSCA Postdoctoral fellowship

Grant ID: 101062916

Abstract

This protocol describes the microscopy-based bead assay.

Materials

Materials

- Glutathione Sepharose 4B beads (GE Healthcare)
- dH₂O
- MgCl₂
- FIP200-GFP
- SINTBAD-mCherry
- TBK1
- SINTBAD-GFP
- mCherry-OPTN
- Zeiss LSM 700 confocal microscope

Bead assay buffer

	A	B
	Tris-HCl pH 7.4	25 mM
	NaCl	150 mM
	DTT	1 mM



Microscopy-based bead assay

6h 30m

- 1 We use Glutathione Sepharose 4B beads (GE Healthcare) to bind GST-tagged bait proteins.
- 2 To this end, wash  20 μL of beads twice with dH_2O and equilibrate with bead assay buffer. 
- 3 Then, resuspend beads in  40 μL bead assay buffer, to which bait proteins are added at a final concentration of  5 micromolar (μM) .
- 4 Incubate the Beads with the bait proteins for  01:00:00 at  4 $^{\circ}\text{C}$ at a horizontal tube roller.  1h
- 5 Wash the beads three times to remove unbound GST-tagged bait proteins and resuspend in  30 μL bead assay buffer. 
- 6 Where needed, also add MgCl_2 and ATP to the buffer to allow the phosphorylation of targets by TBK1 or other kinases. 
- 7 Prepare glass-bottom 384-well microplates (Greiner Bio-One) with  20 μL samples containing prey proteins at the concentrations described below and dilute in bead assay buffer, and add  3 μL of beads per well. 
- 8 For the experiments in which full-length FIP200-GFP is recruited to GST-4xUb beads in presence of NDP52 and/or SINTBAD-mCherry, use NDP52 at a final concentration of  50 nanomolar (nM) , FIP200-GFP, SINTBAD-mCherry, and TBK1 are used at a final concentration of  100 nanomolar (nM) .
- 9 For recruitment of SINTBAD-GFP to GST-LC3/GABARAP beads, use a final concentration of  200 nanomolar (nM) SINTBAD-GFP.
- 10 For the TBK1-binding competition experiment between OPTN and NAP1, use mCherry-OPTN and GFP-TBK1 at a final concentration of  500 nanomolar (nM) , and use NAP1 from  100 nanomolar (nM) to  10 micromolar (μM) .

11 Incubate the beads with the prey proteins for  00:30:00 prior to imaging, with the exception of experiments where full-length FIP200 is recruited, where proteins are co-incubated for  04:00:00 before imaging. For recruitment of autophagy components to GST-LC3/GABARAP beads, proteins are co-incubated for  01:00:00 before imaging.

5h 30m



12 Image samples with a Zeiss LSM 700 confocal microscope equipped with Plan Apochromat 20X/0.8 WD 0.55 mm objective.

**Note**

Three biological replicates were performed for each experimental condition.

13 For the quantification, we employ an artificial intelligence (AI) script that automatically quantifies signal intensities from microscopy images by drawing line profiles across beads and recording the difference between the minimum and maximum grey values along the lines.

Note

The AI was trained to recognize beads employing cellpose [60].

14 Processing is composed of two parts, with the first operating in batch mode. Multichannel input images are split into individual TIFF images and passed to cellpose (running in a Python environment).

15 The labeled images produced by cellpose are re-assembled into multichannel images.

16 Circular regions of interest (ROIs) are fitted to the segmented particles, and a pre-defined number of line profiles (here set to 20) are drawn automatically, starting at the center of the ROI and extending beyond the border of the circular ROI.

17 This results in line profiles from the center of the bead into the inter-bead space of the well, allowing us to quantify the signal intensities at the rim of the beads.

- 18 To prevent line profiles from protruding into adjacent beads, a combined ROI containing all beads is used.
- 19 Inspect the AI-generated results manually for undetected beads, incorrect line profiles, or false-assigned bead structures.
- 20 For each bead, a mean fluorescence and standard deviation are obtained based on the 20 line profiles per bead.
- 21 Either exclude or subject the beads with standard deviations equal to or greater than half the mean value to manual inspection for correction.
- 22 If needed, correct for inter-experiment variability in absolute values by dividing the mean values for each bead by the average bead intensity of the control condition, to obtain relative values.
- 23 These values are then plotted and subjected to statistical significance calculations.

Protocol references

Schneider CA, Rasband WS, Eliceiri KW: **NIH Image to ImageJ: 25 years of image analysis.** *Nat Methods* 2012, **9**(7):671-675.

Stringer C, Wang T, Michaelos M, Pachitariu M: **Cellpose: a generalist algorithm for cellular segmentation.** *Nat Methods* 2021, **18**(1):100-106.