

Dec 09, 2024

Microsomal stability assay for human and mouse liver microsomes - drug metabolism

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvkdb9l4o/v1>

Lina Bortnichuk¹, Anna Pashchenko¹, Yurii Kheilik¹

¹Enamine Ltd

ASAP Discovery



Nick Lynch

Curlew Research, ASAP Discovery Consortium

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvkdb9l4o/v1>

Protocol Citation: Lina Bortnichuk, Anna Pashchenko, Yurii Kheilik 2024. Microsomal stability assay for human and mouse liver microsomes - drug metabolism. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.5qpvkdb9l4o/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: August 14, 2024

Last Modified: December 09, 2024

Protocol Integer ID: 106088

Keywords: Microsomal protein, Glucose-6-phosphate dehydrogenase, NADPH-cofactor , Intrinsic clearance , HPLC-MS/MS , Microsomes, microsomal stability, oral bioavailability, drug metabolism, drug metabolism microsomal stability, liver microsomal stability, mouse liver microsome, microsomal stability, metabolism by liver enzyme, liver microsomal incubation medium, cytochrome p450, mg of liver, liver enzyme, metabolism, microsomal protein, microsomal incubation, mouse microsome, key parameter in drug discovery, microsomal incubation medium, liver, microsome, assay, drug discovery, tandem mass spectrometer

Disclaimer

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Microsomal stability is a key parameter in drug discovery and development to support understanding of a compound's susceptibility to liver metabolism.

This assay assesses a compound's susceptibility to metabolism by liver enzymes, primarily cytochrome P450 (CYP) enzymes. Using either human or mouse microsomes.

The data analysis provides elimination constant, half-life, and intrinsic clearance. These are determined as results using linear regression analysis.

In this protocol, we measure (human or mouse) liver microsomal stability assay for 12 test articles in 96-well plate format. Microsomal incubations are carried out in 5 aliquots of 30 μ L (one for each time point). The liver microsomal incubation medium comprises phosphate buffer (100 mM, pH 7.4), $MgCl_2$ (3.3 mM), NADPH (3 mM), glucose-6-phosphate (5.3 mM), glucose-6-phosphate dehydrogenase (0.67 units/mL) with 0.415 mg of liver microsomal protein per mL. In the negative control reactions, the NADPH-cofactor system is substituted with phosphate buffer. Test compounds (2 μ M, final acetonitrile concentration 1.6 %) are incubated with microsomes at 37°C, shaking at 100 rpm. Five time points over 40 minutes are analyzed (0, 7, 15, 25, and 40 min). The reactions are stopped by adding 5 volumes of acetonitrile with internal standard to incubation aliquots, followed by protein sedimentation by centrifuging at 5500 rpm for 5 minutes. Supernatants are analyzed using the HPLC-MS/MS system coupled with a tandem mass spectrometer.

The elimination constant (k_{el}), half-life ($t_{1/2}$), and intrinsic clearance (Cl_{int}) are determined in a plot of $\ln(AUC)$ versus time, using linear regression analysis:

$$k_{el} = -slope$$

$$t_{1/2} = \frac{0.693}{k}$$

$$Cl_{int} = \frac{0.693}{t_{1/2}} \times \frac{\mu l_{incubation}}{mg_{microsomes}}$$

Materials

Reagents and preparations:

- 100 mM (1x) Potassium phosphate buffer (PPB), pH 7.4
- 33 mM $MgCl_2 \times 6H_2O$ solution
- Glucose-6-phosphate dehydrogenase (G6PDH) solution, 250 U/mL
- Human or mouse liver microsomes, 20 mg/mL
- Compounds stock solutions

Safety warnings

- ! Always wear appropriate PPE for this protocol
Refer to Material Safety Data Sheets for additional safety and handling information.

Reagents and preparations

- 1 Prepare [M] 20 millimolar (mM) stocks of reference and test compounds in DMSO.
 - Then, dissolve DMSO stock in acetonitrile (AcN) to [M] 125 micromolar (μM) , and mix by vortexing. Place AcN stock solutions as shown in **Fig. 1**.
 - **Diclofenac** and **propranolol** are used as reference compounds in the assay with human microsomes, **imipramine** and **propranolol** - with mice microsomes.

Compounds stock solutions in AcN

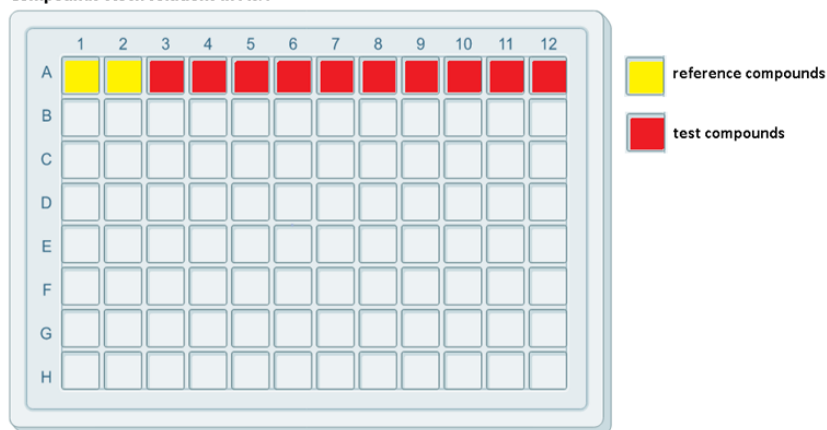


Fig. 1 Plate layout for compounds stock solutions in acetonitrile

2 Quenching solution with internal standards

Dissolve appropriate for LC-MS/MS analysis internal standards in ACN.

3 NADPH tetrasodium salt, D-Glucose-6-phosphate sodium salt

Weigh out 17.88 mg of NADPH tetrasodium salt (MW=833.35 g/mol) and

9.72 mg of D-Glucose-6-phosphate sodium salt (MW=282.12 g/mol) into separate tubes. Place tubes with powders at $-20\text{ }^{\circ}\text{C}$.

Samples preparation

10m

- 4 Set the incubator-shaker at 37°C and 100 rpm .
- 5 Place aliquots of microsomes and G-6-PDH 0°C On ice to thaw.
- 6 Prepare five 96-well plates and label them as «Preincubation», «K+/K-», «0+/0-», «40-», and «Incubation».
 - Place microtubes into plates as shown in **Fig. 2**.

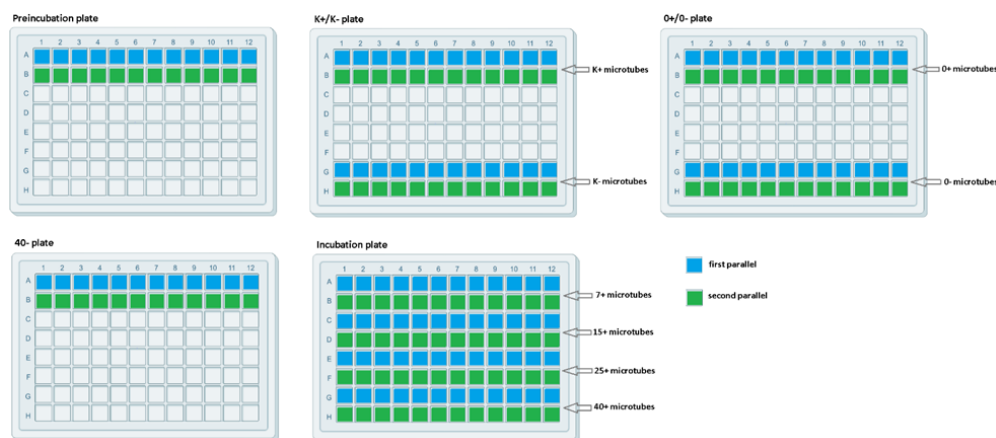


Fig. 2 Preparation of the 96-well plates with 0.75mL microtubes for the assay

- 7 Label tubes as «K-», «K+», and «Microsomes».

Prepare the following solutions:

- «K-». Mix $1021\ \mu\text{L}$ [M] 100 millimolar (mM) PPB and $325\ \mu\text{L}$ [M] 33 millimolar (mM) $\text{MgCl}_2 \times 6\text{H}_2\text{O}$.
- «K+». Mix $1388\ \mu\text{L}$ [M] 100 millimolar (mM) PPB and $650\ \mu\text{L}$ [M] 33 millimolar (mM) $\text{MgCl}_2 \times 6\text{H}_2\text{O}$ solution.
- «Microsomes». Fill the tube with $5785\ \mu\text{L}$ [M] 100 millimolar (mM) PPB.

- 8 Aliquot $52\ \mu\text{L}$ of «K-» solution into the corresponding microtubes of «K+/K-» plate, and then cover the microtubes with mates until starting the experiment.



- 9 Dispense the stop solution into the «0+/0-» plate, and then cover the microtubes with mates until the start of the experiment.
- 10 After thawing, add  216 μL of **human** or **mouse liver microsomes** into the tube labeled «Microsomes», mix manually, and place the tube  On ice . 
- 11 Aliquot  231 μL of «Microsomes» solution into the «Incubation» plate. Then, add  6.5 μL  125 micromolar (μM) AcN-stock solutions of compounds, and mix by pipetting.  10m
- Cover microtubes with mates and place the plate in the incubator-shaker for  00:10:00 to pre-warm the solution before the experiment.
- 12 Within 10 minutes perform the following steps:
- 12.1 Dissolve  17.88 mg of NADPH in  325 μL of  100 millimolar (mM) PPB, mix by pipetting; 
- 12.2 Dissolve  9.72 mg of D-Glucose-6-phosphate in the same way.
- 12.3 Add  325 μL of both NADPH and D-Glucose-6-phosphate solutions into the «K+» tube;
- 12.4 Add  16 μL of G-6-PDH into the «K+» tube, gently mix by pipetting; 
- 12.5 Aliquot  104 μL of «K+» solution into the corresponding microtubes of «K+/K-» plate.
- 12.6 Fill in the «0+/0-» plate with the desired amount of quenching solution.

Incubation

45m

- 13 After 10 minutes of pre-incubation, remove the «Preincubation» plate from the incubator-shaker, and mix samples by pipetting. ✂
- 14 **«K-» samples.** Transfer 🧪 71 µL of pre incubation solution into the first parallel of «K-» microtubes, mix by pipetting, and aliquot solution into the corresponding «40-» and «0-» microtubes. Perform the same operation with another parallel. ✂
- 15 **«K+» samples.** Transfer 🧪 142 µL of preincubation solution into the first parallel of «K+» microtubes, mix by pipetting, aliquot solution into the corresponding «0+» microtubes (start the timer from this moment), and into the microtubes of «Incubation» plate (7-, 15-, 25-, 40-min time points). ✂
 - Perform the same operation with another parallel and record the time shift between the 0-min time points of the first and second incubation parallels.
- 16 Cover «0+/0-» microtubes with caps, vortex plate and place in the refrigerator. Cover «7», «15», «25», «40», and «40-» microtubes with mats, and place in the incubator-shaker.
- 17 At each time point, remove corresponding microtubes from the incubator-shaker and add quenching solution to stop the reaction.
- 18 After the total incubation time, centrifuge the final plates at 🌀 5500 rpm, 00:05:00. 5m
The final plates should look as shown in **Fig. 3**. 🌀

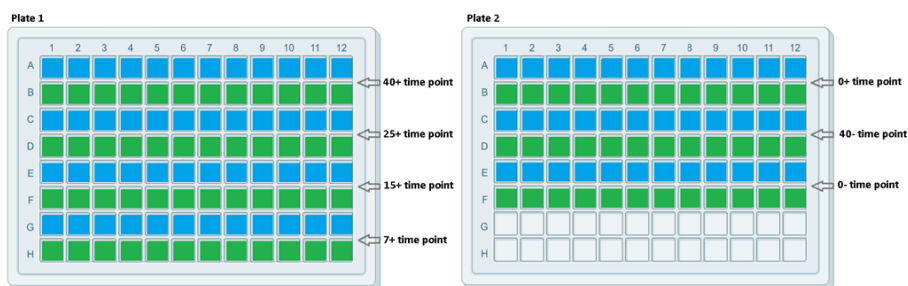


Fig. 3 Sample layout on the final plate

- 19 Analyze the samples using HPLC-MS/MS in the following order (start the analysis from the samples with theoretically lower concentrations of the tested compound): 40m



00:40:00

time point→25 min→15 min→7 min→0 min→40- min →0-.



Data Analysis

- 20 The elimination constant (k_{el}), half-life ($t_{1/2}$), and intrinsic clearance (Cl_{int}) are determined in a plot of $\ln(AUC)$ versus time, using linear regression analysis:

$$k_{el} = -slope$$

$$t_{1/2} = \frac{0.693}{k}$$

$$Cl_{int} = \frac{0.693}{t_{1/2}} \times \frac{\mu l_{incubation}}{mg_{microsomes}}$$

Equations used