

Aug 10, 2023

Miltenyi Biotec: Adipose Tissue Dissociation -- University of Minnesota TMCs

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg39jxqg25/v1>

Miltenyi Biotec¹, Laura J Niedernhofer², David A Bernlohr²

¹Miltenyi Biotec, technicalsupportUS@miltenyi.com; ²University of Minnesota, Minneapolis, MN USA

Cellular Senescence Net...



Mickayla DuFresne-To

University of Minnesota - Twin Cities

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg39jxqg25/v1>

Protocol Citation: Miltenyi Biotec, Laura J Niedernhofer, David A Bernlohr 2023. Miltenyi Biotec: Adipose Tissue Dissociation -- University of Minnesota TMCs. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.kqdg39jxqg25/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 04, 2023

Last Modified: August 10, 2023

Protocol Integer ID: 81368

Keywords: Dissociation , Adipose, Miltenyi, University of Minnesota, adipose tissue dissociation, human adipose tissue, human adipose, adipose, miltenyi biotec, using miltenyi dissociation kit catalog, miltenyi dissociation kit catalog, minnesota tmcs dissociation, university of minnesota tmcs dissociation, dissociation

Funders Acknowledgements:

NIH

Grant ID: 1U54AG07041-01

NIH

Grant ID: 1U54AG079754-01

Abstract

Dissociation of human adipose tissue using Miltenyi dissociation kit Catalog #130-105-808 and Protocol #140-004-708.02



1



Adipose Tissue Dissociation Kit (M...

<https://www.miltenyibiotec.com/US-en/products/adipose-tissue-dissociation-kit-mouse-and-rat.html#130-105-808>