

Sep 23, 2023

Mitochondrial isolation protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg3x4zzg25/v1>

Elias Adriaenssens¹

¹Sascha Martens lab, University of Vienna, Max Perutz Labs - Vienna (AT)



Elias Adriaenssens

Sascha Martens lab, University of Vienna, Max Perutz Labs - ...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg3x4zzg25/v1>

Protocol Citation: Elias Adriaenssens 2023. Mitochondrial isolation protocol. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.kqdg3x4zzg25/v1>

Manuscript citation:

The methodology is based on Wieckowski et al. 2009 Nat Protocols.

Wieckowski, M., Giorgi, C., Lebiezinska, M. et al. Isolation of mitochondria associated membranes and mitochondria from animal tissues and cells. Nat Protoc 4, 1582–1590 (2009).

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 26, 2023

Last Modified: May 31, 2024

Protocol Integer ID: 84020

Keywords: Mitochondrial isolation, HeLa cells, ASAPCRN, crude mitochondrial fractions from hela cell, mitochondrial isolation protocol, mitochondrial isolation protocol this protocol, crude mitochondrial fraction, hela cell

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000350

Marie Skłodowska-Curie MSCA Postdoctoral fellowship

Grant ID: 101062916

Abstract

This protocol describes how to isolate crude mitochondrial fractions from HeLa cells.

Attachments



[752-1920.pdf](#)

50KB

Materials

MATERIALS

Mitochondrial isolation buffer:

	A	B
	HEPES (pH 7.4)	5 mM
	mannitol	250 mM
	EGTA	0.5 mM

make fresh on the day itself!

Oligomycin/Antimycin A cocktail:

	A	B
	Oligomycin	10 μ M
	Antimycin A	4 μ M

In case cells were treated for more than 8 hours, we also added 10 μ M Q-VD-OPh (A1901, ApexBio) to suppress apoptosis.

RIPA buffer:

	A	B
	Tris-HCl pH 8.0	50 mM
	NaCl	150 mM
	sodium deoxycholate	0.50%
	SDS	0.10%
	NP-40	1%

supplemented by cOmplete EDTA-free protease inhibitors (11836170001, Roche) and phosphatase inhibitors (PhosphoSTOP, 4906837001, Roche).

 Oligomycin A **ApexBio Technology Catalog #A5588**

 Antimycin A from Streptomyces sp. **Merck MilliporeSigma (Sigma-Aldrich) Catalog #A8674**



 Q-VD-OPh **ApexBio Technology Catalog #A1901**

 cOmplete™, Mini, EDTA-free (Protease Inhibitor) **Roche Catalog ##11836170001**

 Roche PhosSTOP™ **Merck MilliporeSigma (Sigma-Aldrich) Catalog #4906837001**



Mitochondrial isolation

1h 15m

- 1 Seed HeLa cells in 15 cm dishes and grow until confluence. Treat the cells with DMSO or the mitophagy-inducing cocktail Oligomycin/Antimycin A (O/A) if needed.
- 2 Collect HeLa cells by trypsinization and resuspension in DMEM medium. Centrifuge the cells at  300 x g, 4°C, 00:10:00 .

10m


- 2.1 Resuspend the cell pellet in  1 mL of ice-cold PBS (1x) and spin again for  300 x g, 4°C, 00:05:00 .

5m
- 3 Remove the PBS and resuspend the cell pellet in  1 mL mitochondrial isolation buffer, which was cooled to  4 °C .

Note

Note that the mitochondrial isolation buffer is prepared fresh on the day itself.

- 4 To lyse cells without damaging the mitochondria, pipet the  1 mL cell suspension up and down (15-20x) with a 26.5G needle.

Note

This should lyse the plasma membrane but leave organelles intact.

- 5 Spin the suspension down at  600 x g, 4°C, 00:10:00 .

10m



- 5.1 Keep the supernatant and centrifugation  600 x g, 4°C, 00:10:00 .

10m

Note

The mitochondria are located in the supernatant at this speed, the pellet contains intact cells, cell nuclei, and other large cell debris.



6 After two spins at $600 \times g$, subject the supernatant to $7000 \times g$, 4°C , 00:10:00 .

10m



6.1 The pellet contains the mitochondria, so remove the supernatant and resuspend the pellet in 1 mL mitochondrial isolation buffer.

6.2 Then, centrifugation $7000 \times g$, 4°C , 00:10:00 .

10m

Note

The supernatant from the first spin at $7000 \times g$ can be stored as a cytosolic fraction.

7 After two spins at $7000 \times g$, subject the resuspended pellet to $10000 \times g$, 4°C , 00:10:00 .

10m



7.1 The pellet contains the mitochondria, so remove the supernatant and resuspend the pellet in 1 mL mitochondrial isolation buffer.

7.2 Then, centrifugation $10000 \times g$, 4°C , 00:10:00 .

10m

8 After two spins at $10000 \times g$, the pellet can be resuspended in RIPA lysis buffer or used for further procedures such as mitochondrial import assays, proteinase K protection assays, etc.