

Oct 16, 2025

Mitochondrial ROS Analysis

 [Science Advances](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vz9x98gx1/v1>

Szu-Chi Liao¹, Ken Nakamura¹

¹Gladstone Institutes



Szu-Chi Liao

Gladstone Institutes

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vz9x98gx1/v1>

External link: <https://doi.org/10.1126/sciadv.adu0726>

Protocol Citation: Szu-Chi Liao, Ken Nakamura 2025. Mitochondrial ROS Analysis. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.rm7vz9x98gx1/v1>

Manuscript citation:

Liao S, Kano K, Phanse S, Nguyen M, Margolis E, Fu Y, Meng JX, Moutaoufik MT, Chatterton Z, Saccon T, Broderick K, Aoki H, Simms J, Suteja FX, Sei Y, Huang EJ, McAvoy K, Manfredi G, Halliday G, Babu M, Nakamura K CHCHD2 mutant mice link mitochondrial deficits to PD pathophysiology. Science Advances 11(46). doi: [10.1126/sciadv.adu0726](https://doi.org/10.1126/sciadv.adu0726)



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 19, 2025

Last Modified: October 16, 2025

Protocol Integer ID: 225026

Keywords: ASAPCRN, mitochondrial ros analysis this protocol, mitochondrial ros analysis, levels in isolated mitochondria, mitochondria, isolated mitochondria, ro, h2o2

Abstract

This protocol is used to measure ROS (H₂O₂) levels in isolated mitochondria.

Materials

- Amplex Red Hydrogen Peroxide/Peroxidase Assay Kit (Invitrogen A22188)
- Mitochondrial respiration buffer
- Respiratory substrate solution

Reagent Preparation

- 1 Prepare mitochondrial respiration buffer (100 mL; 120 mM KCl, 5 mM KH_2PO_4 , 10 mM HEPES, 2 mM MgCl_2 , 1 mM EGTA).
 - 894.6 mg KCl
 - 68.0 mg KH_2PO_4
 - 238.8 mg HEPES
 - 19.0 mg MgCl_2
 - 38.0 mg EGTA
 - pH 7.2 (adjusted with KOH)
- 2 Prepare 10 mM Amplex Red reagent stock solution.
 - 2.1 Warm up Amplex Red reagent (blue cap) and DMSO (green cap) at RT.
 - 2.2 Immediately before use, add 60 μL of DMSO into a vial of Amplex Red (each vial of Amplex Red is sufficient for 100 assays of 100 μL).
- 3 Prepare 1X reaction buffer by diluting 4 mL of 5X reaction buffer (white cap) with 16 mL of deionized water.
- 4 Make 10 U/mL horseradish peroxidase (HRP) stock solution.
 - 4.1 Reconstitute a vial of HRP (yellow cap) with 1 mL of 1X reaction buffer.
 - 4.2 After the assay, aliquot unused HRP stock solution into single-use tubes and store at -20°C .
- 5 Prepare 20 mM hydrogen peroxide (H_2O_2) working solution.
 - 5.1 Immediately before use, dilute the ~3% (0.88 M) H_2O_2 (red cap) into appropriate volume of 1X reaction buffer to make a 20 mM working solution; exact concentration is on the label.
- 6 Prepare respiratory substrate solution (10 mL; 50 mM pyruvate and 25 mM malate).



- 55.0 mg pyruvate
- 33.5 mg malate
- 10 mL of mitochondrial respiration buffer

Measure H₂O₂ level

- 7 Dilute 30 µg crude mitochondrial preps to 40 µL with mitochondrial respiration buffer.
- 8 Prepare positive (10, 5, 2, 1 and 0.5 µM H₂O₂) and negative (1X mitochondrial respiration buffer) controls.
- 9 Load samples (40 µL/well).
- 10 Prepare a working solution of 100 µM Amplex Red reagent and 0.2 U/mL HRP by mixing the following:
 - 50 µL of 10 mM Amplex Red reagent stock solution
 - 100 µL of 10 U/mL HRP stock solution
 - 4.85 mL of 1X reaction buffer
- 11 Add 50 µL of the working solution into each well.
- 12 Record fluorescence emission at 590 nm (excitation at 560 nm) for 8 min.
- 13 Add 10 µL of respiratory substrate into each well; final concentration is 5 mM pyruvate and 2.5 mM malate.
- 14 Record fluorescence emission at 590 nm (excitation at 560 nm) for 45 min.