

Jul 11, 2024

MLi-2 treatment of astrocytes

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkwzzwl5r/v1>

Shiyi Wang¹

¹Duke University



Shiyi Wang

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkwzzwl5r/v1>

Document Citation: Shiyi Wang 2024. MLi-2 treatment of astrocytes. protocols.io
<https://dx.doi.org/10.17504/protocols.io.j8nlkwzzwl5r/v1>

License: This is an open access document distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: May 24, 2023

Last Modified: July 11, 2024



Document Integer ID: 82325

Keywords: ASAPCRN, Irrk2 inhibitor

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-020607

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

LRRK2 inhibitor MLI-2 treatment of astrocytes



1. Cortical astrocytes were isolated from P0-P1 rat female pups as described in the “**Cortical astrocyte isolation and culture**” section.
2. Cultured astrocytes were treated with 0, 30, and 100 nM LRRK2 kinase inhibitor, MLI-2 (Tocris; Catalog number: 5756), for 72 hours.
3. Cultured astrocytes were lysed at DIV 11 for protein extraction and Western blot analysis.