



May 22, 2024

Modelling Insulin resistance in vitro

DOI

<https://dx.doi.org/10.17504/protocols.io.kxygxy3mwl8j/v1>

Laura Mahoney-Sanchez¹

¹The Francis Crick Institute, 1 Midland Road, London NW1 1AT, UK.



Laura Mahoney-Sanchez

The Francis Crick Institute, 1 Midland Road, London NW1 1AT,...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygxy3mwl8j/v1>

Protocol Citation: Laura Mahoney-Sanchez 2024. Modelling Insulin resistance in vitro. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.kxygxy3mwl8j/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 22, 2024



Last Modified: May 22, 2024

Protocol Integer ID: 100256

Keywords: ASAPCRN, modelling insulin resistance, insulin resistance in ipsc, minutes insulin treatment the following day, cells of insulin, minutes insulin treatment, insulin, derived neuron, ipsc, neuron

Abstract

This protocol contains the instruction for modelling insulin resistance in iPSC-derived neurons by depriving the cells of insulin overnight followed by a 30 minutes insulin treatment the following day.



- 1 The media is removed from the cells
- 2 Cells are washed 1x with PBS
- 3 Fresh media is added the cells supplemented with B27 without insulin (Thermofisher A1895601).
- 4 Cells are left back in the incubator  Overnight at  37 °C and 5% CO2 18h
- 5 The following day, insulin was added to the media at either  100 nanomolar (nM) ,  250 nanomolar (nM) or  1 micromolar (μM) for  00:30:00 30m