

Mar 14, 2023

Version 2

Modified Promega Wizard Extraction for Barcoding Macrofungi V.2

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzb3p4vx1/v2>


Stephen Douglas Russell¹

¹The Hoosier Mushroom Society



Stephen D Douglas Russell

Mycota Lab, Biodiverse, MycoMap

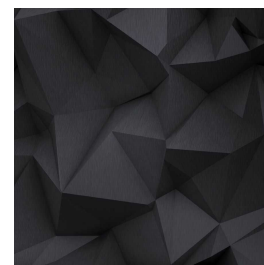

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzb3p4vx1/v2>

Protocol Citation: Stephen Douglas Russell 2023. Modified Promega Wizard Extraction for Barcoding Macrofungi. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.rm7vzb3p4vx1/v2> Version created by **Stephen D Douglas Russell**

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: March 13, 2023

Last Modified: March 14, 2023

Protocol Integer ID: 78721

Keywords: preparing macrofungal specimen, secondary extraction protocol for ont nanopore, macrofungal specimens for sanger, dna extraction method, dna extraction, modified promega wizard extraction, general fungal collection, nanopore barcoding, result with nanopore barcoding, quick extraction protocol, extraction protocol, fungi, secondary extraction protocol, extraction, ont nanopore, important specimen, specimen

Abstract

'This protocol is best used when preparing macrofungal specimens for Sanger sequencing or as a secondary extraction protocol for ONT nanopore barcoding.

The quality of a DNA extraction method is a primary limiting factor in the total number of samples that will return a result with nanopore barcoding of fungi. The "quick" extraction protocol will often yield a positive result for 80-85% of general fungal collections. Utilizing this extraction protocol pushes that number to nearly 100%. It is more time consuming and utilizes more expensive chemicals, but may be worth considering for important specimens that fail with the quick extraction protocol.



Materials

Equipment:

Tube Racks for 1.5uL eppi tubes

Tweezers

Pestles

Heat Block

Vortexer

Centrifuge

Consumables:

1.5uL eppi tubes

Molecular water

70% ethanol

Kimwipes

Reagents:

⊗ Nuclei Lysis Solution, 1000ml **Promega Catalog #A7943**

⊗ Protein Precipitation Solution 350ml **Promega Catalog #A7953**

⊗ Isopropanol **IBI Scientific**

Protocol materials

⊗ Nuclei Lysis Solution, 1000ml **Promega Catalog #A7943**

⊗ Protein Precipitation Solution 350ml **Promega Catalog #A7953**

⊗ Isopropanol **IBI Scientific**



- 1 Add 600uL of  Nuclei Lysis Solution, 1000ml **Promega Catalog #A7943** to 1.5mL eppi tubes. One tube for each specimen you are planning an extraction for.
- 2 Place tissue from your specimens into each tube using tweezers. Utilize a piece about the size of a grain of rice. The tissue can be either fresh or dried. Label the tube with the appropriate number. Wipe the tweezers off with a Kimwipe or paper towel in between each specimen. These tubes can be stored at room temperature until they are ready to be used.
- 3 Grind the tissue in each tube using a sterile pestle.

- 4 Heat the tubes at  65 °C for  00:15:00 .

15m

- 5 Centrifuge the tubes for  00:03:00 .

3m

- 6 Transfer the supernatant (liquid on top) to a new 1.5mL eppi tube.

6m 20s

Add  200 µL of

 Protein Precipitation Solution 350ml **Promega Catalog #A7953** to the tube.

Vortex the tube for  00:00:20 .

Centrifuge the tube for  00:06:00 .

- 7 Transfer the supernatant (liquid on top) to a new 1.5mL eppi tube.

1m

Add  600 µL of 100%  Isopropanol **IBI Scientific** to the tube. This precipitates the DNA.

Centrifuge the tube for  00:01:00 . The DNA will now be in a pellet stuck to the bottom of the tube.

Discard the supernatant. It can just be poured out of the tube into a waste container.



8 Add  600 μ L of 70% ethanol to the tube.

16m

Centrifuge the tube for  00:01:00 .

Discard the supernatant. It can just be poured directly out of the tube into a waste container.

Place the tube upside down on a Kimwipe for at least  00:15:00 , or until all of the ethanol has evaporated from the tube. I usually leave the tube to dry overnight.

9 Add 30uL of molecular water to the tube.

Your DNA template is now ready for amplification.