



MojoSort™ Isolation Kits Column Protocol - 1

DOI

dx.doi.org/10.17504/protocols.io.7bahiie



Sam Li¹

¹BioLegend

BioLegend

Tech. support email: tech@biolegend.com



Sam Li

BioLegend



DOI: dx.doi.org/10.17504/protocols.io.7bahiie

External link: <https://www.biolegend.com/protocols/mojosort-isolation-kits-column-protocol-1/4589/>

Protocol Citation: Sam Li . MojoSort™ Isolation Kits Column Protocol - 1. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.7bahiie>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: September 12, 2019

Last Modified: September 12, 2019

Protocol Integer ID: 27714

Keywords: MojoSort™, nanobeads, column, cell separation

Abstract

BioLegend MojoSort™ nanobeads work in commonly used separation columns, based on our internal research as well as validation by external testing by academic labs. This simple protocol consists of following the MojoSort™ protocol to label the cells with **pre-diluted** MojoSort™ reagents and using the columns as indicated by the manufacturer.

Note: Due to the properties of our beads, it may be possible to use far fewer beads and less antibody cocktail than with other commercial suppliers. We recommend a titration to find the best dilution factor. However, as a general rule, dilutions ranging from 1:2 to 1:10 for the antibody cocktail can be used. Dilutions ranging from 1:5 to 1:20 for the Streptavidin Nanobeads can be used. Please contact BioLegend Technical Service (tech@biolegend.com) if further assistance is needed.

Guidelines

MojoSort™ magnetic particles can be used with other commercially available magnetic separators, both free standing magnets and column-based systems. Because MojoSort™ protocols are optimized for the MojoSort™ separator, the protocols may need to be adjusted for other systems. Please contact BioLegend Technical Service (tech@biolegend.com) for more information and guidance. We do not recommend using MojoSort™ particles for BD's IMag™ or Life Technologies' DynaMag™.

Materials

MATERIALS

 MojoSort™ Buffer **BioLegend Catalog #480017**

Additional reagents:

- commercially available cell separation columns
- 5 mL polypropylene tubes
- 70 µm cell strainer

This protocol works with the following MojoSort™ Kits (cat#):

480005, 480006, 480033, 480007, 480008, 480035, 480009, 480010, 480021, 480022, 480023, 480024, 480031, 480039, 480040, 480041, 480042, 480043, 480044, 480045, 480046, 480051, 480052, 480061, 480062, 480063, 480064, 480067, 480068, 480081, 480082, 480087, 480088

- 1 Prepare cells from your tissue of interest or blood without lysing erythrocytes. Kits for human samples have been optimized for PBMCs, please prepare the cells using a suitable method.
- 2 In the final wash of your sample preparation, resuspend the cells in MojoSort™ Buffer by adding up to 4 mL in a 5 mL (12 × 75 mm) polypropylene tube.
Note: Keep MojoSort™ Buffer on ice throughout the procedure.
- 3 Filter the cells with a 70 µm cell strainer, centrifuge at 300xg for 5 minutes, and resuspend in an appropriate volume of MojoSort™ Buffer. Count and adjust the cell concentration to 1×10^8 cells/mL.
- 4 Aliquot 100 µL (10^7 cells) into a new tube. **Add 10 µL of the pre-diluted Biotin-Antibody Cocktail.** Mix well and **incubate on ice for 15 minutes.** Scale up the volume if separating more cells. For example, add 100 µL of pre-diluted antibody cocktail for separating 1×10^8 cells in 1 mL of MojoSort™ Buffer. When working with less than 10^7 cells, use indicated volumes for 10^7 cells.
- 5 Vortex the Streptavidin conjugated Nanobeads (to resuspend) at max speed, 5 touches, and prepare the dilutions to test. **Add 10 µL of pre-diluted Streptavidin Nanobeads.** Mix well and **incubate on ice for 15 minutes.** Scale up the volume accordingly if separating more cells. For example, add 100 µL of pre-diluted Nanobeads for separating 1×10^8 cells in 1 mL of MojoSort™ Buffer. When working with less than 10^7 cells, use indicated volumes for 10^7 cells.
- 6 Add the appropriate amount of MojoSort™ Buffer and proceed to separation. At least 500 µL is needed for column separation.
Note: There are several types of commercially available columns, depending on your application. Choose the one that fits best your experiment:

	Max. number of labeled cells	Max. number of total cells	Cell suspension volume	Column rinse volume	Cell wash volume	Elution volume
Small Capacity	1×10^7	2×10^8	500µL for up to 10^8 cells	1ml	1 ml	1 ml
Medium Capacity	1×10^8	2×10^9	500µL for up to 10^9 cells	3ml	3 ml	5 ml
Large Capacity	1×10^9	2×10^{10}	500µL for up to 10^{10} cells	20-50ml	30 ml	20 ml

Example of magnetic separation with medium capacity columns:

- 7 Place the column in a magnetic separator that fits the column.



- 8 Rinse the column with 3 mL of cell separation buffer.
- 9 Add the labeled cell suspension in at least 500 μ L of buffer to the column through a 30 μ m filter and collect the fraction containing the unlabeled cells. These are the cells of interest; do not discard.
- 10 Wash the cells in the column **2 times** with 3 mL of buffer and collect the fraction containing the untouched cells. Combine with the collected fraction from step 3.
- 11 If desired, the labeled cells can be collected by taking away the column from the magnet and place it on a tube. Then add 5 mL of buffer and flush out the magnetically labeled fraction with a plunger or supplied device. The labeled cells may be useful as staining controls, to monitor purity/yield, or other purposes.