



MojoSort™ Mouse anti-PE Nanobeads Protocol

DOI

<https://dx.doi.org/10.17504/protocols.io.7xshpne>



Sam Li¹

¹BioLegend

BioLegend

Tech. support email: tech@biolegend.com



Sam Li

BioLegend

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account



DOI: <https://dx.doi.org/10.17504/protocols.io.7xshpne>

External link: <https://www.biolegend.com/protocols/mojosort-mouse-anti-pe-nanobeads-protocol/4864/>

Protocol Citation: Sam Li . MojoSort™ Mouse anti-PE Nanobeads Protocol. [protocols.io](https://dx.doi.org/10.17504/protocols.io.7xshpne)

<https://dx.doi.org/10.17504/protocols.io.7xshpne>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: October 04, 2019



Last Modified: October 04, 2019

Protocol Integer ID: 28370

Keywords: mojosort, anti-PE, nanobeads, selection, isolation, target cell, pe nanobeads protocol product description, cells per tube, nanobead, functional assay, antibody, conjugated antibody, cell, specific cell number, assay

Abstract

Product description and procedure summary: Target cells are positively selected or depleted by incubating the sample with an anti-human PE conjugated antibody, followed by incubation with magnetic anti-PE Nanobeads. The magnetically labeled fraction is retained by the use of a magnetic separator. These are the PE⁺ cells, do not discard them if those are the cells of interest. Some of the downstream applications include functional assays, gene expression, phenotypic characterization, etc.

Note: This procedure is optimized for the isolation of 10^7 to 2×10^8 cells per tube. If working with fewer than 10^7 cells, keep volumes as indicated for 10^7 cells. For best results, optimize the conditions to your specific cell number and tissue. Prepare fresh MojoSort™ Buffer solution by diluting the 5X concentrate with sterile distilled water. *Scale up volumes if using 14mL tubes and Magnet, and place the tube in the magnet for 10 minutes.*

Guidelines

MojoSort™ magnetic particles can be used with other commercially available magnetic separators, both free standing magnets and column-based systems. Because MojoSort™ protocols are optimized for the MojoSort™ separator, the protocols may need to be adjusted for other systems. Please contact BioLegend Technical Service (tech@biolegend.com) for more information and guidance. We do not recommend using MojoSort™ particles for BD's IMag™ or Life Technologies' DynaMag™.

Application notes: To use this product in magnetic separation columns, a titration of the Nanobeads should be performed. Optimal concentration for magnetic separation columns is lot-specific. Please contact BioLegend Technical Service (tech@biolegend.com) for further assistance on how to use MojoSort™ Nanobeads in magnetic separation columns.

Materials

MATERIALS

⊗ MojoSort™ Magnet **BioLegend Catalog #480019**

⊗ MojoSort™ Buffer **BioLegend Catalog #480017**

⊗ MojoSort™ Mouse anti-PE Nanobeads **BioLegend Catalog #480079, 480080**

- Adjustable pipettes
- 70µm filters (one per sample)
- 5mL (12 × 75mm) or 14mL (17 × 100mm) polypropylene tubes
- Reagents for sample preparation
- Reagents and instruments (Flow cytometer) to determine yield and purity



- 1 Prepare cells from your tissue of interest or blood without lysing erythrocytes.
- 2 In the final wash of your sample preparation, resuspend the cells in MojoSort™ Buffer by adding up to 4 mL in a 5 mL (12 × 75 mm) polypropylene tube.
Note: Keep MojoSort™ Buffer on ice throughout the procedure.
- 3 Filter the cells with a 70 µm cell strainer, centrifuge at 300xg for 5 minutes, and resuspend in an appropriate volume of MojoSort™ Buffer. Count and adjust the cell concentration to 1×10^8 cells/mL.
- 4 Aliquot 100µL of cell suspension (10^7 cells) into a new tube. Check the recommended usage for flow cytometric staining of the PE-conjugated antibody indicated in the antibody technical datasheet. Calculate the volume to stain 10^7 cells (or desired amount of cells). **Add the appropriate volume of PE-conjugated antibody** to the cell suspension, mix well and **incubate on ice for 15 minutes**.
Optional: Take an aliquot before adding the cocktail to monitor purity and yield.
- 5 Wash the cells by adding MojoSort™ Buffer up to 4mL. Centrifuge the cells at 300xg for 5 minutes.
- 6 Discard supernatant and resuspend in 100µL of MojoSort™ Buffer.
- 7 Resuspend the beads by vortexing, maximum speed, 5 touches. Add **10µL of PE Nanobeads**. Mix well and **incubate on ice for 15 minutes**. Scale up the volume accordingly if separating more cells. For example, add 100 µL of Nanobeads for separating 1×10^8 cells in 1 ml of MojoSort™ Buffer. When working with less than 10^7 cells, use indicated volumes for 10^7 cells.
- 8 Wash the cells by adding MojoSort™ Buffer up to 4 mL. Centrifuge the cells at 300xg for 5 minutes.
- 9 Discard the supernatant.
- 10 Add 2.5 mL of MojoSort™ Buffer.
Note: If you observe aggregates, filter the suspension. To maximize yield, you can disrupt the aggregates by pipetting the solution up and down.
- 11 Place the tube in the magnet for 5 minutes.
Optional: Take a small aliquot before placing the tube in the magnet to monitor purity and yield. Keep unused cells to be used as control or other applications if needed.

- 12 Pour out the unlabeled fraction. If these are your cells of interest, **DO NOT DISCARD**. Resuspend the labeled cells in 2.5 mL MojoSort™ Buffer.
- 13 Repeat steps 11-12 on the labeled fraction twice more for a total of **3 separations**. Pool the unlabeled fractions and keep the labeled cells. The fraction that is not of interest may be useful as staining controls, to monitor purity/yield, or other purposes.
Optional: Take a small aliquot to monitor purity and yield.

