

Feb 21, 2025

Version 3

Mouse Brain Perfusion and Flash Freezing V.3

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkodr6v5r/v3>

Allen Institute¹

¹Allen Institute

Allen Institute for Brain S...

Allen Institute for Neural...



Allen Institute

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkodr6v5r/v3>

Protocol Citation: Allen Institute 2025. Mouse Brain Perfusion and Flash Freezing . **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.j8nlkodr6v5r/v3>Version created by **Allen Institute**

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: February 21, 2025

Last Modified: February 21, 2025

Protocol Integer ID: 123226

Keywords: transcardial perfusion with hepe, transcardial perfusion, perfusion, freezing, flash freezing, fast brain dissection, sucrose cutting solution, hepes hereafter, hepe

Abstract

This protocol is used for transcardial perfusion with HEPES-Sucrose Cutting Solution (Referred as HEPES hereafter), followed by fast brain dissection and flash freezing.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



AF0138_Mouse_Brain_P.

⌵

664KB

