

Jul 10, 2024

Mouse brain slice electrophysiology

DOI

<https://dx.doi.org/10.17504/protocols.io.n92ldp44nl5b/v1>

Shiyi Wang¹

¹Duke University



Shiyi Wang

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldp44nl5b/v1>

Document Citation: Shiyi Wang 2024. Mouse brain slice electrophysiology. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.n92ldp44nl5b/v1>

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: May 23, 2023

Last Modified: August 13, 2024



Document Integer ID: 82335

Keywords: ASAPCRN, mouse, brain

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-020607

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

The protocols.io team notes that research involving animals and humans must be conducted according to internationally-accepted standards and should always have prior approval from an Institutional Ethics Committee or Board.

Abstract

Mouse brain slice electrophysiology

1. For whole-cell patch-clamp recordings, 3–4 mice of each genotype and condition were used for miniature excitatory postsynaptic current (mEPSC) and miniature inhibitory postsynaptic current (mIPSC) measurements.
2. WT and *LRRK2* G2019S^{ki/ki} mice of both sexes were anesthetized with 200 mg/kg tribromoethanol (avertin) and decapitated.
3. After decapitation, the brains were immersed in ice-cold artificial cerebrospinal fluid (aCSF, in mM): 125 NaCl, 2.5 KCl, 3 mM MgCl₂, 0.1 mM CaCl₂, 10 glucose, 25 NaHCO₃, 1.25 NaHPO₄, 0.4 L-ascorbic acid, and 2 Na-pyruvate, pH 7.3–7.4 (310 mOsmol).
4. 350 µm thick coronal slices containing the ACC were obtained using a vibrating tissue slicer (Leica VT1200; Leica Biosystems).
5. Slices were immediately transferred to standard aCSF (33°C, continuously bubbled with 95% O₂– 5% CO₂) containing the same as the low-calcium aCSF but with 1 mM MgCl₂ and 1–2 mM CaCl₂.
6. After 30-minute incubation at 33°C, slices were transferred to a holding chamber with the same extracellular buffer at room temperature (RT: ~25°C).
7. Brain slices were visualized by an upright microscope (BX61WI, Olympus) through a 40× water-immersion objective equipped with infrared-differential interference contrast optics in combination with a digital camera (ODA-IR2000WCTRL).
8. Patch-clamp recordings were performed by using an EPC 10 patch-clamp amplifier, controlled by Patchmaster Software (HEKA). Data were acquired at a sampling rate of 50 kHz and low-pass filtered at 6 kHz.
9. To measure mEPSCs, the internal solution contained the following (in mM): 125 K-gluconate, 10 NaCl, 10 HEPES, 0.2 EGTA, 4.5 MgATP, 0.3 NaGTP, and 10 Na-phosphocreatine, pH adjusted to 7.2 – 7.4 with KOH and osmolality set to ~ 300 mOsmol.
10. mEPSCs were measured in the aCSF bath solution containing 1 µM tetrodotoxin and 50 µM Picrotoxin at -70 mV in voltage-clamp mode.
11. To measure mIPSCs, the internal solution contained the following (in mM): 77 K-gluconate, 77 KCl, 10 HEPES, 0.2 EGTA, 4.5 MgATP, 0.3 NaGTP, and 10 Na-phosphocreatine, pH adjusted to 7.2 – 7.4 with KOH and osmolality set to ~ 300 mOsmol.
12. mIPSCs were measured in the aCSF bath solution containing 1 µM tetrodotoxin and 10 µM 6-cyano-7-nitroquinoxaline-2,3-dione (CNQX), and 50 µM D-2-amino-5-phosphonopentanoate (D-AP5) at -70 mV in voltage-clamp mode.



13. mEPSCs and mIPSCs recorded at -70 mV were detected using Minhee Analysis software (https://github.com/parkgilbong/Minhee_Analysis_Pack).

14. To analyze the frequency, events were counted over 5 minutes of recording.

15. To obtain the average events for each cell, at least 100 non-overlapping events were detected and averaged. The peak amplitude of the average mEPSCs was measured relative to the baseline current.