

Apr 04, 2025

Version 8

Mouse Cardiac Perfusion Fixation and Brain Collection V.8

DOI

<https://dx.doi.org/10.17504/protocols.io.8epv51bejl1b/v8>

Allen Institute for Brain Science¹

¹Allen Institute

Allen Institute for Brain Science
Tech. support
[Click here to message tech. support](#)



Allen Institute

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8epv51bejl1b/v8>

Protocol Citation: Allen Institute for Brain Science 2025. Mouse Cardiac Perfusion Fixation and Brain Collection. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.8epv51bejl1b/v8> Version created by **Allen Institute**

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 04, 2025

Last Modified: April 04, 2025

Protocol Integer ID: 126182

Keywords: AF0075, perfusion, brain, mouse, cardiac, mouse cardiac perfusion fixation, intracardiac perfusion fixation, procedures for intracardiac perfusion fixation, intracardiac perfusion, postnatal mice, brain removal, fixation, mice, brain collection this protocol, including anesthesia, brain collection,

Abstract

This protocol describes the procedures for intracardiac perfusion fixation of postnatal mice, including anesthesia, exsanguination, fixation, brain removal and post-fixation storage.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



LAAS Manual.pdf

1.9MB



557766_Pump_66,67_Ma

...
1MB



AF0075_Mouse_Cardiac

...
782KB

