


Aug 22, 2023

Mouse colon vasculature labeling combined with immunofluorescence

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvwddd7lmk/v1>

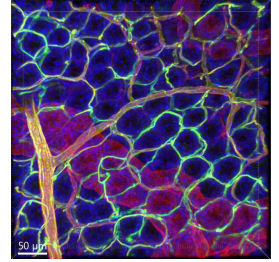

Lixin Wang¹

¹University of California, Los Angeles

Lixin Wang: We thank Ms. Honghui Liang (UCLA) for her excellent technical assistance.

SPARC

Tech. support email: info@neuinfo.org



Manmeet Bains

UCLA

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvwddd7lmk/v1>

Protocol Citation: Lixin Wang 2023. Mouse colon vasculature labeling combined with immunofluorescence. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.e6nvwddd7lmk/v1>

Manuscript citation:

Wang, L., P. Q. Yuan, C. Challis, S. Ravindra Kumar, and Y. Taché. 2022. "Transduction of Systemically Administered Adeno-Associated Virus in the Colonic Enteric Nervous System and c-Kit Cells of Adult Mice." *Front Neuroanat* 16: 884280. <https://doi.org/10.3389/fnana.2022.884280>.

Wang, L., P. Q. Yuan, and Y. Taché. 2023. "Vasculature in the mouse colon and spatial relationships with the enteric nervous system, glia, and immune cells." *Front Neuroanat* 17: 1130169. <https://doi.org/10.3389/fnana.2023.1130169>.

Yang, B., J. B. Treweek, R. P. Kulkarni, B. E. Deverman, C. K. Chen, E. Lubeck, S. Shah, L. Cai, and V. Gradinaru. 2014. "Single-cell phenotyping within transparent intact tissue through whole-body clearing." *Cell* 158 (4): 945-958. <https://doi.org/10.1016/j.cell.2014.07.017>. <https://www.ncbi.nlm.nih.gov/pubmed/25088144>.

Yuan, P. Q., J. P. Bellier, T. Li, M. R. Kwaan, H. Kimura, and Y. Taché. 2021. "Intrinsic cholinergic innervation in the human sigmoid colon revealed using CLARITY, three-dimensional (3D) imaging, and a novel anti-human peripheral choline acetyltransferase (hpChAT) antiserum." *Neurogastroenterol Motil* 33 (4): e14030. <https://doi.org/10.1111/nmo.14030>.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 16, 2023

Last Modified: August 22, 2023

Protocol Integer ID: 86562

Keywords: colon, fluorescent vessel painting, immunohistochemistry, microvessel, mouse, vasculature, wheat germ agglutinin, mouse colon vasculature labeling, different mouse colonic segment, spatial relationships of the vasculature, innervation of vasculature, immunofluorescence, enteric plexus, macrophage, innervation

Funders Acknowledgements:

NIH

Grant ID: OT2OD024899

Abstract

To study the distribution, morphology and innervation of vasculature in different mouse colonic segments and layers, as well as spatial relationships of the vasculature with the enteric plexuses, glia and macrophages

Materials

Animals

Mice, C57BL/6J (000664, Jackson Laboratories, Sacramento, CA), male and female 8-12 weeks old

Reagents and supplies

1. Isoflurane
2. 0.9% saline
3. wheat germ agglutinin-Alexa Fluor (WGA-AF) 448 (W11261, ThermoFisher Scientific)
4. 0.01 M phosphate-buffer saline (PBS), stock solution 10X
5. paraformaldehyde (P6148, Millipore-Sigma)
6. Sylgard™ 184 silicone elastomer (Electron Microscopy Science, Hatfield, PA)
7. Acrylamide
8. VA-044 Initiator
9. A4F0 hydrogel (Yang et al. 2014): 4% acrylamide and 0.25% VA-044 Initiator in 0.01M PBS. (A4F0: 4% acrylamide without paraformaldehyde)
10. Sucrose
11. Primary antibodies (Table 1)
12. Secondary antibodies: donkey anti-rabbit IgG 555 or 647, donkey anti-rat IgG 594 (all purchased from ThermoFisher)
13. Normal donkey serum
14. Triton-X 100
15. OCT Compound (Tissue-Plus™, 23-730-571, Fisher Scientific)
16. DAPI (4', 6-diamidino-2-phenylindole; 62247, ThermoFisher Scientific)
17. RIMS (refractive index matching solution) (Yang et al. 2014)
18. Vectashield anti-fade mounting medium (Vector Laboratories, Burlingame, CA)
19. iSpacer (Sunjin Lab, Hsinchu City, Taiwan, R.O.C.) (Yuan et al. 2021)

Table

1. Primary antibodies

	A	B	C	D	E	F
	Antibody	RRID*	Species	Source	Catalog No.	Dilution
	CD31	AB_393571	Rat	BD Biosciences	550274	1:50
	PGP9.5	AB_10891773	Rabbit	Abcam	ab108986	1:1,000
	αCGRP	AB_518147	Rabbit	Peninsula	T-4032	1:2,000
	VIP	AB_2890602	Rabbit	CURE/UCLA	ab7913	1:1,000
	TH	AB_390204	Rabbit	Millipore	AB152	1:1,000
	GFAP	AB_305808	Rabbit	Abcam	ab7260	1:2,000



	A	B	C	D	E	F
	S100B	AB_882426	Rabbit	Abcam	AB56242	1:1,000
	Iba1	AB_839504	Rabbit	WAKO	019-19741	1:1,000

*: RRID: Research Resource Identifiers

Tissue collection and preparation

1

- 1.1 Remove the whole colon is removed from the ileocecal junction to the end of distal colon at the level of pelvic brim where the iliac artery runs, after the cardiovascular perfusion of WGA.
- 1.2 Open the colon along the mesenteric margin, and pin the colon flat on a dish coated with Sylgard™ 184 silicone elastomer.
- 1.3 Fix the colon in 4% paraformaldehyde in 0.1M phosphate buffer (pH 7.4) overnight at 4°C.
- 1.4 Divide the colon samples to the proximal, mid and distal colon (Wang et al. 2022; Wang et al. 2023).
- 1.5 To improve penetration of antibodies for immunofluorescence, or for visualization of vasculature in different layers, the colonic samples could be prepared as the following approaches:
 - (1) Clear the samples using A4P0 hydrogel. The methodological details as described in previous report (Yang et al. 2014) and the protocol on protocols.io (Yuan et al. <https://dx.doi.org/10.17504/protocols.io.bqi2muge>).
 - (2) Treat the samples with 2% Triton-X 100 and 10% normal donkey serum in 0.01M PBS for 1 h at room temperature (RT) and then overnight at 4°C.
 - (3) Because the layers of the proximal colon with mucosal folds are thick and poorly immunolabeled, the proximal colon samples are prepared as: all layers without the mucosa and separating the wall in two parts at the submucosal layer.
 - (4) Frozen thick sections of colonic walls. The colon of mice perfused with the paraformaldehyde immersed in 20% sucrose for 2 days for cryoprotection, and embedded in Tissue-Plus™ O.C.T. Compound, snap-frozen and sectioned at 200 µm in a cryostat.

Vasculature painting

2

- 2.1 Dilute WGA AF-488 in 0.1M phosphate-buffer saline (PBS) at 30 µg/ml.

2.2 Anesthetized mice deeply with 5% isoflurane.

2.3 Flush the blood system with 0.9% saline via a cardiac cannula

2.4 Perfuse WGA-AF 488 (30 µg/g of body weight) at 1 ml/g of mouse body weight and a rate of 3 ml/min.

For immunostaining in thick colonic wall sections, some mice were perfused with 4% paraformaldehyde in 0.1M phosphate buffer (pH 7.4) following WGA perfusion.

Immunohistochemistry

3 Process samples of mice perfused with WGA-AF 488 by the following steps:

3.1 2% Triton-X 100 and 10% normal donkey serum in 0.01M PBS for 1 h at room temperature (RT) and then overnight at 4°C (the same step of 5(2) in Tissue Preparation)

3.2 Primary antibodies (Table 1) in 0.3% Triton-X 100-PBS for 2 h RT, at 4°C 2-5 days

3.3 Secondary antibodies for 3 h at RT, then overnight at 4°C.

3.4 Counterstaining: DAPI 1:2,000, 30 min at RT.

3.5 Mount the colonic tissues onto glass slides in a frame (iSpacer), and seal by glass coverslip in Vectashield anti-fade mounting medium or RIMS

Imaging and quantitative analysis

4

4.1 Acquire microscopic images were acquired with Z-stacks in Zeiss confocal microscopes (LSM 710 and 880, Carl Zeiss Microscopy GmbH, Germany), using objectives of 10X, 20X



and 63X.

4.2 The optical sections were scanned at intervals of 2 or 2.5 μm in 10X, 1 μm in 20X and 0.5 μm in 63X objectives.

4.3 The image segmentation, quantitation and visualization using Imaris 9.8 and 9.9 for neuroscientists (Bitplane Inc., Concord, MA).

4.4 The image segmentation, quantitation and visualization using Imaris 9.8 and 9.9 for neuroscientists (Bitplane Inc., Concord, MA).