



Aug 02, 2023

Mouse perfusions and brain tissue processing

DOI

<https://dx.doi.org/10.17504/protocols.io.bp2l69kn1lqe/v1>

Ayse Ulusoy¹, Shirley Lee¹, Angela Rollar¹, Michael Helwig¹, Michael Klinkenberg¹, Sinead O'Sullivan¹, Rita Pinto-Costa¹, Donato Di Monte¹

¹DZNE



Ayse Ulusoy

German Center for Neurodegenerative Diseases

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.bp2l69kn1lqe/v1>

Protocol Citation: Ayse Ulusoy, Shirley Lee, Angela Rollar, Michael Helwig, Michael Klinkenberg, Sinead O'Sullivan, Rita Pinto-Costa, Donato Di Monte 2023. Mouse perfusions and brain tissue processing. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.bp2l69kn1lqe/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: August 25, 2022

Last Modified: May 31, 2024

Protocol Integer ID: 69172

Keywords: ASAPCRN, mouse perfusion, harvesting from perfused rodent, brain tissue, perfused rodent, perfusion, brain fixation, processing, tissue, mouse, histological analysis, brain

Abstract

The protocol describes tissue processing for histological analyses and includes perfusions, brain fixation, and tissue harvesting from perfused rodents.



- 1 Inject sodium pentobarbital (600 mg/kg, i.p.) to sacrifice the mouse
- 2 Once the respiration is stopped, open the chest and place the cannula connected to a peristaltic pump into the heart. Make sure that the cannula has access to the ascending aorta.
- 3 Perfuse, first with 20 ml saline solution, kept at room temperature, and then with ice-cold 60 ml 4% (w/v) paraformaldehyde.
- 4 Quickly remove the brain and immerse in 4% paraformaldehyde for 24 hours.
- 5 For cryopreservation, place the brain in a 30% (w/v) sucrose solution.
- 6 Once sunk, section the brain on the coronal plane ( 35 μ L) using a freezing microtome.