

Jul 13, 2019

## Mouse Stellate Immunohistochemistry protocol

DOI

[dx.doi.org/10.17504/protocols.io.2qggdtw](https://dx.doi.org/10.17504/protocols.io.2qggdtw)

Pradeep Rajendran<sup>1</sup>

<sup>1</sup>[University of California, Los Angeles]



Scott John

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.2qggdtw>

**Protocol Citation:** Pradeep Rajendran 2019. Mouse Stellate Immunohistochemistry protocol. **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.2qggdtw>

**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** May 09, 2019



**Last Modified:** July 13, 2019

**Protocol Integer ID:** 23016

**Keywords:** immunohistochemistry protocol, mouse

## Troubleshooting

- 1 See mouse stellate isolation protocol
- 2 Complete/intact stellate ganglia are fixed in 4% paraformaldehyde overnight at 4°C.
- 3 Fixed tissue is rinsed in phosphate buffered saline (PBS), and stored in PBS + 0.02% sodium azide.
- 4 The stellate ganglion whole mounts are permeabilized in 'block' solution. (10% normal donkey serum in 1.0% BSA + 0.4% Triton X-100 + PBS).
- 5 Ganglia are stained with antibodies directed to either tyrosine hydroxylase (TH) or protein gene product 9.5 (PGP 9.5). Secondary staining with streptavidin conjugated to ATTO-647N was used to visualize neurobiotin filling. Stained tissue was rinsed in PBS and mounted on glass slides.