

Jul 10, 2024

Mouse synapse imaging and analysis

DOI

<https://dx.doi.org/10.17504/protocols.io.kxygx9jjog8j/v1>

Shiyi Wang¹

¹Duke University



Shiyi Wang

Duke University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygx9jjog8j/v1>

Document Citation: Shiyi Wang 2024. Mouse synapse imaging and analysis. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.kxygx9jjog8j/v1>

License: This is an open access document distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Created: May 23, 2023

Last Modified: July 10, 2024



Document Integer ID: 82332

Keywords: ASAPCRN, analysis mouse synapse imaging, imaging, mouse

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP) initiative

Grant ID: ASAP-020607

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Mouse synapse imaging and analysis

1. Staining, image acquisition, and analysis of synaptic density were performed as described previously with slight adjustments.
2. Synaptic staining was performed in coronal sections (30 μm thick) containing the ACC and MOp from WT and LRRK2 G2019S^{ki/ki} mice.
3. To label pre and postsynaptic proteins, the following antibody combinations were used: VGluT1 and PSD95 (excitatory, intracortical), VGAT, and GEPHYRIN (inhibitory).
4. High magnification 60 $\times$ objective plus 1.64x optical zoom z-stack images containing 15 optical sections spaced 0.34 μm apart were obtained using an Olympus FV 3000 inverted confocal microscope.
5. Each z-stack was converted into 5 maximum projection images (MPI), each corresponding to a 1 μm section of z plane, by using FIJI.
6. Synapses were identified by the colocalization of pre and postsynaptic puncta. The number of co-localized synaptic puncta of excitatory intracortical (VGluT1/PSD95) and inhibitory (VGAT/GEPHYRIN) synapses were obtained using the FIJI plugin Puncta Analyzer.
7. 15 MPIs were analyzed for each mouse (5 MPI/tissue section, 3 tissue sections/mouse).
8. Between 4 and 5, age and sex-matched mice/genotype/condition were analyzed for each synaptic staining combination, as indicated in the figure legends.
9. All animals appeared healthy at the time of collection. No data were excluded.