

Nov 16, 2018

Version 2

Mouse Tissue Fixation with Paraformaldehyde for Fluorescent Reporter Mice V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.vnge5bw>

Paola Marcovecchio¹, Sara McArdle¹, Zbigniew Mikulski¹, Angela Denn², katarzyna dobaczewska²

¹La Jolla Institute for Immunology; ²La Jolla Institute for Allergy and Immunology

Optical Clearing of Tissue

La Jolla Institute Micros...



Paola Marcovecchio

La Jolla Institute for Immunology

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.vnge5bw>

Protocol Citation: Paola Marcovecchio, Sara McArdle, Zbigniew Mikulski, Angela Denn, katarzyna dobaczewska 2018. Mouse Tissue Fixation with Paraformaldehyde for Fluorescent Reporter Mice. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.vnge5bw>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 15, 2018

Last Modified: November 16, 2018

Protocol Integer ID: 17832

Keywords: Fixation, Mouse, Tissue preparation, Histology, Microscopy, fluorescent reporter mice fixation of mouse tissue, fluorescent reporter mice fixation, mice with transgenic fluorescent reporter, tissue clearing for whole tissue imaging, cryosections for fluorescence microscopy, fluorescent reporter, whole tissue imaging, transgenic fluorescent reporter, fluorescence microscopy, mouse tissue, signal of the fluorescent reporter, tissue histology, tissue preparation, tissue clearing, fluorescence, essential first step in tissue preparation, perfusion of mice, microscopy, important that the tissue, tissue, histology, cryosection, paraformaldehyde, mice

Abstract

Fixation of mouse tissues for either histology or fluorescence microscopy is an essential first step in tissue preparation. In this protocol, we outline a reliable method for fixation and perfusion of mice for tissue histology or microscopy. Tissues prepared by this method can be used downstream as FFPE sections, cryosections for fluorescence microscopy, or tissue clearing for whole tissue imaging. For mice with transgenic fluorescent reporters, it is especially important that the tissues are preserved properly to ensure that the signal of the fluorescent reporter can be imaged.

Materials

MATERIALS

✂ 50 ml conical tubes

✂ PBS 1X

✂ EDTA Invitrogen - Thermo Fisher Catalog #AM9261

✂ 10mL syringe VWR International (Avantor) Catalog #75846-756

✂ 21G needle VWR International (Avantor) Catalog #BD-305165

✂ Forceps (Dumont) Braintree Scientific Catalog #FC-5043

✂ Forcep (Adson) Braintree Scientific Catalog #FC02 8

✂ Surgical scissors (Iris - straight) Braintree Scientific Catalog #SCT-I 528

✂ Paraformaldehyde 32% (methanol free) Electron Microscopy Sciences Catalog #15714



- 1 Prepare 1.6% PFA in PBS (at least  20 mL per mouse, half to be used during perfusion, half to be used as storage buffer overnight) MAKE THIS FRESH, DO NOT USE ANYTHING OLDER THAN 24 HOURS. Also, make sure PFA has been stored away from light, this will degrade it.
- 2 Prepare PBS with 2mM EDTA (at least  10 mL per mouse).
- 3 Sacrifice the mouse with CO₂. As soon as the mouse has stopped breathing, begin the procedure.
- 4 Dissect open the thoracic cavity, taking care not to puncture the heart with the scissors or forceps.
- 5 Fill a 10mL syringe with  10 mL PBS with 2mM EDTA , attach needle to syringe.
- 6 Insert needle into mouse's right ventricle (left side of heart when looking at the dissected mouse from your perspective). Slowly perfuse mouse until lungs turn white and liver has turned brown (instead of red).
- 7 Fill the syringe with  10 mL 1.6% PFA in PBS . Repeat perfusion. Mouse should appear stiff at the end of this perfusion.
- 8 Remove tissue of interest carefully so that marks are not made in the tissue.
- 9 Place tissue in  10 mL remaining 1.6% PFA in PBS in either 15 or 50mL conical (depending on how large the tissue is).
- 10 For larger, thicker tissues, leave at room temperature overnight ( 16:00:00 -  20:00:00) in a dark place or covered well in foil. For smaller tissues (i.e. lymph nodes) or for tissues that require antibody staining, fix at room temperature for a few hours (2-8, depending on downstream application).



- 11 Wash with PBS 1x the next day, 3 times. Use at least  10 mL (or more) to wash the tube with the fixed tissue in it. Let tissue sit in fresh PBS wash about  00:20:00 each time in Place in  4 °C until ready to use.