

Nov 27, 2025

MTT reagent preparation

DOI

<https://dx.doi.org/10.17504/protocols.io.4r3l21wz4g1y/v1>

Paolo Edoardo Ginatta Arosemena¹

¹Microbiology and Parasitology Department, University of Navarra

Parasitolab



Paolo Edoardo Ginatta Arosemena

Microbiology and Parasitology Department, University of Nava...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.4r3l21wz4g1y/v1>

Protocol Citation: Paolo Edoardo Ginatta Arosemena 2025. MTT reagent preparation. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.4r3l21wz4g1y/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 27, 2025

Last Modified: November 27, 2025

Protocol Integer ID: 233691

Keywords: mtt reagent preparation prep for mtt preparation, mtt preparation, reagent preparation prep, mtt, viability assay, preparation

Abstract

Prep for MTT preparation for viability assays

Safety warnings

 The important thing is that the final concentration of the MTT solution [M] 5 mg/mL



- 1 Add the  1000 mg MTT powder to  200 mL of PBS
- 2 Sonicate the solution until dissolved
- 3 Filter with  0.22 μ m filter
- 4 Aliquote and freeze