

Nov 12, 2023

Version 1

Multi-electrode array (MEA) V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.j8nlkonj6v5r/v1>

Leonardo Parra-Rivas¹

¹University of California, San Diego



Leonardo A Parra-Rivas

University of California, San Diego

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.j8nlkonj6v5r/v1>

Protocol Citation: Leonardo Parra-Rivas 2023. Multi-electrode array (MEA). protocols.io

<https://dx.doi.org/10.17504/protocols.io.j8nlkonj6v5r/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: November 03, 2023

Last Modified: November 12, 2023

Protocol Integer ID: 90325

Keywords: mea

Abstract

Multi-electrode array (MEA)

- 1 One day before plating, 6-well MEA plates (Axion Biosystems, M384-tMEA-6W) were coated with poly-D-lysine (Millipore Sigma P6407-5MG) in borate buffer
- 2 ,then washed six times with sterile water and aspirated prior use.
- 3 150,000 cells in a 15 μ l drop were plated on a 6-well MEA plate. Cultures were maintained for 17-21 days at 37 °C in 5% CO₂
- 4 On the day of recording, MEA plates were equilibrated for 10 min on the pre-warmer reader at 37 °C in 5% CO₂
- 5 2 mins of firing activity was recorded at a 12.5 kHz sampling rate and filtered with a bandpass filter from 200 Hz to 2.5 kHz. Each MEA well contained 64 electrodes that detected changes in extracellular field potentials reflecting the spiking activity of neurons. Spike detection was performed with the adaptive threshold crossing algorithm with a threshold of 6 $\times$ standard deviation using the AxIS software version 2.4 (RRID:SCR_016308) <https://www.axionbiosystems.com/products/axis-software>. All spike trains were exported from AxIS software.