



Oct 01, 2025

MULTI-Seq Barcoding and Library Preparation

DOI

<https://dx.doi.org/10.17504/protocols.io.kxygx3xzkg8j/v1>

Annika Martin¹, Hanqin Li^{1,2}, Dirk Hockemeyer^{1,2}

¹University of California, Berkeley, Berkeley, CA 94720, USA;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 2081



Hanqin Li

UC Berkeley

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kxygx3xzkg8j/v1>

Protocol Citation: Annika Martin, Hanqin Li, Dirk Hockemeyer 2025. MULTI-Seq Barcoding and Library Preparation. [protocols.io https://dx.doi.org/10.17504/protocols.io.kxygx3xzkg8j/v1](https://dx.doi.org/10.17504/protocols.io.kxygx3xzkg8j/v1)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 11, 2024

Last Modified: October 01, 2025

Protocol Integer ID: 93408

Keywords: ASAPCRN, seq barcoding, crispr screening, 10x genomics user guide, chromium next gem single cell 3' reagent kit, library preparation this protocol, crispr, feature barcoding technology, library preparation, list of reagent, reagent, barcoding, seq, preparation, sequencing, genomic

Funders Acknowledgements:

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-000486

Aligning Science Across Parkinson's (ASAP)

Grant ID: ASAP-024409

Abstract

This protocol describes MULTI-Seq barcoding of hESCs and library preparation, it is based on McGinnis et. al. 2019. PMID: [31209384](https://pubmed.ncbi.nlm.nih.gov/31209384/) and the 10x Genomics user guide "Chromium Next GEM Single Cell 3' Reagent Kits v3.1 with Feature Barcoding technology for CRISPR Screening"

Protocol overview

A. Oligonucleotides

B. Sequencing

Initial notes

A list of reagents and relevant vendor information can be found in the table listed under the materials tab.

Attachments



[MULTI-Seq Barcoding...](#)

131KB

Materials

Reagents:

	Item	Vendor	Catalog Number
	DPBS – Calcium and Magnesium Free (PBS-CMF)	Corning	MT21031CV
	Bovine Serum Albumin	Millipore/Sigma	A3803-50G
	SPRIselect Bead-Based Reagent	Beckman Coulter	B23318
	Chromium Single Cell 3' Reagent Kits v3 with Feature Barcode technology for CRISPR Screening	10x Genomics	PN-1000075PN-1000079
	Isopropyl alcohol/ Isopropanol(2-propanol)	Fisher Scientific	MFCD00011674
	Kapa HiFi HotStart ReadyMix (2X)	Fisher Scientific	NC0295239
	Qiagen Buffer EB	Qiagen	19086
	SPRIselect Reagent Kit	Beckman Coulter	B23318

Oligonucleotides

1 **Oligonucleotides** (store in -20°C for large term storage):

- 50 µM Anchor and Co-Anchor
- 10 µM Barcode Oligos
- 10 µM MULTI-seq additive primer
- 10 µM Universal I5 primer
- 10 µM TruSeq RPI primers

Anchor LMO: 5'-TGGAATTCTCGGGTGCCAAGGgtaacgatccagctgtcact-{Lipid}-3'

Co-Anchor LMO: 5'-{Lipid}-AGTGACAGCTGGATCGTTAC-3'

Barcode Oligo: 5'-CCTTGGCACCCGAGAATTCCANNNNNNNNA30-3'

MULTI-seq Primer: 5'-CTTGGCACCCGAGAATTCC-3'

TruSeq RPIX:

5'-

CAAGCAGAAGACGGCATAACGAGATNNNNNNGTGACTGGAGTTCCTTGGCACCCGAGAA
TTCCA-3'

Universal I5:

5'-

AATGATACGGCGACCAACGAGATCTACACTCTTCCCTACACGACGCTCTTCCGATCT-3'

Make a 10X stock of anchor and barcode strands by mixing at 1:1 molar ratio in PBS-CMF to 2 µM concentration. Pipette to mix.

A	B
In total Per sample	22 µL
10 µM unique Barcode Oligo	4.4 µL
50 µM Anchor LMO	0.9 µL
PBS-CMF	16.7 µL

2 Make a 10X solution of the Co-Anchor in in PBS-CMF to a final concentration of 2 µM. Pipette to mix.

A	B
In total Per sample	22 µL
50 µM Co-Anchor LMO	0.9 µL
PBS-CMF	21.1 µL



- 3 Prepare 1% BSA in PBS-CMF and place On ice – at least 3mL per sample.
- 4 Wash cells with PBS-CMF twice.
 - 4.1 If adherent – rinse and aspirate on plate.
 - 4.2 If suspension cells – centrifuge for 00:05:00 at 200 x g to 300 x g in a 15mL conical tube, carefully aspirate the supernatant, and resuspend in PBS-CMF.
- 5 Dissociate or lift cells to obtain single cell suspension (will vary depending on sample type).
 - 5.1 It is crucial that cells are properly resuspended in buffer at single cell suspension before labeling. For standard cell types, I recommend straining cells through a 40µm mesh before counting and labeling to prevent heterogeneous labeling.

Note

If using Trypsin or similar which is inactivated by the addition of FBS or BSA, rinse the cell suspension with PBS-CMF twice as described in step 4b.

- 6 Strain single-cell suspension through a 40 µm cell-strainer and count cells.
- 7 Spin down ~500k cells (or fewer) for 00:05:00 at 200 x g to 300 x g in a 15mL conical tube. Carefully aspirate the supernatant.
- 8 Barcoding:
 - 8.1 Suspend cells in 180 µL of PBS-CMF
 - 8.2 Add 20 µL 10X Anchor:Barcode solution and pipette gently to mix.



8.3 Incubate  On ice for  00:05:00 .

5m

8.4 Add  20 μ L Co-Anchor solution and pipette gently to mix.

8.5 Incubate  00:05:00 longer.

5m

8.6 Add  1 mL of 1% BSA in PBS (ice cold).

9 Transfer each cell sample to a microcentrifuge tube on ice. Keep the labeled cells on ice for the remainder of procedure until starting the 10x workflow to prevent loss of barcodes after washing. The labeling step itself can be done on  On ice or up to  37 °C .

10 Centrifuge cells for  00:05:00 at  200 x g to  300 x g at  4 °C .
Remove supernatant and resuspend pellet in additional  1 mL of ice cold 1% BSA in PBS-CMF.

5m

11 Repeat step 10 for a total of two wash steps.

12 Filter cell suspension through a 40 μ m cell-strainer and count cells again.

13 Combine all samples at desired ratio and continue with scRNA-seq procedure according to 10x Genomics instructions for endogenous transcripts.

14 For each lane of 10X, follow the 10X workflow until cDNA amplification (Step 2.2 in the Chromium Single Cell 3' Reagent Kits v3 with Feature Barcode technology for CRISPR Screening protocol).

14.1 During post GEM-RT cleanup the aqueous layer will be cloudy due to the higher BSA concentration we use during our post-barcoding rinses. The BSA helps to quench excess LMO barcodes and limit off-target labeling. This does not cause any issues or negatively affect results.



15 Prepare the following cDNA amplification master mix (volumes per lane):

A	B
10X Amp Mix (PN2000047/2000103)	50 μ L
Feature cDNA Primers 1 (PN2000096)	15 μ L
2.5 μ M MULTI-seq primer	1 μ L

16 Add  65 μ L cDNA Amplification Reaction Mix to  35 μ L sample.

17 Perform cDNA amplification according to 10X workflow.

18 To prevent SPRI-bead saturation, add  100 μ L water to the  100 μ L cDNA amplification mixture. Pipette to mix and split into two tubes of  100 μ L volume each.

19 Vortex to resuspend the SPRIselect reagent. Add  60 μ L SPRIselect reagent (0.6X) to each sample and pipette mix 15x (pipette set to 150 μ L).

20 Incubate  00:05:00 at  Room temperature .

5m

21 Place on the 10x Magnetic Separator in the magnet(High) orientation as described in the 10x workflow until the solution clears.

22 Transfer and save  150 μ L supernatant per tube in a new tube strip without disturbing the pellet.

22.1 For each sample, you should have two  150 μ L supernatant aliquots. One will be used for feature barcode preparation according to the 10X protocol. One will be used for MULTI-seq library preparation.

23 Maintain at room temperature. DO NOT discard the transferred supernatant DO NOT discard the pellet.



- 24 Continue with 2.3A Pellet Cleanup as described in the 10x workflow until step 2.3A.vi.
- 25 Remove the tube from the magnet. Using  40.5 μL buffer EB, resuspend BOTH SPRI bead pellets per sample by pipetting each pellet up and down 15x to mix thoroughly, combining into a single tube and proceed with protocol for endogenous transcripts without change.
- 26 Use one  150 μL supernatant aliquot for step 2.3B Transferred Supernatant Cleanup as described in the 10x workflow and proceed to Feature Barcode Library prep without change.
- 27 Transfer remaining  150 μL supernatant to fresh 1.5 mL microcentrifuge tube. Add  260 μL SPRI beads and  180 μL 100% isopropanol (for a final ratio of 1.8X SPRI). Pipette mix 10 times, incubate at  Room temperature for  00:05:00 .
- 28 Place tube on magnetic rack and wait for solution to clear.
- 29 Remove and discard supernatant.
- 30 Wash beads twice on magnet with  500 μL of 80% ethanol and allow to stand for  00:00:30 between washes.
- 31 After second wash, briefly centrifuge beads and place back on magnetic rack.
- 32 Remove any remaining ethanol with P10 micropipette.
- 33 Air-dry beads on magnet for  00:02:00 . Do NOT exceed 2 minutes.
- 34 Remove from magnet, resuspend beads in  50.5 μL buffer EB and pipette mix thoroughly to resuspend.

5m

30s

2m



35 Incubate at 0 °C for 00:02:00 .

2m

36 Return to magnet and wait for solution to clear.

37 Transfer 50 µL supernatant to PCR strip tube, pipetting carefully to avoid transferring beads.

38 Quantify barcode DNA concentration using Qubit (typical range is 0.5 - 5 ng/µL).

39 For each lane, prepare the following PCR mix:

A	B
Kapa HiFi HotStart ReadyMix (2X)	26.25 µL
10 µM Universal I5 primer	2.5 µL
10 µM unique RPI primer (choose unique RPI for each sample from 10X lane)	2.5 µL
barcode DNA (volume based on concentration from Qubit)	3.5 ng
Nuclease-free water	To 50 µL

40 Perform library preparation PCR:

A	B
95 °C	5 min
98 °C	15 sec
60 °C	30 sec
72 °C	30 sec
Repeat steps 2-5	8-12 times
72 °C	1 min
4 °C	hold



- 41 Add  80 μL (1.6X) SPRI to each PCR product, pipette mix thoroughly.
- 42 Incubate at  Room temperature for  00:05:00 . 5m
- 43 Place tube on 10x Magnetic Separator in the magnet (HIGH) orientation, wait for solution to clear.
- 44 Remove and discard supernatant.
- 45 Wash beads twice on 10x Magnetic Separator in the magnet (HIGH) orientation with  200 μL of 80% ethanol and allow to stand for  00:00:30 between washes. 30s
- 46 After second wash, briefly centrifuge beads, and invert the 10x Magnetic Separator to place the tubes on the 10x Magnetic Separator in the magnet (LOW) orientation.
- 47 Remove any remaining ethanol with P10 micropipette.
- 48 Air-dry beads on magnet for  00:02:00 . Do NOT exceed 2 minutes. 2m
- 49 Remove from magnet, resuspend beads in  25 μL buffer EB and pipette mix thoroughly to resuspend.
- 50 Incubate at  Room temperature for  00:02:00 . 2m
- 51 Return to 10x Magnetic Separator in the magnet (LOW) orientation, wait for solution to clear, and transfer supernatant to PCR strip tube.
- 52 Quantify barcode library concentration by running 1ul diluted 1:5 on an Agilent Bioanalyzer High Sensitivity chip.

Expected result

Expected library size is approximately 175bp.



Sequencing:

- a. Barcodes can be sequenced independently or as fraction of endogenous cDNA library.
- b. Target 3000-5000 barcode reads per cell.

- 53 Deconvolute barcodes for analysis as described at <https://github.com/chris-mcginnis-ucsf/MULTI-seq>

Protocol references

McGinnis CS, Patterson DM, Winkler J, Conrad DN, Hein MY, Srivastava V, Hu JL, Murrow LM, Weissman JS, Werb Z, Chow ED, Gartner ZJ. MULTI-seq: sample multiplexing for single-cell RNA sequencing using lipid-tagged indices. Nat Methods. 2019 Jul;16(7):619-626. doi: 10.1038/s41592-019-0433-8. Epub 2019 Jun 17. PMID: 31209384; PMCID: PMC6837808.