

Feb 27, 2019

Neural Rosette Formation and Selection

 In 1 collection

DOI

dx.doi.org/10.17504/protocols.io.x86frze

Celeste M M. Karch¹, Rita Martinez¹, Jacob Marsh¹

¹Washington University in St Louis

Neurodegeneration Method Development Community
Tech. support email: ndcn-help@chanzuckerberg.com



Celeste M M. Karch

Washington University in St Louis



Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x86frze>

Protocol Citation: Celeste M M. Karch, Rita Martinez, Jacob Marsh 2019. Neural Rosette Formation and Selection. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.x86frze>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 18, 2019

Last Modified: February 27, 2019

Protocol Integer ID: 20478

Keywords: neural rosette formation, selection, formation

Attachments



IPSC CORTICAL

DIFFER...

179KB

Guidelines

This protocol is part of the [IPSC CORTICAL DIFFERENTIATION](#) collection.


This method should be performed using sterile technique.


Materials

Please refer to the attached full manuscript for required materials.





Troubleshooting





Safety warnings

 Please refer to the SDS (Safety Data Sheet) for information about hazards, and to obtain advice on safety precautions.

- 1 On Day 5 of neural aggregate formation, remove media (by pipetting) and carefully wash spheres with  100 μL of pre-warmed DMEM/F12. Repeat 2 times.






Note

Do not break apart spheres. Neural spheres are very delicate at this stage. An alternative approach is to remove  50 μL of spent media and wash with  50 μL DMEM/F12 . Add  50 μL fresh neural induction media . Transfer  100 μL of spheres and media to the new PLO/laminin-coated well. This approach will transfer more dead cells into the new well.

- 2 Remove the last wash and add  50 μL of neural induction media to each well.
- 3 Aspirate laminin from one well of the pre-coated plate. Using 200 μL sterile tips, carefully pipet up spheres from wells using  100 μL volume and transfer thirty-two spheres per well. Repeat above steps for the remaining wells. Incubate cells in  37 $^{\circ}\text{C}$, 5% CO_2 and 95% humidified chamber and distribute evenly by making a “T” motion.
- 4 After  24:00:00 , examine attached aggregates. Remove medium and replace with 2mls/well fresh neural induction medium daily.


Note

If some aggregates have not attached, carefully pipet out all medium and replace with 1ml/well fresh neural induction medium. Once 90-100% of aggregates attach, exchange medium daily with 2mls/well neural induction medium.

- 5 Monitor spheres daily under microscope for formation of neural rosette structures. Neural rosettes are ready to harvest when spheres have completely flattened and clusters are clearly visible (3-7 days after plating, line dependent).
- 6 Harvest neural rosettes by aspirating spent medium. Add  1 mL of pre-warmed DMEM/F12 to each well to remove unattached cells (repeat if necessary).
- 7 Add  1 mL of Neural Rosette Selection reagent to each well and incubate for up to  01:00:00 at  37 $^{\circ}\text{C}$ (check cells at  00:20:00 . Cells are typically collected



after 30–45 min incubation. Look for rosette structure to be rounding up without the disturbance of other surrounding cells).

- 8 Carefully remove Neural Rosette Selection reagent with a pipet, being careful not to disturb rosette clusters. Add  1 mL DMEM/F12 to each well, then using a p1000 detach rosette clusters by rinsing over them.
- 9 Transfer rosette material from 1 well into a 15 mL conical tube for cryopreservation of neural rosettes and from 2 wells into a separate 15 mL conical tube for neural progenitor expansion. Do not triturate clusters.

Note

To maintain a pure culture, it is best to leave some rosettes behind rather than collect all of the rosettes and additional cells.

- 10 Centrifuge rosette clusters at 750 rpm for  00:03:00 .