

Jan 10, 2024

Neuromelanin Processing and Image Acquisition

DOI

dx.doi.org/10.17504/protocols.io.e6nvwd799lmk/v1

Anastasia Filimontseva^{1,2}, Glenda Halliday^{1,2}, YuHong Fu^{1,2}

¹Brain and Mind Centre & Faculty of Medicine and Health School of Medical Sciences, The University of Sydney, Sydney, NSW 2050, Australia;

²Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD, 20815



Anastasia Filimontseva

Brain and Mind Centre & Faculty of Medicine and Health Schoo...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvwd799lmk/v1>

Protocol Citation: Anastasia Filimontseva, Glenda Halliday, YuHong Fu 2024. Neuromelanin Processing and Image Acquisition. protocols.io <https://dx.doi.org/10.17504/protocols.io.e6nvwd799lmk/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: January 10, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 93362

Keywords: ASAPCRN, neuromelanin processing, intracellular nm quantification

Abstract

Protocol for preparing post-mortem tissue for intracellular NM quantification.

Troubleshooting

Tissue Processing

- 1 Formalin-fixed paraffin-embedded tissue blocks from the pontine and midbrain regions were sectioned at 6 μ m, collected onto adhesive microscope slides and allowed to dry in the oven at 37°C for 48 hours.
- 2 Slides were incubated in the oven at 60°C for 1 hour to melt the paraffin.
- 3 To remove the paraffin, sections were submerged in HistoChoice Clearing Agent (Sigma-Aldrich, H2779) for 2 \times 7 minutes, followed by rehydration in decreasing ethanol concentrations (100% ethanol for 2 \times 3 minutes, 95% ethanol for 3 minutes, 70% ethanol for 3 minutes) and distilled H₂O for 3 minutes.
- 4 Slides were submerged in 1xPBS and coverslipped using Fluorescence Mounting Medium (DAKO, S302380-2).

Image Acquisition

- 5 Sections were scanned using 20x objective (NA=0.8) with pre-set focusing and exposure parameters for optimal NM signal quality with an automated Slide Scanner (SLIDEVIEW VS200, Tokyo, Japan).
- 6 Identical parameters were applied to each scanned section to ensure consistency in capturing NM.
- 7 Once the scan was completed, sections were submerged in 1xPBS to remove coverslips and continue with IF processing.