

Nov 07, 2024

Neuronal differentiation and MEA plating using an inducible NGN2 hiPSC-line

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg3q54ev25/v1>

Silke Frahm¹

¹Stem cell technology platform



Silke Frahm

Stem cell technology platform

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg3q54ev25/v1>

Protocol Citation: Silke Frahm 2024. Neuronal differentiation and MEA plating using an inducible NGN2 hiPSC-line. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.kqdg3q54ev25/v1>

**Manuscript citation:**

This protocol is an adaptation of the published Star protocol Peitz et al. 2020.

Citation

Peitz M, Krutenko T, Brüstle O (2020)
. Protocol for the Standardized Generation of Forward Programmed Cryopreservable Excitatory and Inhibitory Forebrain Neurons.

<https://doi.org/10.1016/j.xpro.2020.100038>

LINK

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: October 01, 2024

Last Modified: November 07, 2024


Protocol Integer ID: 111768

Keywords: Neuronal differentiation, Inducible NGN2 hiPSC-line, NGN2, excitatory neurons, MEA, neuronal differentiation, neuron, inducible ngn2, ngn2, mea plate

Abstract

This protocol details the neuronal differentiation and MEA plating using an inducible NGN2 hiPSC-line. Neurons generated using this protocol will start showing electrical activity 1-2 weeks after seeding on MEA plate.

Guidelines

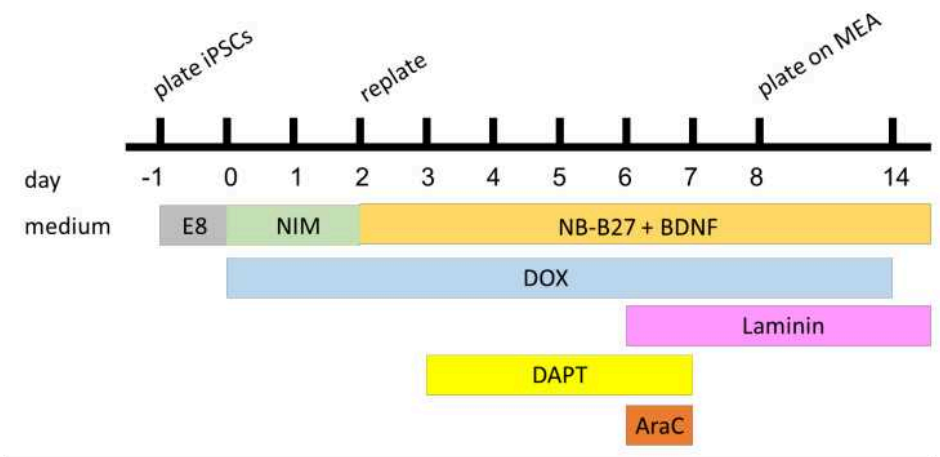

- Adapted from Peitz et al. 2020, Star protocol

Well numbers for plating 1 × 48WP MEA:

- Day -1:** 3 × 6W at 500k/well
- Day 2:** 4 × 6W at 750k/well



Materials

Neuronal Induction Medium (NIM)

- DMEM/F-12 500 mL
- N2 supplement 5 mL
- Doxycycline 2 µg/mL (final concentration)

Neuronal Medium (NB-B27)

- Neurobasal medium 500 mL
 - B-27 supplement (50x) 10 mL stable for one month at  4 °C
 - GlutaMAX (100x) 5 mL
- +
- BDNF 10 ng/mL (final concentration)
 - DAPT 10µM (1:1000)
 - Laminin 1:500 dilution
 - AraC (stock 4mM, use 1:800)
 - Doxycycline 2 µg/mL (final concentration) until day 14

BrainPhys Medium complete

-  BrainPhys™ Neuronal Medium **STEMCELL Technologies Inc. Catalog #05790**
- B-27 supplement 1:50
- Penn/Strep 1:100
- BDNF 10ng/ml
- Laminin 1:1000
- GDNF 10ng/ml
- Ascorbic acid 200µM
- cAMP 500µM

Reagents:

-  Pierce™ Concentrated Buffer Stocks (10X and 20X) **Thermo Scientific Catalog #28341**
-  Poly(ethyleneimine) solution **Merck MilliporeSigma (Sigma-Aldrich) Catalog #03880**
- mouse Laminin (Sigma-Aldrich) Catalog #L2020
- DMEM/F12 (Gibco) Catalog #11320-033
- Neurobasal (Gibco) Catalog #21103-049


Procedure

2h 40m

1 Day -1:

Dissociate 80% confluent hPSCs with Accutase or TrypLE and seed them on Geltrex-coated 6-well plates at a density of 450,000 cells/well in E8 supplemented with

[M] 10 micromolar (μM) RI ( 2 mL /well).

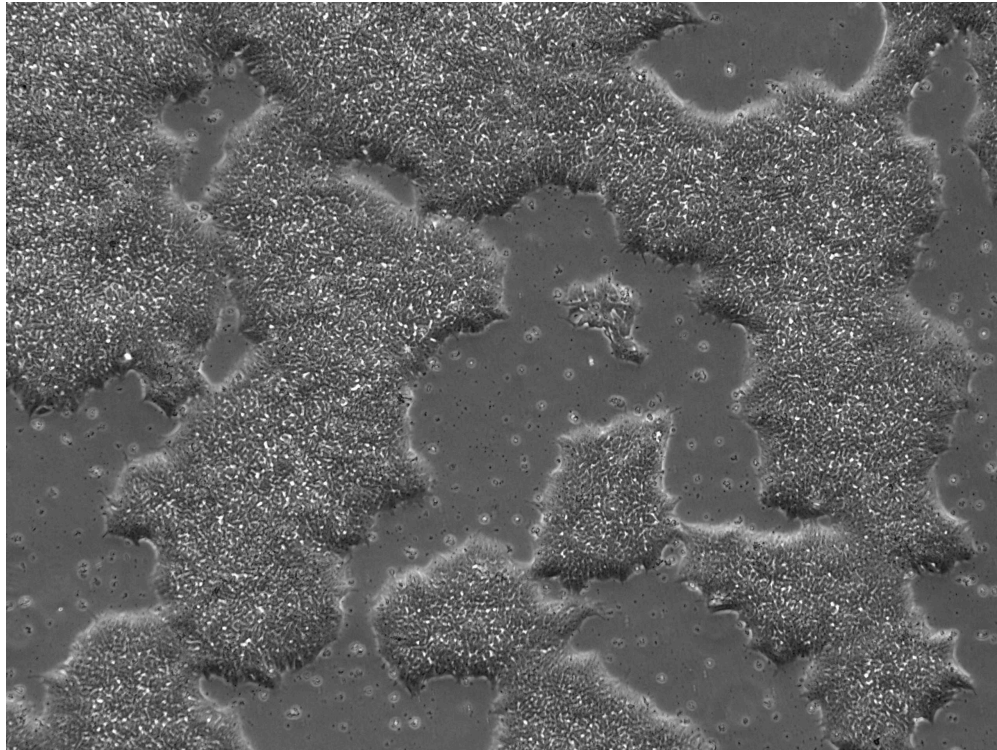

Example of a 80% confluent iPSC culture.

2 Day 0:

Change medium to neuronal induction medium (NIM) supplemented with  2 $\mu\text{g}/\text{ml}$ doxycycline (dox) to induce transgene expression. Dox is added freshly to the medium from now on. Start PO/laminin coating for neural progenitor split on day 2.

**3 Day 1:**

Change medium with NIM supplemented with  2 µg/ml doxycycline (dox).

4 Day 2:

Dissociate cells with Accutase for  00:10:00 at  37 °C . Plate cells on PO/laminin-coated wells at a density 7.5×10^4 cells/cm for the generation of glutamatergic neurons. At this point, change to neuronal medium supplemented with  2 µg/ml dox and  10 micromolar (µM) RI.

10m

**5 Day 3:**

Medium change (dox-supplemented neuronal medium containing  10 micromolar (µM) DAPT; DAPT remains in the medium from day 3 until day 7).

6 Day 5:

Medium change (dox-supplemented neuronal medium containing  10 micromolar (µM) DAPT). From this day on, carefully remove the medium with a serological pipet instead of an aspirator to avoid cell detachment.

**7 Day 6:**

Medium change (dox-supplemented neuronal medium containing  10 micromolar (µM) DAPT,  5 micromolar (µM) Ara-C and 1:500 laminin). Ara-C will induce a growth arrest of remaining proliferative cells.

8 Day 7:

Medium change (dox-supplemented neuronal medium containing 1:500 laminin). Pre-treat MEA plate and dry o/n.

9 Day 8: MEA plating:



- 9.1 Remove complete medium and directly add (no PBS wash) pre-warmed Accutase ( 1 mL /6W) supplemented with  10 micromolar (μM) RI, incubate for  01:00:00 at  37 °C .  1h
- 9.2 Wash off the cells with same amount of medium using a 10ml pipette, gently pipet up and down 5-10times, collect in 50ml Falcon tube.  
- 9.3 Use same amount of medium to wash once more the wells and add to tube (at the end should be double volume medium:accutase). 
- 9.4 Count cell number in suspension before spinning or after resuspending in max.  500 μL medium.
- 9.5 Spin down, resuspend in NB-B27 medium containing Laminin, BDNF, dox and RI at 50k cells per  8 μL (= 6,25Mio/ml).
- 9.6 Seed 8 μL as a drop with the pipet tip in the center of the MEA well covering the electrodes (48WP pretreated with PEI) without touching them and incubate for  00:15:00 -  00:30:00 .   45m
- 9.7 After 30min, add Astrocytes in  10 μL medium at a ratio of 1:5 (astrocyte:neuron) and incubate for another  00:15:00 -  00:30:00 (10.000 cells in  10 μL per well = 1M/ml).   30m
- 9.8 Add  300 μL medium after  00:30:00 .  30m 
- 9.9 Change medium next day (without RI) and half medium every other day thereafter.
- 9.10 After day 14, change gradually to BrainPhys medium complete (half medium changes 3x per week).

PO/Laminin coating

2d



10 Apply  15 µg/ml polyornithine (diluted in ddH₂O and then sterile-filtered) and incubate for  16:00:00 –  24:00:00 at  37 °C .

1d



11 Remove polyornithine and wash three times with PBS.



12 Apply laminin (diluted in DPBS++ 1:100) and incubate  Overnight for  16:00:00 –  24:00:00 at  4 °C .

1d



Note

Use  1.5 mL /6W

MEA coating

1h

13 Dot  8 µL 0.07% PEI solution over the recording electrodes.

14 Incubate  01:00:00 at  37 °C , 5% CO₂.

1h



15 Wash 3-4x with ddH₂O (PEI is toxic).



16 Air dry  Overnight in culture hood (no UV).



Preparing 0.07 % PEI solution

17 Dilute the 20x borate buffer to 1x in water (Thermo Fisher #28341).

18 Prepare a 0.07 % PEI solution in borate buffer using a 50 % PEI stock solution (Sigma: #03880), e.g. by diluting  28 µL of the 50%PEI stock solution in  20 mL 1x





Borate Buffer.

19 Filter the solution through a 0.22 µm filter, store  1 mL aliquots at  -20 °C .



Protocol references

This protocol is an adaptation of the published Star protocol Peitz et al. 2020.

Citations

Peitz M, Krutenko T, Brüstle O. Protocol for the Standardized Generation of Forward Programmed Cryopreservable Excitatory and Inhibitory Forebrain Neurons.

<https://doi.org/10.1016/j.xpro.2020.100038>