

Sep 27, 2019

Version 2

# Neuropathy Phentoying Protocols - Immunohistochemistry for Tissue Sections V.2

DOI

[dx.doi.org/10.17504/protocols.io.7rfhm3n](https://dx.doi.org/10.17504/protocols.io.7rfhm3n)



Eva Feldman<sup>1</sup>

<sup>1</sup>University of Michigan - Ann Arbor

Diabetic Complications Consortium

Tech. support email: [rmcindoe@augusta.edu](mailto:rmcindoe@augusta.edu)



Lili Liang

## Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.7rfhm3n>

External link: <https://www.diacomp.org/shared/document.aspx?id=54&docType=Protocol>



**Protocol Citation:** Eva Feldman 2019. Neuropathy Phentotyping Protocols - Immunohistochemistry for Tissue Sections.

**protocols.io** <https://dx.doi.org/10.17504/protocols.io.7rfhm3n>

**License:** This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working

**We use this protocol and it's working**

**Created:** September 27, 2019

**Last Modified:** September 27, 2019

**Protocol Integer ID:** 28167

**Keywords:** Neuropathy Phentotyping, Diabetic Neuropathy, Immunohistochemistry , diabetic neuropathy at the whole animal, presence of diabetic neuropathy, diabetic neuropathy, development of diabetic neuropathy, detailed measures of neuropathy, assessments of neuropathy, diabetes, neuropathy phentotyping protocol, duration of diabetes, neuropathy, diabetic complication, initial degree of neuropathy, degree of diabetes, month period of diabetes, loss of nerve fiber, tail blood glucose, degree of glycemic control, glycemic control, glycated hemoglobin level, nerve fiber, immunohistochemistry for tissue sections summary, immunohistochemistry, phenotyping of rodent

## Abstract

### Summary:

#### Phenotyping of Rodents for the Presence of Diabetic Neuropathy

In man, the development of diabetic neuropathy is dependent on both the degree of glycemic control and the duration of diabetes. Diabetic neuropathy is a progressive disorder, with signs and symptoms that parallel the loss of nerve fibers over time. Consequently, assessments of neuropathy in mice are not performed at one time point, but are characterized at multiple time points during a 6 month period of diabetes. The degree of diabetes is evaluated in 2 ways: tail blood glucose measured following a 6 hour fast and glycated hemoglobin levels. The initial degree of neuropathy is screened using the methods discussed below. Detailed measures of neuropathy are employed when the initial screening instruments indicate a profound or unique phenotypic difference. This document contains protocols used by the DiaComp staff to examine and measure diabetic neuropathy at the whole animal, tissue and cellular levels.

### Diabetic Complication:



Neuropathy



## Materials

### Reagent List:

Buffers      Phosphate buffer (PB) pH 7.2, 0.1 M  
                 Phosphate buffered Saline (PBS) pH 7.2, 0.1 M, 150 mM NaCl  
                 Tris, pH 7.2- 7.6, 0.1 M  
                 0.1% TX100 in PB

### Fixatives:

4% paraformaldehyde  
2% paraformaldehyde  
4% paraformaldehyde with 0.5% glutaraldehyde (for TEM)  
1:1 acetone: methanol

**Primary Antibodies:** Working dilutions vary

**Secondary Antibodies:** 1:200 to 1:500

Biotinylated  
Avidin HRP kit  
Fluorochrome labeled

### Counterstains:

Hematoxylin  
Eosin  
Methyl Green  
Bis-benzimide (DAPI)

### Other Chemicals:

3'3-diaminobenzidine  
glucose oxidase  
ammonium chloride  
 $\beta$ -D Glucose  
Ethanol 50, 70, 95, 100%  
Hemo-D  
Normal serum  
Non-fat dry milk  
Bleach  
 $H_2O_2$

### Mounting Media:

Permount  
Prolong  
Gelmount



## **Glassware:**

Square Wheaton dishes

Slide racks and handles

Coplin jars

Beakers

Microfuge tubes

## **Troubleshooting**

## 1 **General Concepts:**

- Keep buffers consistent through fix, rinses, antibody dilutions and chromogen development.
- Antibodies are sticky, always use plastic tips and tubes for storing, diluting and pipeting antibodies.
- Check slides after each major step in the protocol to make sure sections are still stuck to the slide. There is no point in continuing if the sections have all been removed due to over zealous washes.

## 2 **Cryosections:**

1. Remove slides from freezer, air dry. (warm plate)
2. Ring sections with PAP pen, if tissue was not fixed prior to embedding and sectioning, fix sections for 10-30 min, 4°C
  - ♦ 4% paraformaldehyde
  - ♦ 2% paraformaldehyde
  - ♦ 1:1 acetone: methanol, ice cold

## 3 **Paraffin Sections:**

1. Deparaffinize in Hemo-D.
2. Hydrate through ethanol 100, 100, 95, 95, 70, 50, distilled water.
3. Ring sections with PAP pen, carefully, do not let sections dry.

## 4 **Prepare Humid Chamber:**

1. Hydrate in buffer 5 min.
2. Rinse in 0.5% H<sub>2</sub>O<sub>2</sub> in buffer (5 ml 30% in 300 ml buffer), 10 min (for immunofluorescence skip and go to blocking)
3. Rinse 3 X 10 min in buffer.
4. Permeabilize and block 10 minutes:
  - ♦ 0.1% TX100 in buffer
  - ♦ 2% non-fat dry milk
  - ♦ 1% serum (serum from the animal in which the secondary antibody is raised e.g. 1° is rabbit, 2° **goat** anti-rabbit, block with normal goat)
5. Diluted primary antibody in buffer containing 0.1% TX100

6. Remove slides from blocking solution, do not rinse, dab away excess fluid with filter paper, place 1° (20-40  $\mu$ l) in PAP pen ring. Dilutions are antibody specific, check antibody book for suggested dilutions.

7. Always wrap parafilm around the edge of the incubation chamber to seal:

- ◆ Incubation times: 1° 's overnight 22°C, parafilm,  
36-48 hours 4°C, parafilm  
2-4 hours 37°C, lean toward 50-60  $\mu$ l

8. Rinse in buffer 3 X 10 minutes.

## 5 **Immunofluorescence:**

1. Turn off overhead fluorescent lights to decrease quenching.

2. 2° antibody labeled with red or green fluorochrome.

3. Incubate 1 hour 22°C.

4. Rinse 3 X 10 min buffer.

5. Counterstain with Bis-benzimide 1 $\mu$ g/ml buffer, 5 min.

6. Rinse in buffer.

7. Cover with Prolong (see Prolong SOP)

## 6 **Immunoperoxidase:**

1. 2° antibody labeled with biotin.

2. Incubate 1 hour 22°C.

3. Rinse 3 X 10 min buffer.

4. Incubate in Avidin HRP 1 hour 22°C \*\*Prepare this solution 1 hour prior to use, it needs to equilibrate\*\*

5. Rinse 3 X 10 min buffer.

6. DAB

7. Rinse 3 X 10 min in buffer, collect first rinse and add bleach to denature DAB (DO NOT PUT BLEACH ON THE SLIDES)



8. Dehydrate through graded ethanol 50, 70, 95, 95, 100, 100% and 2 changes of Hemo-D Counterstain.
9. Mount coverslips to slides with Permount.