



Jul 06, 2024

Nuclei Isolation

 [eLife](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.14egn6wryl5d/v1>

amanda schneeweis¹

¹Northwestern



amanda schneeweis

Northwestern

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn6wryl5d/v1>

External link: <https://doi.org/10.7554/eLife.101035>

Protocol Citation: amanda schneeweis 2024. Nuclei Isolation. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.14egn6wryl5d/v1>

**Manuscript citation:**

Gaertner Z, Oram C, Schneeweis A, Schonfeld E, Bolduc C, Chen C, Dombeck D, Parisiadou L, Poulin J, Awatramani R Molecular and spatial transcriptomic classification of midbrain dopamine neurons and their alterations in a LRRK2^{G2019S} model of Parkinson's disease. eLife 13(). doi: [10.7554/eLife.101035](https://doi.org/10.7554/eLife.101035)

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 06, 2024

Last Modified: July 06, 2024

Protocol Integer ID: 102927

Keywords: nuclei isolation nuclei isolation for snrna, nuclei isolation, snrna, seq

Funders Acknowledgements:

ASAP

Abstract

nuclei isolation for snRNA-seq

- 1 Prepare buffers and filter sterilize, add RNase inhibitor fresh NP40 Lysis Buffer (NST):
0.1% NP-40 Alternative (or NP-40), 10 millimolar (mM) Tris,
146 millimolar (mM) NaCl, 1 millimolar (mM) CaCl₂, 21mM MgCl₂, 40U/mL of Protector RNase inhibitor (add fresh day of) ST Wash Buffer: (10mM Tris,
146 Mass Percent NaCl, 1 millimolar (mM) CaCl₂, 21mM MgCl₂), 0.01% BSA, 40U/mL of Protector RNase inhibitor (add fresh day of) ST Staining buffer (ST-SB):
2%BSA, 0.02%Tween-20, 10 millimolar (mM) Tris, 146 millimolar (mM) NaCl, 1mM CaCl₂, 21mM MgCl₂, 40U/mL of Protector RNase inhibitor (add fresh day of) Note: Keep tissues/homogenate and buffers on ice throughout the protocol. Pre-cool the centrifuge to 4 °C and keep at 4 °C for all steps.
- 2 Tissue collection a) Sacrifice and rapidly decapitate mice. Using chilled brain matrix (Ted Pella), cut a thick coronal section from 5mm back from start of cortex until start of cerebellum. Store section in ice-cold Hibernate-A media in 10cm plate on ice until all brains have been dissected. b) Use a razor blade to dissect out midbrain and collect tissue, placing directly into dounce homogenizer.
- 3 Tissue lysis and homogenizing a) For each sample to barcode and pool: prepare a separate homogenizer and douncing pestles (loose and tight). Add 1 mL NST buffer to the tube with tissue and transfer to dounce homogenizer and keep on ice. b) with a total volume of 1 mL, dounce 20 times with the loose pestle followed by 20 times with the tight pestle. c) Add 1 mL of ST wash buffer, filter through 30µm filters (Milentyi Biotec 130-041-407) and transfer filtered homogenate to a 15mL tube. d) Rinse the homogenizer with 3x 1 mL of ST wash buffer, filter through 30µm filters and add to the filtered homogenate to add up to a final volume of 5 mL. e) Immediately spin down at 500 g for 00:05:00 at 4 °C to pellet the nuclei in swing bucket rotor f) Remove supernatant g) Resuspend nuclei in 200µl of ST-SB and transfer to lo-bind 1.5 mL tube and wash out the original tube with an additional 1 mL and transfer to the same 1.5ml tube. h) Wash by spinning down for 00:05:00 at 500 g at 4 °C and resuspending in 1.2 mL ST-SB twice, for a total of 3 washes i) resuspend in 1 mL ST-SB

10m



4 Proceed with sorting nuclei for GFP+ nuclei via MACS Tyto