

Jan 08, 2020

## Nuclei Isolation for SNARE-seq2

DOI

[dx.doi.org/10.17504/protocols.io.8tvhwn6](https://dx.doi.org/10.17504/protocols.io.8tvhwn6)

Carter Palmer<sup>1</sup>, Christine Liu<sup>1</sup>, Jerold Chun<sup>2</sup>

<sup>1</sup>University of California, San Diego, and Sanford Burnham Prebys Medical Discovery Institute;

<sup>2</sup>Sanford Burnham Prebys Medical Discovery Institute

BICCN / BICAN



Carter Palmer

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.8tvhwn6>

**Protocol Citation:** Carter Palmer, Christine Liu, Jerold Chun 2020. Nuclei Isolation for SNARE-seq2 . **protocols.io**  
<https://dx.doi.org/10.17504/protocols.io.8tvhwn6>

**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** October 28, 2019

**Last Modified:** January 08, 2020

**Protocol Integer ID:** 29269

**Keywords:** Brain, Nuclei, Isolation, Single-Cell, isolation of nuclei, nucleus chromatin accessibility, snare, frozen brain tissue in preparation, nuclei, seq2, frozen brain tissue, chromatin, nucleus, mrna, tissues from mouse, seq2 this protocol, mrna expression, seq, isolation, sequencing

## Abstract

This protocol is intended to be used for the isolation of nuclei from fresh-frozen brain tissue in preparation for analysis by Single-Nucleus Chromatin Accessibility and mRNA Expression sequencing (SNARE-seq). It has been applied to tissues from mouse, marmoset, and human.

## Guidelines

This protocol is designed specifically for isolating tissue for SNARE-seq. RNA stability is considered at every step. Samples are kept on ice throughout the process, all centrifuges are pre-chilled to 4 °C before use, and RNase Away is used to spray down all surfaces and pipets before use.

## Materials

### MATERIALS

- ☒ Sucrose **Fisher Scientific Catalog #S3-212**
- ☒ Protease Inhibitor Tablets cOmplete Mini EDTA free **Roche Catalog #11836170001**
- ☒ RNase Inhibitor **Takara Bio Inc. Catalog #2313A**
- ☒ Bovine Serum Albumin **Gemini Bio-Products Catalog #700-107P**
- ☒ Magnesium acetate tetrahydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #M5661**
- ☒ Calcium chloride dihydrate **Merck MilliporeSigma (Sigma-Aldrich) Catalog #C5080**
- ☒ Ethylenediaminetetraacetic acid (EDTA) **Merck MilliporeSigma (Sigma-Aldrich) Catalog #EDS**
- ☒ Triton X-100 **Merck MilliporeSigma (Sigma-Aldrich) Catalog #T8787**
- ☒ 16% Formaldehyde (w/v) Methanol-free **Thermo Fisher Scientific Catalog #28906**
- ☒ DAPI **Merck MilliporeSigma (Sigma-Aldrich) Catalog #10236276001**
- ☒ 50um filters **Sysmex Catalog #04-0042-2317**
- ☒ Tissue Homogenizer **Catalog #358005**

## Troubleshooting



## Buffer Preparation

1

### **Nuclei Extraction Buffer (NEB):**

		Final Concent ration		Chemica l to Add	Stock Concent ration
	Sucrose	320 mM		5.48g	Solid
	Magnesium Acetate	3 mM		32.16mg	Solid
	CaCl <sub>2</sub>	5 mM		50µl	5M Stock
	EDTA	100 µM		10µl	0.5M Stock
	Tris-HCl pH 8	10 mM		500µl	1M Stock
	Triton X-100	0.1 %		500µl	10% Stock
	MilliQ Water			Bring to 50ml	
	RNase Inhibitor	80 U/ml		100µl	40,000U /ml
	Protease Inhibitor Tablet			1 Tablet	

Prepare NEB as outlined above. 50ml is enough for 3 samples. Combine reagents 2-8 the night before nuclei isolation and chill at 4 °C. Add RNase inhibitor and Protease Inhibitor tablet the morning of isolation.

2

### **PBSE + Sucrose:**

		Final Concent ration		Chemica l to Add	Stock Concent ration
	Sucrose	320 mM		5.48g	Solid



	EGTA	250 $\mu$ M		25 $\mu$ l	0.5M
	1x PBS pH 7.4			Bring to 50ml	
	RNase Inhibitor	40 U/ml		50 $\mu$ l	40,000U/ml
	Protease Inhibitor Tablet			1 Tablet	

Prepare PBSE + Sucrose as outlined above. 50ml is enough for 3 samples. Combine reagents 2-4 the night before nuclei isolation and chill at 4 °C. Add RNase inhibitor and Protease inhibitor tablet the morning of isolation.

### 3 **PBSE + BSA:**





		Final Concentration		Chemical to Add	Stock Concentration
	BSA	1%		500mg	Solid
	EGTA	250 $\mu$ M		25 $\mu$ l	0.5M
	1x PBS pH 7.4			Bring to 50ml	
	RNase Inhibitor	40 U/ml		50 $\mu$ l	40,000U/ml
	Protease Inhibitor Tablet			1 Tablet	

Prepare PBSE + BSA as outlined above. 50ml is enough for 3 samples. Combine reagents 2-4 the night before nuclei isolation and chill at 4 °C. Add RNase inhibitor and Protease inhibitor tablet the morning of isolation.

## Tissue Dissociation

- 4 Remove tissue from -80°C storage and place on ice.
- 5 Image the tissue rapidly in front of a ruler







- 6 Add 1ml of ice cold NEB to the tissue and incubate on ice for  00:15:00
- 7 During the incubation, wash the homogenizer with MilliQ water, 10% bleach, 70% EtOH, and MilliQ water again
- 8 Rinse homogenizer with 1ml of ice cold NEB
- 9 Once the incubation is complete, add 1 ml of fresh NEB to the homogenizer and transfer the tissue and the 1ml of NEB it is in to the homogenizer
- 10 Homogenize the tissue using ~20 compressions with the pestle, or until the tissue is entirely dissociated, this step can be variable
- 11 Pass nuclei suspension through 50 micron filter into a 15 ml conical tube, wash filter with another 4ml of NEB
- 12 Wait for  00:05:00
- 13 Centrifuge the samples  820 x g , 5 min
- 14 Carefully aspirate the supernatant and slowly resuspend the pellet in 1ml of NEB
- 15 Gently add another 9 ml of NEB to the sample
- 16 Centrifuge the samples  820 x g , 5 min


## Nuclei Fixation

- 17 Aspirate the supernatant and slowly resuspend the samples in 1ml of PBSE + Sucrose
- 18 Gently add another 5 ml of PBSE + Sucrose



- 19 Centrifuge the samples  820 x g , 5 min
- 20 Aspirate all but approximately  100  $\mu$ L of the PBSE+Sucrose, and gently resuspend the nuclei pellet in the remaining volume
- 21 While gently vortexing, add 5ml of 0.5% Formaldehyde diluted in 1X PBS
- 22 Fix for 10 min on ice
- 23 Centrifuge the samples  820 x g , 5 min
- 24 Aspirate the supernatant and slowly resuspend the samples in 1ml of PBSE + Sucrose
- 25 Centrifuge the samples  820 x g , 5 min

## DAPI Staining and Sorting

- 26 Aspirate the supernatant and slowly resuspend the samples in 1ml of PBSE + BSA
- 27 Gently add another 10ml of PBSE + BSA
- 28 Centrifuge the samples  820 x g , 5 min
- 29 Make 1ml/sample of PBSE + BSA + DAPI by adding DAPI at a final concentration of 1.25  $\mu$ g/ml to PBSE + BSA. This is a 1:4000 dilution from a DAPI stock at 5mg/ml
- 30 Aspirate the supernatant and slowly resuspend the samples in 1ml of PBSE + BSA + DAPI



- 31 Pass nuclei suspension through 50 micron filter into a FACS tube, gently tap on table to get all nuclei through the filter
- 32 Use the DAPI singlet peak to sort singlet nuclei events into a 2ml low binding eppendorf tube.
- 33 Store all samples on ice until SNARE-seq processing begins. Do not store for longer 4 hours before processing.