

Jul 02, 2025

Nuclei isolation from brain tissue for single-cell ATAC

DOI

<https://dx.doi.org/10.17504/protocols.io.14egnr7ql5d/v1>

Koen Theunis^{1,2,3,4}, Suresh Poovathingal^{5,6}

¹Laboratory of Computational Biology, VIB Center for AI & Computational Biology, Leuven, Belgium;

²VIB-KU Leuven Center for Brain & Disease Research, Leuven, Belgium;

³Department of Human Genetics, KU Leuven, Leuven, Belgium;

⁴Aligning Science Across Parkinson's (ASAP) Collaborative Research Network, Chevy Chase, MD 20815, United States;

⁵VIB KU Leuven Center for Brain & Disease Research;

⁶VIB Center for AI & Computational Biology (VIB.AI), Leuven, Belgium

Aertslab



Koen Theunis

Laboratory of Computational Biology, VIB Center for AI & Com...

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egnr7ql5d/v1>



Protocol Citation: Koen Theunis, Suresh Poovathingal 2025. Nuclei isolation from brain tissue for single-cell ATAC. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.14egnr7q15d/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 30, 2025

Last Modified: July 02, 2025

Protocol Integer ID: 221299

Keywords: ASAPCRN, nuclei isolation from brain tissue, cell atac this protocol, nuclei isolation, cell atac assay, nuclei from brain tissue, isolating nuclei, cell atac, brain tissue, atac, single cell, nuclei, isolation

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-025179

Abstract

This protocol is for isolating nuclei from brain tissue and prepare them for single-cell ATAC assays like the 10X Genomics Chromium Next GEM Single Cell ATAC Kit, HyDrop-ATAC, etc.

Attachments



Nuclei_Isolation_ATA...

18KB

Materials

	A	B	C	D	E	F
	2X HB- (Homogenization Buffer)	Stock conc		Final conc		V (µl)
	Sucrose	2000	mM	640	mM	2245, 6
	NaCl	5000	mM	20	mM	28,1
	MgCl ₂	1000	mM	6	mM	42,1
	Tris pH7,5	1000	mM	20	mM	140,4
	EDTA	500	mM	0,2	mM	2,8
	BSA	7,5	%	0,08	%	74,9
	Tween20	10	%	0,02	%	14,0
	Proteinase inhibitor	50	X	2	X	280,7
	DTT	1000	mM	2	mM	14,0
	Water					4175, 0
	Total Volume					7017, 6

	A	B	C	D	E	F
	1X HB+ (Homogenization Buffer)	Stock conc		Final conc		V (µl)
	2X HB-	2	X	1	X	2200,0
	Igepal	10	%	0,05	%	22,0
	Digitonin	2	%	0,0025	%	5,5
	Water					2178,0
	Total Volume					4400,0



A	B	C	D	E	F
1X HB- (Homogenization Buffer)	Stock conc		Final conc		V (μl)
2X HB- (Homogenization Buffer)	2	X	1	X	3648, 0
Water					3648, 0
Total Volume					7296, 0

A	B	C	D	E	F
GM (Gradient Medium)	Stock conc		Final conc		V(μl)
CaCl ₂	1000	mM	5	mM	12,5
Optiprep	60	%	50	%	2080,0
MgCl ₂	1000	mM	3	mM	7,5
Tris pH 7.5	1000	mM	10	mM	25,0
BSA	7,5	%	0,04	%	13,3
Proteinase inhibitor	50	X	0,2	X	10,0
DTT	1000	mM	1	mM	2,5
Water					345,3
Total Volume					2496,0

A	B	C	D	E	F
ODM (Optiprep Diluent Medium)	Stock conc		Final conc		V(μl)
KCl	3000	mM	150	mM	114,6



	A	B	C	D	E	F
	MgCl ₂	1000	mM	30	mM	68,7
	Tris pH 8	1000	mM	60	mM	137,5
	Sucrose	2000	mM	250	mM	286,4
	Water					1684,3
	Total Volume					2291,5

	A	B	C	D	E	F
	29% Cushion	Stock conc		Final conc		V(μl)
	Optiprep	60	%	29	%	1786,4
	ODM					1909,6
	Total Volume					3696,0

	A	B	C	D	E	F
	Resuspension Buffer	Stock conc		Final conc		V(μl)
	20X Nuclei buffer	20	X	1	X	22
	BSA	7,5	%	1	%	58,7
	Water					359,3
	Total Volume					440

A. Homogenization

- 1 Cut brain piece on dry ice, then transfer immediately in homogenizer containing 1 mL of ice-cold HB+. If tissue is still frozen, leave it in HB for 3 min before starting homogenization to prevent the nuclei from breaking.
- 2 Homogenize tissue with 10 manual gentle strokes (pestle A) + 10 manual gentle strokes (pestle B).
- 3 Put homogenate through 70 μ m cell strainer over a 50 mL falcon (or immediately a 2 mL protein lo-bind tube).
- 4 Wash homogenizer & strainer with 1 mL HB- to a final volume of 2 mL.
- 5 Transfer the contents from the 50 mL falcon to a 2 mL protein lo-bind tube.
- 6 Spin nuclei @500g for 5 min in a swinging bucket and remove supernatant.
- 7 Resuspend pellet with HB- to get a total volume of 520 μ L and transfer to a clean DNA lo-bind tube.
- 8 Add 520 μ L of Gradient medium (GM) to sample (V_f = 1040 μ L).

B. Isolation by centrifugation

- 9 Layer 770 μ L of 29% Cushion in a 2 mL DNA LoBind tube.
- 10 Layer the sample on top of cushion using a P1000, without disturbing the cushion.
- 11 Centrifuge at least 3,000 rcf (you can go as high as 9,000) in a swinging bucket, at 4°C for 20 minutes with brake off.
- 12 Remove supernatant, remove the lower supernatant with P200, leaving about max 30 μ L.



C. Resuspension and counting

- 13 Disrupt nuclei on bottom by pipetting with P200. And bring X μ L to a fresh 1.5 mL LoBind tube.
- 14 Add X μ L of 1X Resuspension buffer to tube to retain extra nuclei, and add them to the 1.5 mL tube from C.1.
- 15 Spin 5 min @ 350-450xg followed by resuspension in 1X Diluted nuclei buffer. Optional filter and count.