

Sep 08, 2017

Oil Red O Staining

7

Forked from Oil Red O Staining Drosophila Larval and Prepupal Tissues

PLOS One

DOI

dx.doi.org/10.17504/protocols.io.iemcbc6

Yuejun Wang¹

¹Department of Prosthodontics, Peking University School and Hospital of Stomatology



Yuejun Wang

OPEN ACCESS



DOI: dx.doi.org/10.17504/protocols.io.iemcbc6

External link: https://doi.org/10.1371/journal.pone.0184182

Protocol Citation: Yuejun Wang 2017. Oil Red O Staining . protocols.io https://dx.doi.org/10.17504/protocols.io.iemcbc6

Manuscript citation:

Wang Y, Liu Y, Fan Z, Liu D, Wang F, Zhou Y (2017) IGFBP2 enhances adipogenic differentiation potentials of mesenchymal stem cells from Wharton's jelly of the umbilical cord *via* JNK and Akt signaling pathways. PLoS ONE 12(8): e0184182. doi: 10.1371/journal.pone.0184182

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

Created: June 12, 2017

Last Modified: March 17, 2018

Protocol Integer ID: 6317

Keywords: stempro adipogenesis differentiation kit, adipogenic differentiation, adipose, stained cell, isopropanol, total protein concentration, protein



Abstract

Adipogenic differentiation was induced by using the StemPro adipogenesis differentiation kit (Invitrogen, Carlsbad, CA, USA). WJCMSCs were grown in the adipose-inducing medium for 3 weeks. For Oil Red O staining, after induction, cells were fixed with 10% formalin for at least 1 h at room temperature. Next, cells were stained with the 60% Oil Red O in isopropanol as working solution for 10 min. The proportion of Oil Red O-positive cells was determined by counting stained cells under a light microscope. The final OD value in each group was normalized with the total protein concentrations prepared from a duplicate plate.

- 1 Cells were grown in the adipose-inducing medium for 2-3 weeks.
- 2 After induction, cells were fixed with 10% formalin for at least 1 h at room temperature.
- 3 Cells were stained with the 60% Oil Red O in isopropanol as working solution for 10 min.
- 4 The proportion of Oil Red O-positive cells was determined by counting stained cells under a light microscope.
- 5 The final OD value in each group was normalized with the total protein concentrations prepared from a duplicate plate.