

Oct 19, 2024

Olfactory Organoids

DOI

<https://dx.doi.org/10.17504/protocols.io.14egn33zzl5d/v1>

Alp Ozgun¹, Priya Suman¹, Josée oulombe¹, Julianna Tomlinson¹, Earl G. Brown¹, John M. Woulfe¹,
Michael G. Schlossmacher¹

¹Ottawa Hospital Research Institute



Alp Ozgun

Ottawa Hospital Research Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.14egn33zzl5d/v1>

Protocol Citation: Alp Ozgun, Priya Suman, Josée oulombe, Julianna Tomlinson, Earl G. Brown, John M. Woulfe, Michael G. Schlossmacher 2024. Olfactory Organoids. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.14egn33zzl5d/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working



Created: August 14, 2023

Last Modified: October 19, 2024

Protocol Integer ID: 86475

Keywords: ASAPCRN, olfactory organoids from krt5, embedded olfactory organoid, olfactory organoids this protocol enable, olfactory organoid, hollow spheres of olfactory epithelium, olfactory horizontal basal cell, olfactory epithelium, hydrogel matrix dome, organoid, obtaining 3d matrix, cell, containing hollow sphere, 3d matrix

Funders Acknowledgements:

Aligning Science Across Parkinson's

Grant ID: ASAP-020625

Abstract

This protocol enables obtaining 3D matrix-embedded olfactory organoids from KRT5+ olfactory horizontal basal cells. Resulting samples consist of hydrogel matrix domes with a volume of  40 µL containing hollow spheres of olfactory epithelium. The amounts were given per sample and should be calculated according to how many samples will be made.



Materials

Organoid initiation medium:

- 1:1 mixture of Advanced DMEM and Neurabasal A media
- 1x B27 supplement
- 1x N2 supplement
- 1% Glutamax
- 1% penicillin/streptomycin
- 3 μ M CHIR99021
- 500 ng/mL R-Spondin 1
- 1.5 μ M DMH-1
- 1.5 μ M A83-01

Maintenance medium:

- 1:1 mixture of DMEM/F12 and Neurabasal A media
- 1x B27 supplement
- 1x N2 supplement
- 1% Glutamax
- 1% penicillin/streptomycin
- 3 μ M CHIR99021
- 1 μ M DMH-1
- 1 μ M A83-01

HBC medium:

- Pneumacult-Ex medium
- 1x Glutamax
- 1x B27 supplement without vitamin A
- 1x N2 supplement
- 1% pen/strep
- 10ng/mL TGF alpha
- 1 μ M A-83-01
- 1 μ M DMH-1



- 1 Thaw reduced growth factor matrix precursor aliquots at  4 °C overnight.  20 µL matrix precursor is required per sample. Geltrex and Matrigel were tested and shown to work equally well with this protocol. Keep these products at  4 °C or on ice throughout the protocol.
- 2 Place 24-well plates into tissue culture incubator at  37 °C to equilibrate the temperature of the plastic. This step is important to prevent sinking of the cells to surface inside the matrix precursor.
- 3 Lift olfactory basal cells from the culture surface by rinsing with dPBS and adding cold accutase. Incubate at  37 °C for up to 15 minutes.
- 4 Add HBC medium equal to the volume of accutase and transfer the cells into a falcon tube. Centrifuge at  350 x g for 4 minutes and discard the supernatant. Resuspend cells in  2 mL HBC medium and perform cell count.
- 5 Transfer 4×10^4 cells per sample into a new falcon tube and centrifuge at  350 rpm . Discard supernatant. Add  15 µL organoid initiation medium per sample and resuspend the cells. Perform the following steps as quickly as possible.
- 6 Add  30 µL matrix precursor per sample to the cell suspension and briefly triturate to homogenize. Retrieve 24-well plates from the incubator and pipette  40 µL of the cell suspension into the center of each well, forming a liquid dome. Immediately return the well plates to the incubator. Aim to finish pipetting and return the plates to the incubator within 3-4 minutes after adding the matrix precursor to the cells.
- 7 Allow the matrix solutions to set in the incubator for 30 minutes. Confirm that the cells did not settle to the bottom using an inverted microscope. Gently fill the wells with  500 µL organoid initiation medium by pointing the pipette tip to the well wall.
- 8 Return the well-plates to the incubator. This day is denoted as 0 days-in-vitro (0 DIV).



- 9 At the end of 4 DIV, carefully aspirate the organoid initiation medium and add  500 μ L maintenance medium. At this stage, cell clumps and hollowing spheres should be observed under the microscope.

- 10 Over the following days, spheres should start to hollow out. At the end of Day 7, supplement the maintenance medium with  3 micromolar (μ M) retinoic acid. The sphere walls will then start to thicken and differentiate into pseudostratified architecture of the olfactory epithelium. The samples are ready for further testing by 20 DIV.