

Aug 21, 2025

Open field test in rats

 Forked from [Open field test in rats](#)

 In 1 collection

DOI

<https://dx.doi.org/10.17504/protocols.io.rm7vzqw78vx1/v1>

Eduard Bentea¹, María Sanchiz Calvo¹, Veerle Baekelandt¹

¹Katholieke Universiteit Leuven



María Sanchiz Calvo

KU Leuven

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.rm7vzqw78vx1/v1>

Protocol Citation: Eduard Bentea, María Sanchiz Calvo, Veerle Baekelandt 2025. Open field test in rats. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.rm7vzqw78vx1/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: June 24, 2025

Last Modified: August 21, 2025

Protocol Integer ID: 220874

Keywords: ASAPCRN, motor function, behavioral test, rats, open field test in rat, open field test in rats protocol, mobility during the open field test, open field test, nigrostriatal dopaminergic neurotransmission show, decreased nigrostriatal dopaminergic neurotransmission show, parkinson, spontaneous locomotion, min free exploration of an open field arena, open field arena, rats protocol, mobility, test

Funders Acknowledgements:

ASAP (Aligning Science Across Parkinson's)

Abstract

Protocol for performing the open field test in rats. The open field test measures spontaneous locomotion during 5-min free exploration of an open field arena (1m x 1m), and can be used to measure hypolocomotion in animal models of Parkinson's disease. Rats with decreased nigrostriatal dopaminergic neurotransmission show decreased mobility during the open field test.



Test

1h 5m

- 1 Bring cages to the behavioral room for at least  01:00:00 prior to the test for habituation 
- 2 Place rat individually in a clean open field arena (1m x 1m)
- 3 Video tape from above for  00:05:00 and record parameters of spontaneous activity (distance traveled, movement time, immobility) using an activity tracking software. 
- 4 Clean arena with mild soap solution