



Mar 30, 2024

Optical Density Analysis for DAB Staining of Mouse Brain Sections

 [eLife](#)

DOI

<https://dx.doi.org/10.17504/protocols.io.81wgbxo2nlpk/v1>

Katerina Rademacher¹, Ken Nakamura¹

¹Gladstone Institute of Neurological Disease



Haru Yamamoto

UCSF

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.81wgbxo2nlpk/v1>

External link: <https://doi.org/10.7554/eLife.98775>

Protocol Citation: Katerina Rademacher, Ken Nakamura 2024. Optical Density Analysis for DAB Staining of Mouse Brain Sections. [protocols.io](#) <https://dx.doi.org/10.17504/protocols.io.81wgbxo2nlpk/v1>

Manuscript citation:

Rademacher K, Doric Z, Haddad D, Mamaligas A, Liao S, Creed R, Kano K, Chatterton Z, Fu Y, Garcia JH, Vance VM, Sei YJ, Kreitzer A, Halliday G, Nelson AB, Margolis E, Nakamura K Chronic hyperactivation of midbrain dopamine neurons causes preferential dopamine neuron degeneration. eLife 13(). doi: [10.7554/eLife.98775](https://doi.org/10.7554/eLife.98775)



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 07, 2024

Last Modified: May 31, 2024

Protocol Integer ID: 93033

Keywords: ASAPCRN, optical density analysis for dab staining, staining of mouse brain section, optical density analysis, steps for optical density analysis, mouse brain section, dab staining, mouse brain sections this protocol, dab

Funders Acknowledgements:

ASAP

Grant ID: 020529

Abstract

This protocol describes steps for optical density analysis of immunohistochemical 3,3'-Diaminobenzidine (DAB) staining of mouse brain sections using ImageJ.

Materials

- FIJI/ImageJ Software (RRID:SCR_002285)



- 1 Open ImageJ, then calibrate for analysis.
 - 1.1 See <https://imagej.net/ij/docs/examples/calibration/> for calibration setup steps and files.
 - 1.2 Select Analyze > Calibrate, then select the following:
 - i. Rodbard (NIH Image)
 - ii. Unit: O.D.
 - iii. Open the calibration.txt file
 - iv. Uncheck 'show plot'
 - v. Check 'Global calibration'
- 2 Open an image and outline the regions of interest, including a background area for subtraction.
- 3 Select 'multi measure' and copy the results to an Excel sheet for analysis. Repeat for all images.
- 4 Take the 'mean OD' value for each region and subtract the background area 'mean OD' value. Values can be plotted directly or normalized to control groups.