


Feb 05, 2024

🌐 Optimising sample multiplexing oligos by flow cytometry

DOI

<https://dx.doi.org/10.17504/protocols.io.81wgbyjpovpk/v1>

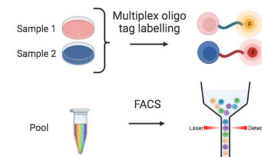

Daniel V Brown¹

¹Walter and Eliza Hall Institute



Daniel V Brown

Walter and Eliza Hall Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.81wgbyjpovpk/v1>

Protocol Citation: Daniel V Brown 2024. Optimising sample multiplexing oligos by flow cytometry. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.81wgbyjpovpk/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 14, 2023

Last Modified: February 05, 2024

Protocol Integer ID: 81857

Keywords: scRNA-Seq, facs, sample multiplexing, hashtag, sample multiplexing oligos by flow cytometry optimisation, sample multiplexing oligos by scrna, sample multiplexing oligo, optimising sample multiplexing oligo, flow cytometry read, flow cytometry optimisation, fluorescent detection, out with fluorescent detection, labelled oligo, scrna, signal swapping

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to [protocols.io](https://www.protocols.io) is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with [protocols.io](https://www.protocols.io), can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Optimisation of sample multiplexing oligos by scRNA-Seq is costly and time consuming. A cheaper and faster method is to use a flow cytometry read-out with fluorescent detection oligonucleotides.

This method can also be used to mix samples with different fluorescently labelled oligos and investigate signal swapping.

Image Attribution

Made with Biorender.com



Guidelines

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Materials

- Phosphate Buffered Saline, without magnesium and calcium
- Sample multiplexing oligos (MULTI-Seq, CellPlex or Hashtag Antibody)
- 10ug/mL DAPI
- 30% BSA stock solution
- Fluorescent detection oligonucleotides

A	B	C
Name	Sequence	Modification
A647_FB2_detect	/5Alex647N/CCTTAGCCGC TAATAGGTGAGC	5' Alexa 647 modification
A594_FB2_detect	/5Alex594N/TTGCTAGGAC CGGCCTTAAAGC	5' Alexa 594 modification
A594_FB1_detect	/5Alex594N/TTGCTAGGAC CGGCCTTAAAGC	5' Alexa 594 modification
A647_Total- SeqC_detect	/5Alex647N/CTGTCTCTTA TACACATCTCCG	5' Alexa 647 modification
AF647_oligo_dT_detect	/5Alex647N/TTTTTTTTTTTT TTTTTTTTTTTTTTTTTTTT	5' Alexa 647 modification

Fluorescent detection oligonucleotides. Order with HPLC modification

MULTI-Seq barcoding oligos. I substituted the poly-A tail for 10x Genomics feature barcode 2 sequence.

A	B	C
Name	Sequence	Barco de
multiSeq_FB2_B C2	CCTTGGCACCCGAGAATT CCACCACAATGGCTCACC TATTAGCGGCTAAGG	CCAC AATG
multiSeq_FB2_B C3	CCTTGGCACCCGAGAATT CCATGAGACCTGCTCACCT ATTAGCGGCTAAGG	TGAGA CCT
multiSeq_FB2_B C4	CCTTGGCACCCGAGAATT CCAGCACACGCGCTCACC TATTAGCGGCTAAGG	GCAC ACGC
multiSeq_FB2_B C5	CCTTGGCACCCGAGAATT CCAAGAGAGAGGCTCACC TATTAGCGGCTAAGG	AGAG AGAG

	A	B	C
	multiSeq_FB2_B C6	CCTTGGCACCCGAGAATT CCATCACAGCAGCTCACC TATTAGCGGCTAAGG	TCACA GCA
	multiSeq_FB2_B C7	CCTTGGCACCCGAGAATT CCAGAAAAGGGGCTCACC TATTAGCGGCTAAGG	GAAA AGGG
	multiSeq_FB2_B C8	CCTTGGCACCCGAGAATT CCACGAGATTTCGCTCACC TATTAGCGGCTAAGG	CGAG ATTC
	multiSeq_FB2_B C9	CCTTGGCACCCGAGAATT CCAGTAGCACTGCTCACCT ATTAGCGGCTAAGG	GTAGC ACT
	multiSeq_FB2_B C10	CCTTGGCACCCGAGAATT CCACGACCAGCGCTCACC TATTAGCGGCTAAGG	CGAC CAGC
	multiSeq_FB2_B C11	CCTTGGCACCCGAGAATT CCATTAGCCAGGCTCACCT ATTAGCGGCTAAGG	TTAGC CAG
	multiSeq_FB2_B C12	CCTTGGCACCCGAGAATT CCAGGACCCCAGCTCACC TATTAGCGGCTAAGG	GGAC CCCA
	multiSeq_FB2_B C13	CCTTGGCACCCGAGAATT CCACCAACCGGGCTCACC TATTAGCGGCTAAGG	CCAA CCGG

Safety warnings

 Please follow all Manufacturer safety warnings and recommendations.



Prepare multiplexing reagent

1 For CellPlex and Hashtag antibody the reagent comes ready to use.

1.1 Prepare a dilution series for titration if desired

MULTI-Seq oligo preparation

2 Mix anchor and barcode strands in 1:1 molar ratio in PBS (without FBS or BSA at 2 μ M concentration (10X stock).

2.1 This is 6uL 50uM anchor LMO, and 15uL 10uM barcode oligo to 129uL plain PBS.

2.2 Make one unique barcode solution per sample.

2.3 Total 25 μ L per sample.

3 Make a 10X solution of the Co-Anchor in PBS.

3.1 Add 3uL 50uM co-anchor to 141uL plain PBS.

3.2 Add 6uL 100uM fluorescent detection oligo. For example:
Alexa 647 feature barcode 2 detection oligo.

3.3 It is highly recommended to label in at least duplicate. On these occasions I label the other replicate with a Alexa 594 detection oligo and mix immediately before FACS analysis.

Sample preparation

4 I use suspension cell lines to titrate cell multiplexing oligos so the sample preparation is easy.



- 4.1 Prepare a single-cell suspension of the sample to be tested.
- 4.2 Wash cells once in plain PBS without additives. I centrifuge suspension cell lines at 400xg.
Resuspend in PBS.
- 4.3 Count cells and transfer 100k to 1M cells (preferably 500k) into a 1.5mL tube for labelling

Labelling samples with multiplexing oligos

- 5 This is largely based on original protocols but with the addition of a fluorescent secondary oligo.
- 5.1 Resuspend cells in 180uL of plain PBS.
- 5.2 Add 20 μ L 10X Anchor:Barcode solution and pipette gently to mix 10 – 15 times.
- 5.3 Incubate on ice for 5 minutes.
- 5.4 Add 20 μ L Co-Anchor solution and pipette gently to mix.
- 5.5 Incubate 5 minutes longer on ice.
- 5.6 Add 1.5mL of 1% BSA in PBS (ice cold) to quench.
- 5.7 Add 1.5mL of 1% BSA in PBS (ice cold) to quench.
- 5.8 Add 1mL of 1% BSA in PBS (ice cold) to quench.
- 5.9 Centrifuge cells at 4°C 400xg 5min.



- 5.10 Resuspend cells in the remaining 100uL of supernatant then transfer to a new 1.5mL tube.
- 5.11 Add 1.9mL 1% BSA in PBS and spin 5 minutes 400g at 4°C.
- 5.12 Repeat wash 2 more times for a total of 3 washes.
- 5.13 Resuspend cells in 500 µL of PBS + 1% BSA and transfer to 5mL polystyrene FACS tube

Flow cytometry

- 6 Use a FACS analyser to compare the signal of labelled samples to a control sample where only the fluorescent detection oligo has been added.
If you have replicate labelled samples with different fluorescent detection oligos, you may combine them at this step.
- 6.1 Add DAPI to a final concentration of 0.1ug/mL (1/100 stock tube).
- 6.2 Gate for debris, single cells and viable cells.
- 6.3 Acquire at least 10,000 viable cells per test.
- 6.4 Analyse with relevant software.



Protocol references

Marlon Stoeckius, Shiwei Zheng, Brian Houck-Loomis, Stephanie Hao, Bertrand Z Yeung, William M Mauck, Peter Smibert, and Rahul Satija. Cell hashing with barcoded antibodies enables multiplexing and doublet detection for single cell genomics. *Genome biology*, 19(1):1–12, 2018.

Christopher S McGinnis, David M Patterson, Juliane Winkler, Daniel N Conrad, Marco Y Hein, Vasudha Srivastava, Jennifer L Hu, Lyndsay M Murrow, Jonathan S Weissman, Zena Werb, et al. Multi-seq: sample multiplexing for single-cell rna sequencing using lipid-tagged indices. *Nature methods*, 16(7):619–626, 2019