

Oct 28, 2024

p-S65-Ub sandwich ELISA

DOI

<https://dx.doi.org/10.17504/protocols.io.n2bvjng3bgk5/v1>

Jens Watzlawik¹, Fabienne Fiesel¹, Wolfdieter Springer¹

¹Mayo Clinic



Francesca Tonelli

MRC-PPU at The University of Dundee

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvjng3bgk5/v1>

Protocol Citation: Jens Watzlawik, Fabienne Fiesel, Wolfdieter Springer 2024. p-S65-Ub sandwich ELISA. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.n2bvjng3bgk5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 17, 2024



Last Modified: October 28, 2024

Protocol Integer ID: 104415

Keywords: ASAPCRN, ELISA, PINK1, Ubiquitin, Serine-65, ub sandwich elisa, ubiquitin, linked immunosorbent assay, immunosorbent assay, levels of phosphorylation, phosphorylation, enzyme, assay, vivo by enzyme, elisa

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to **protocols.io** is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with **protocols.io**, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

Here we describe an assay to measure the levels of phosphorylation of Ubiquitin at Serine-65 in vivo by enzyme linked immunosorbent assay (ELISA). This assay was previously described by Watzlawik and colleagues [<https://doi.org/10.1080/15548627.2020.1834712>].

Materials

-  Phospho-Ubiquitin (Ser65) (E2J6T) Rabbit mAb **Cell Signaling Technology Catalog #62802**
-  Ubiquitin Monoclonal Antibody (eBioP4D1 (P4D1)), eBioscience™ **Thermo Fisher Scientific Catalog #14-6078-37**
- 200 mM sodium carbonate buffer pH 9.7
-  MULTI-ARRAY 96 Plate Pack, SECTOR Plate **MESO SCALE DIAGNOSTICS, LLC. Catalog #L15XA-3**
- **ELISA washing buffer**

	A	B
	Tris, pH 7.4	150 mM
	NaCl	150 mM
	Tween-20	0.1% [v:v]

- **ELISA blocking buffer**

	A	B
	Tris, pH 7.4	150 mM
	NaCl	150 mM
	Tween-20	0.1% [v:v]
	BSA	1% [w:v]

-  Anti Mouse Antibody Goat SULFO-TAG Labeled **MESO SCALE DIAGNOSTICS, LLC. Catalog #R32AC-1**
-  MSD GOLD Read Buffer A **MESO SCALE DIAGNOSTICS, LLC. Catalog #R92TG**
- microplate mixer (USA Scientific, 8182–2019)



Protocols:

3h

- 1 Prepare the sodium carbonate buffer, ELISA washing buffer and ELISA blocking buffer.
- 2 Use rabbit monoclonal p-S65-Ub antibody as capture antibody and used in a final concentration of 1 µg/ml in 200 millimolar (mM) sodium carbonate buffer 9.7 and coated Overnight at 4 °C with 30 µL per well in a 96-well MSD plate (MSD, MULTI-ARRAY 96-well Plate; L15XA-3).
- 3 Wash the next morning MSD plates twice with 0.22-micron filtered ELISA washing buffer using a 1200 µL multi-(12)-channel pipettor with 300 µL per well. Perform the washing by plate inversion and gentle tapping on paper towels (not by pipette aspiration) and with no incubation time between washes.



- ELISA washing buffer

	A	B
	Tris, pH 7.4	50 mM
	NaCl	150 mM
	Tween-20	0.1% [v:v]

- 4 Block the plates with 0.22-micron filtered ELISA blocking buffer for 01:00:00 at 22 °C without shaking.

1h



- ELISA blocking buffer

	A	B
	Tris, pH 7.4	50 mM
	NaCl	150 mM
	Tween-20	0.1% [v:v]
	BSA	1% [w:v]



- 5 Run all samples in duplicates and diluted in blocking buffer. Load  30 μg of total protein per well for all mouse tissues in a total volume of  30 μL per well. Adjust the detergent volumes across all samples.
- 6 Incubate the antigens for  02:00:00 at  22 $^{\circ}\text{C}$ on a microplate mixer (USA Scientific, 8182–2019) at  600 rpm and perform the three washing steps as described before.  2h
- 7 Subsequently add mouse total Ub antibody (clone P4D1; Thermo Fisher #14-6078-37) as detecting antibody in a final concentration of  5 $\mu\text{g}/\text{ml}$ in blocking buffer in  30 μL total volume per well. Incubate detecting antibody for  02:00:00 at  22 $^{\circ}\text{C}$ on a microplate mixer at  600 rpm.  2h
- 8 After three washing steps, add the SULFO-TAG labeled goat anti-mouse antibody (MSD, R32AC-1) in blocking buffer in a final concentration of  1 $\mu\text{g}/\text{ml}$ using  50 μL per well and incubate for  01:00:00 at  22 $^{\circ}\text{C}$ on a microplate mixer at  500 rpm.  1h
- 9 After another three washing steps,  150 μL MSD GOLD Read Buffer (MSD, R92TG-2) were finally added to each well and the plate being read on a MESO QuickPlex SQ 120 reader.
- 10 Perform the data analysis in GraphPad prism 9 or similar software.