



Oct 06, 2015

pGGDestTol2LC-sgRNA ligation mixture

DOI

<https://dx.doi.org/10.17504/protocols.io.drw57d>

Linlin Yin¹, Lisette A. Maddison¹, Mingyu Li¹, Nergis Kara¹, Matthew C. LaFave¹, Gaurav K. Varshney¹, Shawn M. Burgess¹, James G. Patton¹, and Wenbiao Chen¹

¹Genetics

Genetics



Tracey Depellegrin

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.drw57d>

Protocol Citation: Linlin Yin, Lisette A. Maddison, Mingyu Li, Nergis Kara, Matthew C. LaFave, Gaurav K. Varshney, Shawn M. Burgess, James G. Patton, and Wenbiao Chen 2015. pGGDestTol2LC-sgRNA ligation mixture. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.drw57d>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

Created: September 08, 2015

Last Modified: November 09, 2017

Protocol Integer ID: 1558

Keywords: sgrna ligation mixture, pggdesttol2lc



- 1 2µl 10x CutSmart buffer
- 2 2 µl T4 DNA ligase buffer
- 3 100ng of each pU6x-sgRNA plasmid
- 4 50ng empty pGGDestTol2LC-sgRNA vector of choice (see plasmid list and diagrams in Guidelines),
- 5 1µl BsaI
- 6 1µl T4 DNA ligase
- 7 adjust volume to 20µl with H₂O