

Mar 29, 2023

Version 3

PAS Staining of Fresh Frozen or Paraffin Embedded Human Kidney Tissue V.3

DOI

<https://dx.doi.org/10.17504/protocols.io.x54v98nkzl3e/v3>

Elizabeth Neumann¹, Jamie Allen¹, Angela Kruse¹, Jennifer Harvey¹, Maya Brewer², Carrie Romer¹, Mark De Caestecker², Jeff Spraggins¹

¹Vanderbilt University; ²Division of Nephrology, Vanderbilt University Medical Center

VU Biomolecular Multim...

Human BioMolecular Atl...



Jamie Allen

Vanderbilt University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.x54v98nkzl3e/v3>

Protocol Citation: Elizabeth Neumann, Jamie Allen, Angela Kruse, Jennifer Harvey, Maya Brewer, Carrie Romer, Mark De Caestecker, Jeff Spraggins 2023. PAS Staining of Fresh Frozen or Paraffin Embedded Human Kidney Tissue. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.x54v98nkzl3e/v3> Version created by **Angela R.S. Kruse**

**Manuscript citation:**

References: 1. Luna, Lee (ed.). Manual of histological staining methods of the Armed Forces Institute of Pathology. 2. Dubowitz, B. Muscle Biopsy. A practical approach, 2nd edition, Bailliere, Tindall, London, 1985. 3. Dr. Fogo Clinical lab PAS protocol.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: February 23, 2023

Last Modified: October 18, 2023

Protocol Integer ID: 77523

Keywords: paraffin embedded human kidney tissue scope, kidney pathology, pas staining of fresh frozen, glycoprotein, such as glycoprotein, stain, tissue, mucins in tissue, staining, pas staining, polysaccharides such as glycogen, glycogen, pas stain, polysaccharide

Abstract

Scope:

The PAS stain is used to demonstrate polysaccharides such as glycogen, and mucosubstances such as glycoproteins, glycolipids and mucins in tissues. It is used as a replacement for the H&E in kidney pathology.

Expected Outcome:

Intermyofibrillar Network.....	Pink to Rose
Glycogen.....	Pink to Rose
Myofibrils.....	Unstained
Type I fibers.....	Lighter
Type II fibers.....	Darker
Blood Vessel walls & connective tissue.....	Faintly Stained

Guidelines

1. Leftover matrix will change the pH of the solutions and prevent adequate staining.
2. Bubbles left on tissue will prevent acquisition of high-quality microscopy images.



Materials

Reagents:

1. Commercial kit AbCam 150680
2. Ethanol, Fisher BP2818500
3. Hematoxylin Solution, Mayer's, Sigma MHS32-1L
4. Xylenes, Histological Grade, Sigma 534056

Materials:

1. Easy Dip Slide Staining Jars, Mercedes Medical SIM M90012AS
2. Coplin Dish Staining Dish, Fisher S17495
3. Microscope Cover Slips, Creative Waste Solutions

Solutions:

1. 0.5% Periodic Acid:
Periodic Acid.....5g
Distilled Water....1000mL
Pour out what is needed and reuse for ~1 month
Store at room temperature for 6 months
2. Schiff's Reagent – Commercially Prepared
Richard Allan Scientific Catalog number: 88017
Solution may be reused for ~1 month.
***Check for effectiveness by placing 3 drops of formalin in 2mL Schiff's reagent.
The solution should immediately turn purple. Follow manufacturer expiration date on bottle.
3. Stock of 0.5% Ammonium Hydroxide
2.5mL of Ammonium Hydroxide
497.5mL Milli-Q H₂O
4. Hematoxylin (filtered at least 1x/week)
5. Ethanol solutions
70% - 350mL EtOH + 150mL Milli-Q H₂O
95% - 350mL EtOH + 25mL Milli-Q H₂O

Safety warnings

1. Safety glasses or goggles, proper gloves, and a lab coat required. The area should be adequately vented and a lab mat placed underneath all solutions.
2. Xylenes should be used in the fume hood.

Before start

Allow PAS reagents to come to room temperature for ~30 minutes.



Start with FFPE here:

- 1 Allow PAS "kit" to come to room temperature on the bench.
- 2 For paraffin sections, deparaffinize in xylene, two changes,  00:03:00 each.
- 3 Hydrate through graded alcohols,  00:01:00 each:
100%, 100%, 95%, 70%, water
- 4 Rinse well in distilled water by holding finger over slides, pouring water into sink and adding water. Do this 5 times.

Start Frozen samples here:

5m

- 5 Remove frozen slides from freezer and let equilibrate to room temperature, and then place in 10% Formalin for  00:05:00 . Proceed to step 7.
- 6 If staining is performed following MALDI analysis and samples have matrix on them, remove matrix in 90% ethanol (~2-3 min or until matrix is gone)
 00:03:00 Until Matrix is removed then in 70% ethanol for  00:03:00 .
Proceed to step 7.
- 7 Rinse with distilled water (follow Step 4).
- 8 Place in 0.5% Periodic acid for  00:10:00 (to oxidize).
- 9 Rinse well in distilled water (follow Step 4).

5m



- 10 Pour Schiff's reagent into Coplin jar containing slides. Allow to sit 15-30 minutes at room temperature (kidney samples ~30 minutes).
- 11 Rinse in running warm (~40 °C) distilled or tap water (follow Step 4).

The tissue should appear pink after rinsing with warm water.
- 12 Place slides in movable gray slide holder.
- 13 Counterstain in Hematoxylin (staining line) for  00:01:00 .
- 14 Rinse well in distilled water, starting with blue container next to hematoxylin (follow Step 4).
- 15 Quickly dip slides into "Bluing" agent (0.5% ammonium hydroxide).
- 16 Rinse 1 minute in distilled water until pink color is visible (follow Step 4).
- 17 Dehydrate through graded alcohols, 10 short dips in each:
95%, 95%, 100%, 100%

If the last 100% ethanol rinse is colored by residual stain, additional washing in new 100% ethanol solution should be performed.
- 18 Fix in 2 rounds of xylenes,  00:01:00 each (in the hood).
- 19 Coverslip slides:
 1. Place coverslip on paper towel
 2. Add 2-3 drops of cyto seal to edge (depending on size)
 3. Dip slide into xylenes, take out and roll the xylene lengthwise on slide
 4. Line up slide and cover slip and slowly place the slide on the coverslip
 5. If needed, use dissecting tool to remove bubbles
- 20 Image slides using a Leica brightfield scanner, AxioScan Z1 slide scanner (or equivalent) at 20x (~1 µm spatial resolution) and save as .tiff or .jpg file.

