



May 03, 2024

Patient PBMC collection and cryopreservation and cryorecovery

DOI

dx.doi.org/10.17504/protocols.io.n92ld895ov5b/v1

Rebecca Wallings¹

¹University of Florida



Rebecca Wallings

University of Florida

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ld895ov5b/v1>

Protocol Citation: Rebecca Wallings 2024. Patient PBMC collection and cryopreservation and cryorecovery. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.n92ld895ov5b/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: April 24, 2024



Last Modified: May 03, 2024

Protocol Integer ID: 98735

Keywords: cryorecovery isolation of pbmc, patient pbmc collection, cryorecovery of cell, cryopreservation, cryorecovery isolation, pbmc, cryorecovery, using cpt tube, human blood, cpt tube

Abstract

Isolation of PBMC from human blood using CPT tubes, cryopreservation and cryorecovery of cells

Troubleshooting

PBMC collection and cryopreservation

- 1 Collect 8 mL of blood per BD Vacutainer CPT Cell Preparation Tube with Sodium Citrate (BD Biosciences, 362761).

Invert tubes 8-10 times and centrifuged at room temperature at 1500x *g* for 20 minutes with no brake.

Using a 10mL stripette, transfer the PBMC enriched layer to a new 50-mL conical tube, being sure not to touch the ficoll layer, and add MACS buffer (PBS, 0.5% bovine serum albumin, 20 mM EDTA, pH 7.2) to a final volume of 50 mL, followed by centrifugation at 1800x *g* for 10 minutes at room temperature.

Following removal of the supernatant, resuspend PBMCs in 10mL MACS buffer and count using a hemocytometer with Trypan blue at a 1:20 dilution (10uL of cells mixed with 10uL of 1:20 Trypan blue solution).

Centrifuge PBMCs for 10 min at 300x *g* at 4°C.

Aspirate supernatant and gently resuspend cell pellet in cryopreservation media (RPMI 1640, with 20% FBS and 10% DMSO) at a final concentration of 1×10^7 cells/mL in cryovials (Simport, T311-2).

Place cryovials in a room-temperature Mr. Frosty freezing container with isopropanol as per manufacturer's instructions and store at -80°C overnight.

Remove cryovials from freezing containers and immediately placed into liquid nitrogen for long-term storage.

Cryorecovery

- 2 For cryorecovery, retrieve PBMCs from liquid nitrogen, and thaw at 37°C in water bath.

Once thawed (roughly 1 minute with agitation) slowly add contents of one cryovial to 25mL 37°C filter sterilized complete culture media (RPMI 1640 media, 10% heat-inactivated FBS, 1mM Penicillin-Streptomycin) in 50mL falcon and pellet via centrifugation at 300x *g* for 10 min at room temperature.

Remove supernatant and resuspend cells in 37°C complete culture media for cell counting using Trypan blue.