

Jun 09, 2018

Version 4

PBMC Isolation from apheresis collars with SepMate tubes V.4

DOI

dx.doi.org/10.17504/protocols.io.qpcdviw

Girija Goyal¹

¹Wyss Institute, Harvard University



Girija Goyal

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.qpcdviw>

Protocol Citation: Girija Goyal 2018. PBMC Isolation from apheresis collars with SepMate tubes. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.qpcdviw>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: June 03, 2018

Last Modified: June 09, 2018

Protocol Integer ID: 12740

Keywords: PBMC Isolation , pbmc isolation from apheresis collar, pbmc isolation, apheresis collar, peripheral blood mononuclear cells from whole human blood, peripheral blood mononuclear cell, apheresis pack, peripheral blood, pbmc, mononuclear cell, whole human blood, sepmate tube, isolation, blood

Abstract

Commonly used protocol to isolate peripheral blood mononuclear cells from whole human blood or apheresis packs

Guidelines

Objective: Isolate peripheral blood mononuclear cells from fresh whole blood or apheresis packs, also referred to as Leukopaks or collars. In our case, these are platelet-depleted samples of human blood given from a donor. These can vary in volume and cell composition.

Materials

STEP MATERIALS

⊗ SepMate™-50 (IVD) 100 Tubes **STEMCELL Technologies Inc. Catalog #85450**

⊗ Lymphoprep™ 250 mL **STEMCELL Technologies Inc. Catalog #7801**

⊗ SepMate™-50 (IVD) 100 Tubes **STEMCELL Technologies Inc. Catalog #85450**

⊗ Lymphoprep™ 250 mL **STEMCELL Technologies Inc. Catalog #7801**

Protocol materials

⊗ SepMate™-50 (IVD) 100 Tubes **STEMCELL Technologies Inc. Catalog #85450**

⊗ Lymphoprep™ 250 mL **STEMCELL Technologies Inc. Catalog #7801**

⊗ SepMate™-50 (IVD) 100 Tubes **STEMCELL Technologies Inc. Catalog #85450**

⊗ Lymphoprep™ 250 mL **STEMCELL Technologies Inc. Catalog #7801**


⊗ SepMate™-50 (IVD) 100 Tubes **STEMCELL Technologies Inc. Catalog #85450**

⊗ Lymphoprep™ 250 mL **STEMCELL Technologies Inc. Catalog #7801**

Troubleshooting



Safety warnings

 Any materials that come into contact with blood should be sterilized with 10% bleach before discarding








Before start

- Make sure to repeatedly label sample with donor number, especially if working with multiple donors
- The protocol here is optimized for 10ml of material from platelet apheresis collars. Variations for other sources have been described.



- 1 Cut cone from collar and place on 50mL conical tube. Cut the top end to release blood and allow it to drip into the tube
- 2 Dilute blood collected from apheresis collar with equal volume RPMI or PBS. Mix well.

Isolation of PBMCs

- 3 Use a 50ml SepMate tube and follow manufacturer's directions for density centrifugation
 -  SepMate™-50 (IVD) 100 Tubes **STEMCELL Technologies Inc. Catalog #85450**
 -  Lymphoprep™ 250 mL **STEMCELL Technologies Inc. Catalog #7801**
- 4 Collect the white layer of cells just above the plastic insert into a final volume of 20 ml of media or buffer. Limit the amount of Lymphoprep that is harvested. Mix well.
- 5 Centrifuge at 300 g for 5 minutes.
 -  00:05:00 Centrifugation
- 6 Aspirate media and resuspend cells in 20 mL media/buffer depending on next steps.
 -  20 mL Warm media
- 7 Dilute 100x by adding 10ul cell solutions to 890 ul media/buffer and 100ul of Trypan blue
 -  10 µL Cell solution
 -  890 µL Media/buffer
 -  100 µL Trypan blue

Cell counting

- 8 Count cells using a hemocytometer. Count the number of cells in each of the four quadrants. Use the following formula to find the total number of cells.

$$\text{Cells/ml} = \text{cells counted} * \text{dilution factor} * 10^4 \text{ cells/4}$$
- 9 Cells can be kept in solution in the refrigerator for up to two hours.