




Jun 10, 2018

Version 1

🌐 PBMC isolation from buffy coat V.1

DOI

dx.doi.org/10.17504/protocols.io.quzdw6

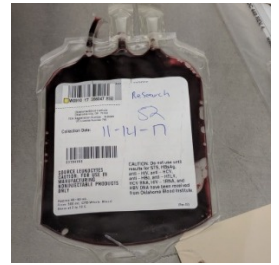

Bulent Arman Aksoy¹, Pinar Aksoy¹, Megan Wyatt¹, Chrystal M. Paulos¹, Jeff Hammerbacher¹

¹Medical University of South Carolina



Bulent Arman Aksoy

Medical University of South Carolina

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.quzdw6>

Protocol Citation: Bulent Arman Aksoy, Pinar Aksoy, Megan Wyatt, Chrystal M. Paulos, Jeff Hammerbacher 2018. PBMC isolation from buffy coat. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.quzdw6>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: June 10, 2018

Last Modified: June 10, 2018

Protocol Integer ID: 12921

Keywords: pbmc isolation from buffy coat, pbmc isolation, buffy coat, isolation

Materials

MATERIALS


✕ Corning™ cellgro™ Lymphocyte Separation Medium **Fisher Scientific Catalog #MT25072CV**

✕ Countess™ Cell Counting Chamber Slides **Catalog #C10314**

Troubleshooting



- 1 Fill 4× 50-ml falcons with 15 ml isolation buffer.
- 2 Cut the buffy coat and fill another 50 ml falcon tube.
- 3 Split the 50 ml blood between 4X 50-ml falcons that have the isolation buffer.
- 4 All tubes should have ~30 ml blood/buffer mix at this point.
- 5 Using a 10 ml pipet, gently underlay 14 ml of Corning™ cellgro™ Lymphocyte Separation Medium below the blood.
- 6 Centrifuge at 800 X g for 25 min at RT with soft deceleration.
- 7 In the meantime, get 4 new 50 ml falcons.
- 8 After the centrifuge, use a 10 ml pipet to transfer the cloudy buffy layer to a fresh tube.
- 9 Wait for 1-2 mins for more of the buffy layer to form and transfer that to the fresh tube as well.

 00:02:00
- 10 Add cold isolation buffer to each tube so that the final volume is now 40 ml.
- 11 Mix by gently inverting tubes for a few times and centrifuge at 500 X g for 10 min at 4°C.
- 12 Discard the supernatant (leave ~5ml liquid) into a bleach filled container.
- 13 Combine the pellet from all tubes into one tube and increase the volume to 40 ml by adding isolation buffer.



Count the cells (1:10 dilution)

- 14 Take 10 ul from the sample and mix with 90 ul of PBS (1:10)
- 15 Mix with 10 ul of Trypan Blue with 10 ul of the sample and load 10 ul to the cell counter chamber for counting.

Equipment

new equipment	NAME
Thermo Fisher Scientific	BRAND
AMQAX1000	SKU
Countess™ II Automated Cell Counter	SPECIFICATIONS

- 16 Continue with the desired T cell isolation kit or freeze PBMC cells.