

Sep 02, 2024

Version 1

PELSA protocol V.1

DOI

dx.doi.org/10.17504/protocols.io.q26g717x1gwz/v1

Kejia Li^{1,2}, Mingliang Ye¹

¹Dalian Institute of Chemical Physics, Chinese Academy of Sciences; ²European Molecular Biology Laboratory



Kejia Li

European Molecular Biology Laboratory

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.q26g717x1gwz/v1>

Protocol Citation: Kejia Li, Mingliang Ye 2024. PELSA protocol. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.q26g717x1gwz/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: September 01, 2024

Last Modified: September 02, 2024

Protocol Integer ID: 106807

Keywords: functional proteomics, drug target discovery, ligand-protein interactions, proteome-wide profiling, pelsa protocol, pelsa method, step procedure of pelsa method, pelsa protocol this protocol detail, unprecedented sensitivity in proteome, affinities of the analyte ligand, analyte ligand, binding protein, proteome, protein

Funders Acknowledgements:

National Key Research and Development Program of China

Grant ID: 2021YFA1302601

the National Natural Science Foundation of China

Grant ID: 92153302

National Natural Science Foundation of China

Grant ID: 22437007

Abstract

This protocol details a step-by-step procedure of PELSA method, a modification-free approach that achieves unprecedented sensitivity in proteome-wide target identification and binding-region determination. Following this procedure, one can determine the binding proteins, binding regions, and binding affinities of the analyte ligand.

Attachments



[PELSA_protocol.pdf](#)

252KB

Troubleshooting

