

Sep 26, 2024

pGEM-T Cloning

 Journal of Virology

DOI

<https://dx.doi.org/10.17504/protocols.io.yxmvme389g3p/v1>

Carolina Lopez¹

¹Washington University



Cecilia Escudero

wustl

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.yxmvme389g3p/v1>

External link: <https://doi.org/10.1128/jvi.00894-25>

Protocol Citation: Carolina Lopez 2024. pGEM-T Cloning. protocols.io
<https://dx.doi.org/10.17504/protocols.io.yxmvme389g3p/v1>

Manuscript citation:

Pye SE, Achouri E, Yang Y, Musa A, López CB Validation of diverse and previously untraceable Sendai virus copyback viral genomes by direct RNA sequencing. Journal of Virology 99(8). doi: [10.1128/jvi.00894-25](https://doi.org/10.1128/jvi.00894-25)



License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: May 29, 2024

Last Modified: September 26, 2024

Protocol Integer ID: 100760

Keywords: cloning, pgem, protocol

Abstract

Protocol for cloning in pGEM-T vector

Materials

Reagents:

- pGEM-T Easy Kit (Promega, A1360)
- Competent E. coli cells (NEB, C3040H)
- LB plates with necessary antibiotic

Before start

PRINCIPLES BEHIND THE PROCEDURE MUST BE UNDERSTOOD. PLEASE CONSULT WITH EXPERIENCED LAB MEMBER THE FIRST TIME YOU USE THIS PROCEDURE. UPDATE AS A GENERAL PROCEDURE AS NECESSARY BUT DO NOT MODIFY WITH SPECIFICS TO YOUR PROJECT, INSTEAD DOWNLOAD AND PASTE A MODIFIED COPY IN YOUR NOTEBOOK. Updated by MH, April 2023



Insert Generation

1h 2m

- 1 1. Perform PCR to generate bands of interest in total volume of  50 μL .

1h 2m

	Volume (μL)
NF H2O	37
Fwd Primer (10 μM)	1
Rev Primer (10 μM)	1
dNTPs	1
Plasmid (20ng)	1
Buffer	5
MgCl ₂	3
Taq Polymerase	1

2. Make 1% low melt-agarose gel.

a) Mix 1 g of Agar with  100 mL of TAE Buffer.

b) Microwave to boil agarose and let cool until you can touch bottle, but gel is not solid.

c) Add  1.5 μL of EtBr to agarose and pour into DNA gel mold with 10 well comb.

d) Let gel solidify.

3. Add  10 μL of loading buffer to PCR reaction then Load  60 μL of reaction into well of gel

4. Run gel for  00:45:00 at 120V.

5. Visualize band with UV light and cut out section of gel with band with new razor blade and place in  1.5 mL tube.

6. Purify Band from gel with Promega Wizard SV Gel and PCR Purification Kit (A9282)

a) <https://www.promega.com/products/nucleic-acid-extraction/clean-up-and-concentration/wizard-sv-gel-and-pcr-clean-up-system/?catNum=A9281>

b) Weigh DNA gel fragment and add  10 μL of Membrane Binding Solution per 10 mg of gel slice.

c) Incubate mixture at  65 $^{\circ}\text{C}$ for  00:10:00 or until gel is completely melted.

d) Add melted gel mixture to SV minicolumn in Collection Tube and incubate at room temperature for  00:01:00 .

e) Centrifuge at max speed for  00:01:00 . Discard flowthrough and reinsert column into tube.



f) Add  700 μL of Membrane Wash Solution. Centrifuge at max speed for

 00:01:00 . Discard flowthrough and reinsert column into tube.

g) Add  500 μL of membrane wash solution. Centrifuge at max speed for

 00:01:00 . Discard flowthrough and reinsert column into tube.

h) Spin empty column for  00:01:00 at max speed to remove excess ethanol.

i) Transfer column to labelled  1.5 mL tube and add  35 μL of NF H_2O .

Incubate for  00:01:00 at room temperature.

j) Centrifuge at max speed for  00:01:00 .

k) Keep eluate and store at  $-20\text{ }^\circ\text{C}$.

pGEM Ligation Protocol

2h 24m

2 1. Set up Ligation Reaction in  0.2 mL PCR tube

2h 24m

Component	Volume (μL)
2x Rapid Ligation Buffer	5
pGEM-T Vector (50ng)	1
PCR Product	X*
H2O	3-X
T4 DNA Ligase	1

a) *Volume of PCR product should be 2 insert:1 pGEM vector molar ratio

b) For example:

	A	B	C	D	E	F	G
	Component	Length of DNA (bp)	Molar ratio	ng of DNA	Volume of 50ng/ μL solution		
	pGEM-T Vector	3015	1	50	1 μL		
	PCR Fragment	500	2	16.58	0.33 μL		
	H2O				2.67 μL		

2. Incubate for  01:00:00 at room temperature

3. Transform Product into E. coli



- a) Add  2 μL of product to  50 μL of TOP10 cells and mix up and down slowly with a pipette.
- b) Incubate for  00:20:00 on ice.
- c) Heat shock bacteria in  42 °C waterbath for  00:01:00 .
- d) Incubate on Ice for  00:03:00 .
- e) Add  100 μL of SOC media and shake in warm room for  01:00:00 .
- f) Plate bacteria onto LB-Antibiotic/Xgal Plate and incubate overnight at  37 °C
- i. Add  40 μL of Xgal to plate and spread with plate spreader. Let dry
 - ii. Then plate  100 μL of bacteria onto plate.
 - iii. Spread cells with plate spreader to get individual colonies
- g) Pick white colonies for miniprep growth and sequencing.

Protocol references

pGEM-T Cloning:

<https://www.promega.com/products/pcr/pcr-cloning/pgem-t-easy-vector-systems/?catNum=A1360>