



Oct 01, 2019

Phage lambda genome concatemerization

DOI

dx.doi.org/10.17504/protocols.io.7k3hkyn

Marijn Ceelen¹

¹Wageningen University

iGEM Wageningen 2019



Marijn Ceelen

Wageningen University

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.7k3hkyn

Protocol Citation: Marijn Ceelen 2019. Phage lambda genome concatemerization. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.7k3hkyn>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: September 23, 2019

Last Modified: October 01, 2019

Protocol Integer ID: 28027

Abstract

This protocol is used to create concatemers of phage genomes. This is done by ligating phage lambda genomes using the overhangs at the cos site with T4 ligase.



- 1 Add the following ingredients in a microcentrifuge tube.

Ingredient	Amount
Ligation buffer 10X	2 μ L
DNA	6 μ g
PEG 4000 (40% W/W)	2.5 μ L
T4 DNA ligase	1 μ L (400 Weiss units)
Water	To 20 μ L

- 2 Incubate at  15 °C overnight