

Mar 26, 2024

## Phenol-chloroform DNA extraction from *Sporosarcina pasteurii*

DOI

<https://dx.doi.org/10.17504/protocols.io.kqdg324eev25/v1>

Michael S. Carter<sup>1</sup>, Matthew J. Tuttle<sup>1,2</sup>, Maneesh K. Gupta<sup>1</sup>

<sup>1</sup>Materials and Manufacturing Directorate, Air Force Research Laboratory, Wright-Patterson Air Force Base, Dayton, Ohio, USA;

<sup>2</sup>Biological and Nanoscale Technologies Division, UES Inc., a BlueHalo Company, Dayton, Ohio, USA



**Matthew J. Tuttle**

Air Force Research Laboratory

### Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.kqdg324eev25/v1>

**Protocol Citation:** Michael S. Carter, Matthew J. Tuttle, Maneesh K. Gupta 2024. Phenol-chloroform DNA extraction from *Sporosarcina pasteurii*. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.kqdg324eev25/v1>

**License:** This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



**Protocol status:** Working

**We use this protocol and it's working**

**Created:** March 25, 2024

**Last Modified:** March 26, 2024

**Protocol Integer ID:** 97344

**Keywords:** DNA extraction, *Sporosarcina pasteurii*, chloroform dna extraction from *sporosarcina pasteurii*, chloroform dna extraction method, chloroform dna extraction, *sporosarcina pasteurii*, dna extraction, extraction of genomic dna, genomic dna, dna, phenol, standard phenol, extraction

## Abstract

This protocol is for extraction of genomic DNA from *Sporosarcina pasteurii*. It is based on a standard phenol-chloroform DNA extraction method.

## Attachments



Phenol-chloroform DN...

179KB

## Guidelines

All steps should be performed in a chemical fume hood.

## Materials

- Brain-heart infusion (BHI) broth supplemented with 330 mM urea
- Resuspension buffer (50 mM Tris pH 8.0, 10% sucrose)
- Lysis buffer (Tris pH 8.0, 10 mg/mL lysozyme)
- 10% sodium dodecyl sulfate (SDS)
- 100 mg/mL RNase A
- 50 mg/mL proteinase K
- 3.0 M sodium acetate, pH 5.5
- 100% ethanol
- TE buffer (10 mM Tris, 1 mM EDTA, pH 8.0)
- Phenol:chloroform:isoamyl alcohol (25:24:1, saturated with 10mM Tris, pH 8.0, 1mM EDTA)
- Chloroform
- 70% ethanol

## DNA extraction



- 1 Grow *Sporosarcina pasteurii* cells in 150 mL of BHI/330 mM urea at 30°C with 200 rpm shaking to an OD600 of 3.5.
- 2 Concentrate cells by centrifugation at 15,000 x g for 10 min and pour off the supernatant.
- 3 Resuspend cells in 10 mL of Resuspension Buffer (50 mM Tris pH 8.0, 10% sucrose).
- 4 Add 2.5 mL Lysis Buffer (Tris pH 8.0, 10 mg/mL lysozyme), 1.5 mL 10% SDS, 5 µL 100 mg/mL RNase A, and 25 µL 50 mg/mL proteinase K to lyse cells and stabilize DNA.
- 5 Incubate for 1 h at 37°C.
- 6 Add 1.3 mL of 3.0 M sodium acetate (pH 5.5) and 30 mL of 100% ethanol to separate DNA from other biomacromolecules by precipitation.
- 7 Spool DNA onto a glass hook and transfer to 20 mL of TE buffer.
- 8 Dissolve DNA in TE buffer for 1 h at 37°C.
- 9 Add a 5 mL aliquot of DNA to 7 mL of phenol:chloroform:isoamyl alcohol (25:24:1, Saturated with 10mM Tris, pH 8.0, 1mM EDTA) and mix by inversion.
- 10 Separate phases by centrifugation at 10,000 x g for 10 min.
- 11 Pipette off the aqueous phase and add to 5 mL of chloroform. Mix by vortexing.
- 12 Separate phases by centrifugation at 10,000 x g for 10 min.
- 13 Pipette off the aqueous phase and add to 30 mL of 100 % ethanol. Mix by inversion.



- 14 Collect precipitated DNA by centrifugation at 10,000 x g for 10 min.
- 15 Wash pelleted DNA twice by adding 5 mL of 70% ethanol, mixing by inversion, centrifuging at 10,000 x g for 5 min, and removing the supernatant.
- 16 Allow the pellet to dry.
- 17 Dissolve the final pellet in 500 µL of TE buffer at 37°C for 1 h.

## Protocol references

Distribution A. Approved for public release: distribution unlimited. AFRL-2024-1482.