

May 19, 2020

Version 2

Phosphate Buffered Saline (PBS) V.2

DOI

<https://dx.doi.org/10.17504/protocols.io.be2hjgb6>

Allen Institute for Brain Science¹

¹Allen Institute

BICCN / BICAN

Allen Institute for Brain S...



Allen Institute

Allen Institute

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.be2hjgb6>

Protocol Citation: Allen Institute for Brain Science 2020. Phosphate Buffered Saline (PBS). **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.be2hjgb6>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: April 14, 2020

Last Modified: March 17, 2021

Protocol Integer ID: 35625

Keywords: RP0001, buffer, ISH, In Situ Hybridization, Immunohistochemistry, IHC, fixation, wash, solution, phosphate buffered saline, neutral wash solution, immunohistochemistry, neutral wash solutions in many process,

Abstract

This protocol is used to prepare Phosphate Buffered Saline (PBS). PBS (-) and PBS (+) which includes Tween are used as biologically neutral wash solutions in many processes including In Situ Hybridization (ISH), Immunohistochemistry (IHC) and Fixation.

Note: Research reported in this publication was supported by the National Institute Of Mental Health of the National Institutes of Health under Award Number U19MH114830. The content is solely the responsibility of the authors and does not necessarily represent the official views of the National Institutes of Health.

Attachments



RP0001_Phosphate_Buf.



39KB

