

Dec 02, 2021 Version 2

Pilot Exercise: Generating the Illumina SampleSheet and sharing data via BaseSpace V.2

DOI

dx.doi.org/10.17504/protocols.io.b2i3qcgn

Ruth Timme¹, Maria Balkey², Tunc Kayikcioglu², Candace Bias², Cameron Boerner², James Pettengill²

¹US Food and Drug Administration; ²US Food And Drug Administration

GenomeTrakr

Tech. support email: genomeTrakr@fda.hhs.gov



Maria Balkey

US Food and Drug Administration

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.b2i3qcgn>

Protocol Citation: Ruth Timme, Maria Balkey, Tunc Kayikcioglu, Candace Bias, Cameron Boerner, James Pettengill 2021. Pilot Exercise: Generating the Illumina SampleSheet and sharing data via BaseSpace. **protocols.io**

<https://dx.doi.org/10.17504/protocols.io.b2i3qcgn> Version created by **Maria Balkey**



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: December 02, 2021

Last Modified: December 02, 2021

Protocol Integer ID: 55611

Keywords: illumina samplesheet, data transfer through basespace, data via basespace, project with fda, samplesheet example, replaced samplesheet example, sharing data, basespace, regarding data transfer, data, fda, pilot exercise, illumina

Abstract

This protocol describes the required steps to generate Illumina SampleSheet and provides guidelines for sharing data via BaseSpace with FDA for the 2021 Pilot Exercise

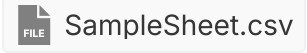
V2:

- added more details regarding data transfer through BaseSpace, including steps for sharing a project with FDA-CFSAN.
- replaced samplesheet example.

Troubleshooting



Generating Sequencing SampleSheet

- 1 Create your SampleSheet using Excel or a text editor. Name your sample sheet according to your internal protocols, and use a *.csv extension. The sample sheet is organized in sections titled Header, Reads, Settings and Data. Section headings are case-sensitive and shown in brackets []. Details for each section are described below.

1.1 Header Section.

A	B
Parameter	Description
Investigator Name	First Name and Last Name for laboratory scientist
Experiment Name	Pilot Exercise SARS-Cov-2
Date	YYYY-MM-DD
Workflow	GenerateFASTQ
Application	FASTQ Only
Instrument Type	MiSeq
Library Prep Kit	Swift Normalase Amplicon NGS Panels (SNAP)
Index Kit	Swift Normalase Amplicon NGS Panels (SNAP)
Description	FDA Wastewater Protocol Pilot Exercise
Chemistry	Amplicon

Table 1: Header Section in Sequencing SampleSheet

1.2 Reads Section

****Enter the length of reads you are generating (could be different from 151)**

A	B
Parameter	Description
Number of cycles in Read 1	151
Number of cycles in Read 2	151

Table 2: Reader Section in Sequencing SampleSheet

1.3 Settings Section

A	B
ReverseComplement	0

Table 3: Settings Section in Sequencing SampleSheet

1.4 Data Section: Include the machine ID in the Sample_Name column. e.g. sample_B01_M01234 where M01234 corresponds to the machine ID. This is the ID that will get populated to the fastq file names.

A	B	C	D	E	F	G	H	I	J
Sample_ID	Sample_Name	Sample_Plate	Sample_Well	I7_Index_ID	index	I5_Index_ID	index 2	Sample_Project	Description
WPP_sample_B.01	WPP_sample_B.01_M01234	1	A1	SU001	TTGTCCTTG	I5_1	CCG AAC AAC A	Lab Abbreviation-WPP	
WPP_sample_C.01	WPP_sample_C.01_WPP-sample_B.01	1	B1	SU002	AGGTGTGTT	I5_2	AAG AATC GGC	Lab Abbreviation-WPP	



	A	B	C	D	E	F	G	H	I	J
	WPP_sample_SA-1.01	WPP_sample_SA-1.01_WPP-sample_B.01	1	C1	SU003	TATTGGTTGG	I5_3	CAC TTC GCT T	Lab Abbreviation-WPP	
	WPP_sample_SA-2.01	WPP_sample_SA-2.01_WPP-sample_B.01	1	D1	SU004	CACTTCGCTT	I5_4	TATTGGT TGG	Lab Abbreviation-WPP	

Table 4: Data Section in Sequencing SampleSheet (dashes converted to underscores for Illumina requirements)

- 1.5 If a new library or sequencing protocol is utilized for processing the pilot exercise samples, make the changes in the corresponding parameters for the sequencing sample sheet.

2 Starting your MiSeq run and connecting to BaseSpace:

When you set up the run on the MiSeq, select the option to log in to BaseSpace.

Transfer sequencing RUN through BaseSpace

- 3 Click the **Runs** tab in the Illumina BaseSpace website .
 - 3.1 Select the **run** that you would like to share with the FDA LFFM wastewater project team.
 - 3.2 Go to the summary tab and click at the share button.
 - 3.3 Enter the email address for the FDA team (**covidtrkr@fda.hhs.gov**), and then click Add Collaborator.
 - 3.4 Click Save Settings. Your run will be automatically shared with the FDA LFFM wastewater project team.

Share this run

210720 M04822 0084 0000000000-JRT3L

Invite a collaborator
covidtrakr@fda.hhs.gov

Optional Message

Add a personal message (optional)

Remaining characters : 140/140

☒ Share the associated Project(s) as well (**Recommended**)
Unchecking this box will limit what the receiptient will have access to.

ADD COLLABORATOR

Collaborators

GenomeTrakr@fda.h...	Pending	Read Only
----------------------	---------	-----------

CANCEL **SAVE SETTINGS**

Figure 1: Sharing data via BaseSpace by Email Invitation

Transfer sequencing PROJECT through BaseSpace

- 4 Click the **Projects** tab in the Illumina BaseSpace website .
- 4.1 Select the **project** that you would like to share with the FDA LFFM wastewater project team.



- 4.2 Click at the share project button.
- 4.3 Enter the email address for the FDA team (**covidtracr@fda.hhs.gov**), and then click Add Collaborator.
- 4.4 Click Save Settings. Your project will be automatically shared with the FDA LFFM wastewater project team.