

Mar 03, 2025

PKH67 Extracellular Vesicle Labeling

DOI

<https://dx.doi.org/10.17504/protocols.io.n2bvj9kmwlk5/v1>

Emily Tabaie¹

¹University of California, Riverside



Emily Tabaie

University of California, Irvine

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n2bvj9kmwlk5/v1>

Protocol Citation: Emily Tabaie 2025. PKH67 Extracellular Vesicle Labeling. protocols.io

<https://dx.doi.org/10.17504/protocols.io.n2bvj9kmwlk5/v1>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: March 03, 2025



Last Modified: March 03, 2025

Protocol Integer ID: 123715

Keywords: pkh67 extracellular vesicle labeling, pkh67 extracellular vesicle labeling this protocol, extracellular vesicles with the lipid dye, extracellular vesicle, lipid dye, pkh67

Abstract

This protocol is designed to visualize extracellular vesicles with the lipid dye, PKH67.



PKH67 Extracellular Vesicle Labeling

- 1 In a microcentrifuge tube add 100 ug (100ul) of EVs in 1 mL of PBS
- 2 Centrifuge at 16,100 g for 10 minutes to form a pellet
- 3 Discard supernatant without disturbing pellet and dissolve the pellet in 100 ul of diluent C and pipette up and down
- 4 In a new microcentrifuge tube add 4 ul of PKH67 to 100 ul of Diluent C and Vortex
- 5 Combine 100 ul of 2X EV with an equal volume of the pKH67 solution
- 6 Pipette the mixture 10 times
- 7 Incubate solution at 37°C for 5 minutes at 300 rpm away from light
- 8 To stop staining add 400 ul of 1% BSA to the mixture and incubate for a minute for excess dye to bind
- 9 Wash stained EVs with PBS 3 times and centrifuge the sample at 16,100 g for 5 minutes
- 10 Discard the supernatant and dissolve the pellet in 100 ul of PBS to get a final concentration of 1 ug/ul
- 11 Two days before transfer endothelial cells to a poly-L-lysine coated coverslip
- 12 Allow the cells to grow a monolayer on the coverslip



- 13 After the monolayer is formed removed the medium carefully and add 1 mL of DMEM to wash the cells twice
- 14 Add 50 ul of PKH67 stained EVs to the coverslip and incubate for 2 hours
- 15 After incubation aspirate the media and wash with PBS 3 times
- 16 Add 1 mL of 4% PFA (made in PBS) to coverslips and fix cells for 15 minutes covered in foil away from light shaking
- 17 Wash coverslips with 1X PBS 3X for 1 min each wash
- 18 Add 1 mL to each well of Perm buffer (Triton X, BSA, PBS) to permeabilize the membrane on the shaker for 10 minutes shaking
- 19 Aspirate the Perm Buffer and wash coverslips with 1X PBS 3X for 1 min each wash
- 20 Block with 1 mL of 5% Donkey Serum (DS) diluted in 1X PBS and incubate on the shaker for 30 min at room temperature shaking
- 21 Aspirate and add 800 ul Primary Antibody Staining in Blocking Buffer and incubate for 1 hour at room temperature shaking
- 22 Wash off unbound antibodies 3X with 1X PBS for 1 min each wash
- 23 Aspirate and add 800 ul of Secondary Antibody Staining in Blocking Buffer and incubate for 1 hour at room temperature shaking covered in foil
- 24 Wash off unbound antibodies with 1X PBS 3 × 10 min each wash shaking
- 25 Grab microscope slides and label them with the experiment, Ab used, initials, and date
- 26 Add (a small) drop of DAPI mounting media for every coverslip, and mount on glass slide



27 Cure slides overnight and image